

Micro Analytical Balances

BMM

Series



BM-20

22 g × 0.001 mg (1 µg)

BM-22 (Smart range)

22 g × 0.01 mg

5.1 g × 0.001 mg (1 µg)

BM-252

250 g × 0.01 mg

BM-200

220 g × 0.1 mg

BM-300

320 g × 0.1 mg

BM-500

520 g × 0.1 mg

Built-in Fanless Ionizer!



...Clearly a Better Value

A&D Company, Limited


How to Eliminate the Greatest Enemy of Precision Weighing

Static electricity, while often overlooked, can be a fatal source of error in precise weight measurements. With the BM Series, A&D seek in various ways to help you perform sometimes-tricky analytical weighing with simplicity and accuracy, including the elimination of static electricity. The world of micro grams (μg) is now accessible to all researchers!

Errors due to static electricity

An electrostatically charged object induces the opposite charge in nearby objects. The attraction between the two charges causes an inaccurate value.


1

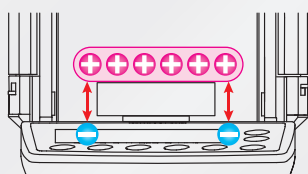

Error when weighing a charged object

The value will be shifted in the positive direction. It then gradually changes as static electricity is discharged into the air or via the weighing pan.

2

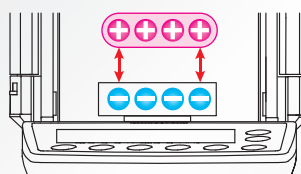

Error when a charged object approaches

The value will be shifted in the negative direction. Static induction from a charged object like clothing (frequent in dry winters) can be a cause of a drifting value.

Built-in Fanless Ionizerⁱ

*i Patent Pending

The internal ionizer neutralizes an electrostatically charged object instantly. Simply hold the sample in the neutralizing chamber for a second before placing it in the weighing chamber.

Direct-current (DC) method

The DC ionizer is fanless so you can even neutralize fine powders without having them blown away.

Long life of approximately 10,000 hours

The discharge electrode pins of the ionizer can be used continuously for approximately 10,000 hours and are safely replaceable.

Please also visit our video library at www.aandd.jp to see how you can easily eliminate static electricity from the sample!

Anti-static Breeze Break

The thin evaporated metal coating of the breeze break blocks outside static electricity.


IBM Series



Neutralizing chamber

Separation plate

Weighing Chamber

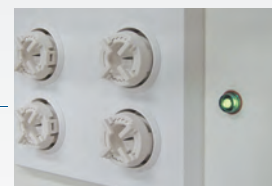

The green LED is lit while the ionizer is on.



Just press the ION key to switch on the ionizer. The "ION" mark blinks on the display.

Motor-driven Internal Calibration Mass

The internal mass of the BM Series allows you to perform calibration quickly and correctly even if you are not skilled at calibrating balances using external masses.

Automatic Self Calibration

The balance can self-calibrate automatically when it detects a change in ambient temperature.

One Touch Automatic Calibration

The accuracy of your balance is assured with one key press.

Automatic Response Adjustment

The balance automatically adjusts its response characteristics (FAST, MID, or SLOW) by analyzing the influence of drafts or vibrations on the stability of the display value. (Manual adjustment is also possible.)

Special Pans and Holder

The BM-20/22 includes three sizes (10 pieces each) of aluminum analytical pans, two pans for weighing PM filtersⁱⁱ, and a microfuge tube holder as standard.

*ii Particulate matter



Aluminum analytical pans

ø8 mm (0.05 ml) / ø12 mm (0.3 ml) / ø15 mm (0.8 ml)



Weighing pans for PM filters

ø50 mm / ø95 mm



Microfuge tube holder

Large Weighing Chamber for Improved Usabilityⁱⁱⁱ

*iii Patent Pending

To block drafts, A&D adopted a double-ring structure around the weighing pan and two vertical shields under the separation plate. This offers more space and easier operation than conventional 1 µg balances with a secondary breeze break in the chamber.



Double-ring structure
(for BM-20/22)

A large flask or cylinder can be weighed by removing the separation plate (for a resolution of 10 µg or higher).



Clear, Reverse Backlit LCD

The contrast of black and white display provides excellent readability.

Interlockable Sliding Doors

It is possible to slide a door open using the handle on the opposite side. Effective use of both hands lets you perform weighing more efficiently.



The "door" sign is displayed until the two side doors are completely closed.)

Built-in Environmental Sensors

The balance displays the temperature, humidity, and atmospheric pressure for users who need to record these environmental conditions.

Twelve Weighing Units

You can easily select from g, mg, oz, ozt, ct, mom, dwt, GN, tael, tola, in addition to PC (counting mode) and % (percent mode).

Density Determination Function

For the BM-252/200/300/500, an optional density determination kit (AD-1653) is available for easy calculations of density.

Diverse Data Management Solutions

Data Memory Function

The balance can store up to 200 sets of weighing data (100 sets with date & time) or 50 sets of calibration history data.

GLP/GMP/GCP/ISO Compliant Output

In accordance with the Good Laboratory Practice (GLP), the balance can output the balance manufacturer, model, serial number, ID number, date, time, calibration mass used, space for signature, etc.

Two Types of Interfaces – RS-232C and Quick USB^{iv} (standard)

Two different devices (e.g. PC and remote controller) can be connected simultaneously to a balance. An external key plug for the optional foot switch (AX-SW128) is also provided as standard.



^{iv}Weighing data transfer from the balance to a PC only. Please use the AX-USB-9P for two-way communication or GLP output via USB.

LAN-Ethernet Interface (BM-08, optional) with WinCT-Plus Software

You can send commands for controlling and acquiring data from multiple balances.



AD-1688 Weighing Data Logger (Standard Accessory)

The AD-1688 is a handheld device (55 × 103 × 16.5 mm) that stores up to 5,000 weighing results with date and time. You can retrieve the saved data later by connecting the AD-1688 to a USB port of your computer (no driver software is necessary).



Save



Carry



and Retrieve the Data!

Specifications

| | BM-20 | BM-22 | BM-252 | BM-200 | BM-300 | BM-500 |
|---|---|------------------------------|----------------------------------|--------------------|-----------------------------|------------------------------------|
| Weighing capacity | 22 g | 5.1 g / 22 g ^v | 250 g | 220 g | 320 g | 520 g |
| Minimum weighing value | 0.001 mg | 0.001 mg / 0.01 mg | 0.01 mg | 0.1 mg | | |
| Repeatability (standard deviation) | 0.0025 mg (for 1 g) | 0.004 mg (for 1 g) / 0.01 mg | 0.03 mg (for 100 g) | 0.1 mg | 0.2 mg | |
| Linearity | ±0.010 mg | ±0.010 mg / ±0.02 mg | ±0.10 mg | ±0.2 mg | ±0.3 mg | ±0.5 mg |
| Stabilization time (when set to FAST) | Approx. 10 secs | Approx. 10 secs / 8 secs | Approx. 8 secs | Approx. 3.5 secs | | |
| Sensitivity drift (10 °C to 30 °C, when automatic self calibration is not used) | ±2 ppm / °C | | | | | |
| Operating temperature | 5 °C to 40 °C (41 °F to 104 °F), 85%RH or less (no condensation) | | | | | |
| Display refresh rate | 5 times / second or 10 times / second | | | | | |
| Display mode | g, mg, oz, ozt, ct, mom, dwt, GN, tael, tola, PC (counting mode), % (percent mode), Specific gravity | | | | | |
| Counting mode Minimum unit mass | 0.1 mg | | | | | |
| Number of samples | 10, 25, 50 or 100 pieces | | | | | |
| Percent mode Minimum 100% reference mass | 10.0 mg | | | | | |
| Minimum 100% display | 0.01%, 0.1%, 1% (depends on the reference mass stored) | | | | | |
| Standard interface | RS-232C, Quick USB | | | | | |
| Internal calibration | Motor-driven internal calibration | | | | | |
| Applicable masses for external calibration | 20 g, 10 g, 5 g, 2 g, 1 g | | 200 g, 100 g 50 g, 20 g, 10 g | 200 g, 100 g, 50 g | 300 g, 200 g 100 g, 50 g | 500 g, 300 g 200 g, 100 g, 50 g |
| Discharge electrode pins of the ionizer | Life approx. 10,000 hours (continuous use) | | | | | |
| Weighing pan (diameter) | 25 mm (50 mm, 95 mm for weighing filters) | | 90 mm | | | |
| Power consumption | Approx. 30 VA (supplied to the AC adaptor) | | | | | |
| External dimensions | 259 (W) × 466 (D) × 326 (H) mm | | | | | |
| Net weight | Approx. 10 kg | | | | | |
| Standard accessories | AC adapter, Instruction manual, WinCT (CD-ROM), Weighing data logger (AD-1688), Tweezers for calibration mass (AD-1689) Tweezers and two weighing pans for PM filters ^{vi} ; Microfuge tube holder ^{vi} ; Three sizes of aluminum analytical pans (10 pcs each) ^{vi} | | | | | |

* v : Smart range function: The minimum weighing value will switch to 0.01 mg automatically when the display value exceeds 5.1 g but returns to 0.001 mg by pressing the RE-ZERO (tare) key.

* vi : For BM-20/22 only.

Options

BM-08^{vii} Ethernet interface with WinCT-Plus software

BM-014 Pipette accuracy testing kit


* vii : BM-08 and the standard Quick USB interface cannot be used at the same time.

External Dimensions (mm/inches)

| Pan size | |
|--------------------|------------|
| Model | A |
| BM-20/22 | Ø25 / 0.98 |
| BM-252/200/300/500 | Ø90 / 3.54 |

Accessories

| | |
|------------------------|--|
| AD-1653 | Density determination kit (for BM-252/200/300/500 only) |
| AD-1671 | Anti-vibration table for balances |
| AD-1672 | Tabletop breeze break |
| AD-1682 | Rechargeable battery unit |
| AD-1684 | Electrostatic fieldmeter |
| AD-1687 | Weighing environment logger |
| AD-1688 | Weighing data logger |
| AD-1689 | Tweezers for calibration mass |
| AD-8121B | Compact printer |
| AD-8526 | Serial/Ethernet converter |
| AD-8920A | Remote display |
| AD-8922A | Remote controller |
| AX-BM-031 | Display covers (5 pcs) |
| AX-BM-032 | Microfuge tube holder (for BM-20/22 only) |
| AX-BM-NEEDLESET | Electrode units |
| AX-KO2466-200 | RS-232C cable (9P-9P, 2 m) |
| AX-ROUND-PAN-L | Aluminum analytical pans (Ø15 mm, 100 pcs) |
| AX-ROUND-PAN-M | Aluminum analytical pans (Ø12 mm, 100 pcs) |
| AX-ROUND-PAN-S | Aluminum analytical pans (Ø8 mm, 100 pcs) |
| AX-SW128 | Foot switch |
| AX-T-314A-S | External key plug |
| AX-USB-9P | Serial/USB converter |

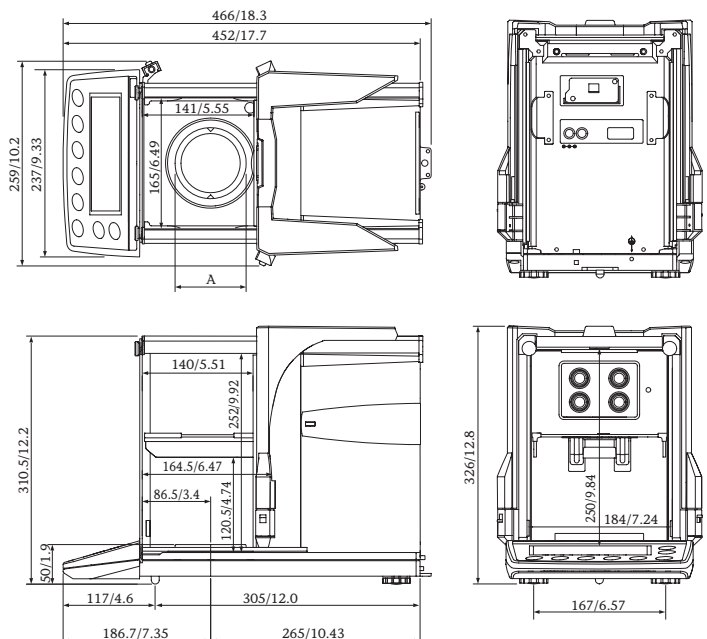



WolfLabs

Pricing on any accessories shown can be found by keying the part number into the search box on our website.

The specifications listed in this brochure are subject to change by the manufacturer and therefore cannot be guaranteed to be correct. If there are aspects of the specification that must be guaranteed, please provide these to our sales team so that details can be confirmed.

www.wolflabs.co.uk

Tel : 01759 301142

Fax : 01759 301143

sales@wolflabs.co.uk

Please contact us if this literature doesn't answer all your questions.