Vortamix Mini Vortexer

Operator's Manual
Manuel de l'opérateur
Bedienungshandbuch
Manual del operador
Manuale d'uso
オペレーターマニュアル

ARGOS
TECHNOLOGIES *manufacturing solutions for science*

Argos Technologies, Inc.
625 Bunker Court
Vernon Hills, IL 60061

Toll Free: 800-886-8675
International: 847-622-0456
www.argos-tech.com

CE  **RoHS**
2011/65/EU

Introduction

Thank you for your purchase of the Argos Vortamix Mini Vortexer. This unit has been designed to rapidly vortex centrifuge tubes with one hand. The small, compact base requires little bench space, yet is stable enough to remain stationary while in use.

Note: Please be sure to read the safety and operating instructions before use. Also, when performing the vortexing function, observe all laboratory safety guidelines including the use of protective clothing, safety goggles and gloves when handling biological samples or chemicals.

Package Contents

Upon receipt of the Vortamix Mini Vortexer, please carefully unpack your unit to ensure it is complete and in good working order. Should you find the unit damaged upon inspection, notify your freight carrier at the time of receipt as well as your distributor.

Should any parts listed below be missing or you discover concealed damage, contact Argos Technologies immediately for resolution (see page 4 for contact information).

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Materials and Specifications

Model #	V0000 (US)	V1000 (International)
Speed	3,000rpm	
Orbital Diameter	Ø4.5mm	
Capacity	Max. 80mL	
Tube Diameter	Max. Ø30mm	
Ambient	5-40°C, ≤80%RH	
Dimensions	Ø95 x 85mm	
Weight	0.55kg	
Power Supply	AC 100-240V 50/60Hz	
Power	5 W	
IP rating (DIN EN 60529)	IP40	
Working Gasket Color	Red	

Warranty Information

Vortamix Mini Vortexer 2 Year Warranty

Argos Vortamix Mini Vortexer is warranted to be free of defects in materials and workmanship for a period of 24 months from the date of purchase, providing that:

- The unit has not been opened or attempted to be opened by the end user.
- The user ensures that it has been carefully cleaned and decontaminated.
- The unit has been operated according to the instructions and not in any improper way.

In order to qualify for warranty repair; the customer must return the warranty card supplied with the unit that includes all of the information required to validate proof of purchase. Should warranty service be required, repair or replacement will be provided by Argos without charge. Freight shipping the unit to Argos will be the responsibility of the purchaser. There are no other warranties or guarantees, implied or otherwise expressed including any warranties of application or merchantability.

Repair Service

Should your Vortamix Mini Vortexer require service or repair, please contact Argos Technologies customer service. They will issue a Return Authorization Number and advise on the proper packaging and method of shipment.

Customer Service: in US: 800-886-8675 outside US: 847-622-0456

Argos Technologies, Inc.

625 Bunker Court
Vernon Hills, IL 60061

Be sure to list your RA number on the shipping label.

Please use the same numbers listed above, should you require any technical assistance or support.

Getting Started

CAUTION

English

- Only use parts and accessories approved to work with the Vortamix Mini Vortexer.
- All power cords are not the same! Only use the power cord adapter supplied with the unit.
- Make sure you cover your tubes prior to vortexing.
- Do not vortex flammable or hazardous liquids.
- The Vortamix Mini Vortexer will respond to light pressure on the top gasket. Avoid applying excess pressure to the tube holder.

Operating Instructions

- Place the vortexer on a flat, sturdy surface to avoid excessive vibration.
- Connect the power supply to the outlet socket.
- Place the tube on the gasket cap and press down lightly to begin vortexing action.
- The unit will automatically stop once the tube has been removed from the gasket cap.

Maintenance, Care and Safety Precautions

WARNING!!

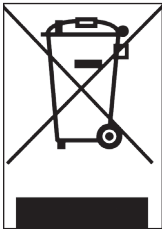
Before cleaning disconnect the Vortamix Mini Vortexer from the power supply. Never clean the power supply or immerse the vortexer in water.

Safety Precautions

- Never use with highly flammable or corrosive liquids.
- Ensure the current of your electrical source matches the voltage of the charger unit. Only use the adapter supplied with the unit.
- Do not open the Vortamix Mini Vortexer as it will void warranty.

Cleaning & Care

- Clean the unit with a soft cloth and a non-corrosive liquid.
- Do not spray the unit with any liquids.
- Make sure the unit is completely dry before use.

This equipment is labeled with the crossed out wheeled bin symbol to indicate that the equipment must not be disposed of with unsorted waste. It is your responsibility to correctly dispose of the equipment by handing it over to an authorized facility for separate collection and recycling. It is also your responsibility to decontaminate the equipment in case of biological, chemical or radioactive contamination. So as to protect from health hazards the persons involved in the disposal or recycling of products.

For more information about where you can drop off your waste equipment, please contact your local dealer from whom you originally purchased the product.

By doing so, you will help to conserve our natural resources and you will ensure that your waste products are recycled in a matter that protects human health and protects the environment.

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