

CAPP's **EXPell Pipette Tips Reload System** is available for 10µL standard and extended length as well as for 200µL and 1000-1250µL tips. Depending on tip volume, it consists of towers with 5, 8 or 10 insert plates containing 96 tips each.

The system is extremely easy to use. Take off the tower cover and pick up the insert plate with tips. Place it into the empty rack. The rack has been refilled with 96 tips.

The **EXPell Reload System** enables multiple savings: Shipping costs are significantly reduced by up to 75% of the packing volume. The tips in an **EXPell Reload System** are packed more compactly compared to standard racked or bulk tips.


In the **EXPell Reload System** the tips from the upper insert plate are partially placed into the tips of the lower plate, resulting in a slim tower, taking much less space than a single box with 10 racked tips.

CAPP's **EXPell Reload System** reduces the plastic waste. Each rack can be reused multiple times and, apart from the insert plates with tips, there is no further plastic component in the **EXPell Reload System** – only the recyclable cardboard packaging.

Further saving with **EXPell Reload System** concerns the space. Thanks to the minimal footprint, the towers of the **EXPell Reload System** take up as less space on the benchtop and in the storage room as possible.

The system is compatible with all **CAPP** racks of corresponding tip volumes.

Spare, empty racks are available as accessories.

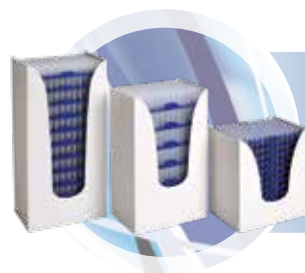


1 Open the refill by sliding it towards the top.

2 Prepare an empty rack. Remove its original grid.



3 Pick up the first grid of refill tips and insert into the rack. The racks can be reused.

4 Once the tips are properly installed in the rack, you may close the hinged lid.

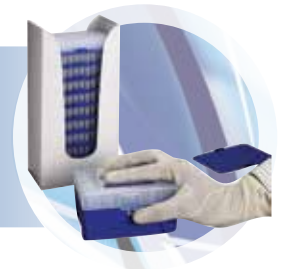

Ordering Information

Cat. No.	Description
5030035C	Expell 10µL, reload tips, 10x10x96 pcs.
5030034C	Empty rack for 10µL, reload tips, 50 pcs.
5030055C	Expell 10/20µL XL, extra narrow, 10x10x96 pcs.
5030054C	Empty Rack ro 10/20µL XL, extra narrow, reload tips, 50 pcs.
5030095C	Expell 200µL, reload tips, 10x10x96 pcs.
5130111C	Expell 300µL, reload tips, 10x8x96 pcs
5030094C	Empty rack for 200µL/300µl, reload tips, 50 pcs.
5130155C	Expell 1000µL (1250µL), reload tips, 10x5x96 pcs.
5130125C	Empty rack for 1000µL (1250µL), reload tips, 32 pcs.



WolfLabs

Pricing on any accessories shown can be found by keying the part number into the search box on our website.

The specifications listed in this brochure are subject to change by the manufacturer and therefore cannot be guaranteed to be correct. If there are aspects of the specification that must be guaranteed, please provide these to our sales team so that details can be confirmed.

www.wolflabs.co.uk

Tel : 01759 301142

Fax : 01759 301143

sales@wolflabs.co.uk

Please contact us if this literature doesn't answer all your questions.