

## ClearLine Mini Vortex Mixer



### Introduction

The ClearLine Mini Vortex Mixer is a compact yet powerful vortex mixer designed to accommodate tubes of varying sizes, from 0.4ml to 50ml. Dynamic balance technology provides smooth speed control, quiet operation and minimal excess vibration.

Please read this manual thoroughly before attempting to operate the ClearLine Vortex Mixer.

### Specifications

Speed	Fixed, 2800 rpm
Operating mode	TOUCH mix
Head	Rubber cup for single tube mixing
Orbit	4mm
Operating environment	+4°C to 65°C
Dimensions (WxDxH)	9.1x9.9x6.6cm/3.7x3.9x2.6"
Weight	0.4kg/1lb
Electrical	Universal, 100 to 240V, 50-60 Hz
Warranty	5 years, parts and labor

### Product Set-Up

Remove the vortexer from the box. Keep all packing materials until the unit is shown to be in good working order. Check that the appropriate power cord has been received. If the included power cord does not match the local requirements, contact your ClearLine dealer.

Place the vortexer on a clean, stable surface and connect the power cord to the appropriate electrical outlet.



# WolfLabs

**Pricing on any accessories shown can be found by keying the part number into the search box on our website.**

The specifications listed in this brochure are subject to change by the manufacturer and therefore cannot be guaranteed to be correct. If there are aspects of the specification that must be guaranteed, please provide these to our sales team so that details can be confirmed.

**[www.wolflabs.co.uk](http://www.wolflabs.co.uk)**

**Tel : 01759 301142**

**Fax : 01759 301143**

**[sales@wolflabs.co.uk](mailto:sales@wolflabs.co.uk)**

Please contact us if this literature doesn't answer all your questions.