



Airstream®

*Airstream Class III, Biological Safety Cabinet,
Model: AC3-4B_*

Class III, Biological Safety Cabinets The Premier Solution for High Containment Laboratories

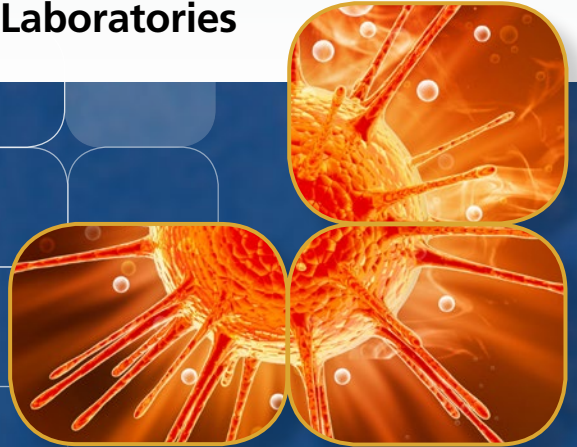

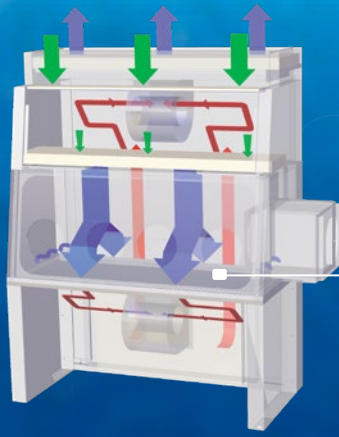

Cabinet Filtration System

Dynamic air barrier, inflow and forward-directed downflow air converge

- Ambient air is pulled into the cabinet through a disposable pre-filter, trapping larger particles and prolonging the life of main filter. Air is then passed through the downflow ULPA filter into the work zone as a vertical laminar flow air stream, bathing the work surface in clean air.
 - The uniform non-turbulent air stream protects against cross contamination within and throughout the work area.
 - Near the work surface, the downflow air stream splits and exits the chamber through perforations around the perimeter of the work zone.
- Air from work zone passes through the first ULPA exhaust filter below the work surface, before traveling through internal ducting and exhausting to the laboratory environment after passing through the second ULPA exhaust filter located above the cabinet's main chamber.

- ULPA-filtered air
- Unfiltered / potentially contaminated air
- Room air / inflow air



Sentinel™ Silver Microprocessor Controller

- Supervises all cabinet functions
- Easy-to-use soft keys for menu, blower, light, UV, electrical outlet, etc
- Large display which monitors the operational parameters



Magnehelic* pressure gauge

- Mounted in the rear of the work zone for at-a-glance monitoring of work zone negative pressure

*Registered trademark of Dwyer Instruments, Inc.



Pass Box

- Equipped with interlocking doors that permits materials transfer without risk of environmental contamination



Glove ports

- For easy and safe glove changing




Neoprene™ Synthetic Arm-length Gauntlets

- Single piece, leak-tested assembly guarantees maximum protection and flame and abrasion resistant

Glove Sizes

- To find out your glove size, measure (in inches) around your hand with a tape measure at the place indicated on the diagram. This measurement, in inches, is the closest to your glove size, i.e. 8" is equal to a glove size of 8.
- You should use your dominant hand, the right if you are righthanded, and the left if you are left-handed.
- Standard glove sizes available are 7, 8 (standard size) and 9. Specify when ordering.

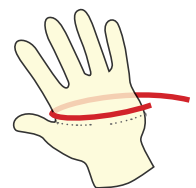

Available in 1.2, 1.5, and 1.8 meter models (4', 5', and 6'). Shown with optional telescoping stand.

ISOCIDE™

ULPA filter

CE
EN12469


| | Biosafety Cabinets | For Air Quality | For Filtration | For Electrical Safety |
|----------------------|--------------------|--|---|--|
| Standards Compliance | EN 12469, Europe | ISO 14644.1 Class 3, Worldwide IEST-G-CC1001, USA IEST-G-CC1002, USA | IEST-RP-CC034.1, Worldwide IEST-RP-CC007.1, Worldwide IEST-RP-CC001.3, Worldwide EN 1822 (H14), Europe | IEC 61010-1, Worldwide EN 61010-1, Europe UL 61010-1, USA CAN/CSA-22.2, No. 61010-1 |

Air-tight Built

- Completely welded construction to ensure absolute containment for operator and environmental safety

Dual Exhaust ULPA FILTER

- 10x Filtration efficiency of HEPA filter
- Creates ISO Class 3 work zone instead of industry-standard ISO Class 5

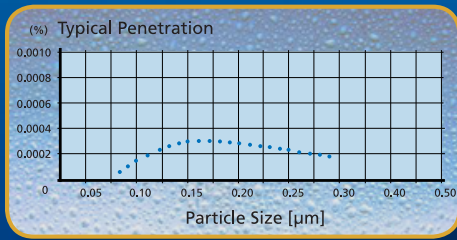

Angled Front

- To optimize user comfort, reduce glare and maximize reach into the work area

ISOCIDE™ Powder Coat

- Silver-ion impregnated powder coat
- Inhibit microbial growth to improve safety



General Specifications, Airstream Class III, Biological Safety Cabinets

Note to customer: Insert electrical voltage number into last model number digits _ when ordering

| Model | AC3-4B_ | AC3-5B_ | AC3-6B_ | |
|---|---|--|---|---------|
| Nominal Size | 1.2 meter (4') | 1.5 meter (5') | 1.8 meter (6') | |
| External Dimensions (W x D x H) | 1665 x 850 x 2250 mm 65.6" x 33.5" x 88.6" | 1970 x 850 x 2250 mm 77.6" x 33.5" x 88.6" | 2275 x 850 x 2250 mm 89.6" x 33.5" x 88.6" | |
| Internal Work Area, Dimensions (W x D x H) | 1340 x 560 x 650 mm 52.8" x 22.0" x 25.6" | 1645 x 560 x 650 mm 64.8" x 22.0" x 25.6" | 1950 x 560 x 650 mm 76.8" x 22.0" x 25.6" | |
| Internal Work Area, Space | 0.75 m ² (8.1 sq.ft) | 0.92 m ² (9.9 sq.ft) | 1.09 m ² (11.7 sq.ft) | |
| Number of Gloves Ports | 2 ports | 4 ports | 4 ports | |
| Glove Type & Sizes Available | Neoprene™ polychloroprene synthetic rubber gauntlets. Available in sizes 7, 8 (standard size) and 9 | | | |
| Initial Airflow Volume | 603 m ³ /h (355 cfm) | 756 m ³ /h (445 cfm) | 902 m ³ /h (531 cfm) | |
| Negative Work Zone Pressure | -275 Pa (-1.1" Wg) | | | |
| Pre-Filter | Disposable and non-washable polyester fibers with 85% arrestance / EU3 rated | | | |
| ULPA Filter Typical Efficiency (Downflow, 1st Exhaust, 2nd Exhaust) | Typical: 99.999% at 0.1 to 0.3 µm and MPPS | | | |
| Sound Emission (Typical)* | NSF 49 | <54 dBA | <55 dBA | |
| | EN 12469 | <51 dBA | <52 dBA | |
| Fluorescent Light Intensity At Zero Ambient | >2000 Lux (>186 foot candles) | >1800 Lux (>167 foot candles) | >2000 Lux (>186 foot candles) | |
| Cabinet Construction | Main Body | 1.5 mm (0.06") 16 gauge electro-galvanized steel with white oven-baked epoxy-polyester ISOCIDE™ antimicrobial powder coated finish | | |
| | Work Zone | 1.2 mm (0.05") 18 gauge stainless steel, type 304, with 4B finish | | |
| Electrical** | 220-240V, AC, 50Hz, 1Ø | AC3-4B1 | AC3-5B1 | AC3-6B1 |
| | Cabinet Full Load Amps (FLA) | 3 A | 3 A | 3 A |
| | Optional Outlets FLA | 5 A | 5 A | 5 A |
| | Cabinet Nominal Power | 361 W | 430 W | 455 W |
| | Cabinet BTU | 1232 | 1467 | 1553 |
| | 110-120V, AC, 60Hz, 1Ø | AC3-4B2 | AC3-5B2 | AC3-6B2 |
| | Cabinet Full Load Amps (FLA) | 10 A | 11.5 A | 11.5 A |
| | Optional Outlets FLA | 5 A | 5 A | 5 A |
| | Cabinet Nominal Power | 536 W | 586 W | 620.5 W |
| | Cabinet BTU | 1829 | 2000 | 2117 |
| Net Weight | 498 kg (1096 lbs) | 598 kg (1316 lbs) | 676 kg (1487 lbs) | |
| Shipping Weight | 606.5 kg / 1337 lbs | 615 kg / 1356 lbs | 720 kg / 1587 lbs | |
| Shipping Dimensions, Maximum (W x D x H) | 2600 x 1950 x 1320 mm 102.4" x 76.8" x 52.0" | 2600 x 2150 x 1320 mm 102.4" x 84.6" x 52.0" | 2600 x 2150 x 1320 mm 102.4" x 84.6" x 52.0" | |
| Shipping Volume, Maximum | 6.69 m ³ (236 cu.ft.) | 7.38 m ³ (261 cu.ft.) | 7.38 m ³ (261 cu.ft.) | |

Accessories for AC3 Biological Safety Cabinets


| Cabinet | Stainless Steel Side Wall | AC3-4B1 (2010204) | AC3-5B1 (2010207) | AC3-6B1 (2010210) |
|-------------------|---------------------------|---------------------|---------------------|---------------------|
| Exhaust Ducting | Anti-blowback Valve | ABBV-10P (5170352) | | |
| | Exhaust Collar 10' dm | ECO-AC34B (5170086) | ECO-AC35B (5170088) | ECO-AC36B (5170139) |
| | Air Tight Damper | B2-Damper (5170104) | | |
| Work Zone | Ultraviolet (UV) Lamp | UV-15A (5170251) | UV-30A (5170255) | |
| Electrical Outlet | Direct Mounted | EO-HD (5170036) | | |

* Noise reading in open field condition/ anechoic chamber.

** Additional voltages may be available; contact Esco for ordering information.

ESCO GLOBAL NETWORK

43 LOCATIONS IN 23 COUNTRIES ALL OVER THE WORLD

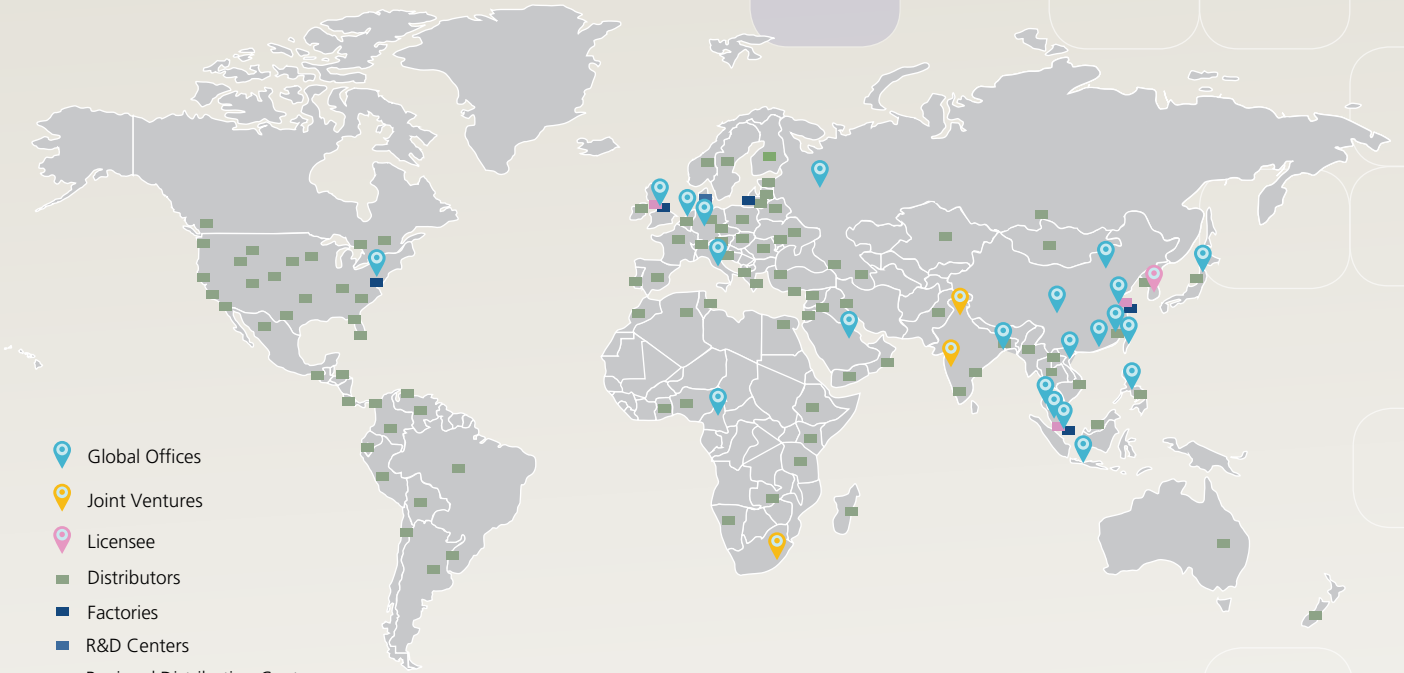

- ART Equipment
- Biological Safety Cabinets
- CO₂ Incubators
- Compounding Pharmacy Equipment
- Containment / Pharma Products
- Ductless Fume Hoods
- Lab Animal Research Products
- Laboratory Centrifuges
- Laboratory Fume Hoods
- Laboratory Ovens and Incubators
- Laboratory Shakers
- Laminar Flow Clean Benches
- PCR Cabinets
- PCR Thermal Cyclers
- Powder Weighing Balance Enclosures
- Ultra-low Temperature Freezers

The Esco Group of Companies is a global life sciences tools provider with sales in over 100 countries. The group is active in lab equipment, pharma equipment and medical devices. Manufacturing facilities are located in Asia and Europe. R&D is conducted worldwide spanning the US, Europe and Asia. Sales, service and marketing subsidiaries are located in 12 major markets including the US, UK, Singapore, Japan, China and India. Regional distribution centers are located in the US, UK, and Singapore.

Life Science • Chemical Research • Assisted Reproductive Technology (ART) • Pharmaceutical Equipment • General Equipment

ESCO

WORLD CLASS. WORLDWIDE.



WolfLabs

Pricing on any accessories shown can be found by keying the part number into the search box on our website.

The specifications listed in this brochure are subject to change by the manufacturer and therefore cannot be guaranteed to be correct. If there are aspects of the specification that must be guaranteed, please provide these to our sales team so that details can be confirmed.

www.wolflabs.co.uk

Tel : 01759 301142

Fax : 01759 301143

sales@wolflabs.co.uk

Please contact us if this literature doesn't answer all your questions.