XR600 & 1300 Cabinets XR2 & 3 Counters with FD2-10 Controller & LCD55 Display

English

Français

Deutsch

Nederlands

A Division of ITW Ltd
Foster Refrigerator
Oldmedow Road
King's Lynn
Norfolk, PE30 4JU
United Kingdom

00-570033 January 2015 Issue 2

Call: +44 (0)843 216 8800

Fax: +44 (0)843 216 4700
www.fosterrefrigerator.co.uk

Email: support@foster-uk.com

Important Note to the Installer:

Please ensure that this document is passed to the user as it contains important guidelines on operation, loading, cleaning and general maintenance and should be kept for reference.

General Electrical Safety

Foster Refrigerator recommends that the equipment is electrically connected via a Residual Current Device; such as a Residual Current Circuit Breaker (RCCB) type socket, or through a Residual Current Circuit Breaker with Overload Protection (RCBO) supplied circuit.

Disposal Requirements

If not disposed of properly all refrigerators have components that can be harmful to the environment. All old refrigerators must be disposed of by appropriately registered and licensed waste contractors, and in accordance with national laws and regulations.

Start-Up and Test Sequence

After unpacking, clean and allow the cabinet to stand for 2 hours before turning on (cleaning directions supplied within this manual). Ensure the cabinet is situated away from both hot and cold air sources, as this will affect its performance. Make sure that a minimum clearance of 150mm above the cabinet is available for ventilation and effective operation. Connect the unit to a suitable mains power outlet and turn the supply on. Do not plug or unplug the unit with wet hands.

Cabinets & counters are supplied ready for operation.

After connecting the unit to the mains the display will briefly show a dash in the centre of the screen. This will then show *OFF*

Press and hold the  button for 3 seconds then release. (If pressed and held for 5 seconds then released will start the 'Test Sequence').

To cancel the 'Test Sequence' press and release the  button during the test.

If not pressed the test will continue and when complete the controller will wait 1 minute, then resume normal operation.

User Adjustments

To Check Storage Temperature Set Point

Press the  button for at least half a second to display the set point.

Manual Defrost

To initiate a manual defrost press and hold the defrost button for 2 seconds 

Set Unit to Standby

When pressing the  button for 3 seconds, the unit will enter the standby mode and display *OFF*

This indication will be displayed while the unit is not operating but with mains power applied to the unit. This mode may be used for interval cleaning regimes and short periods when the unit is not required.

For extended periods of inactivity the mains supply should be isolated.

Alarms and Warnings

During normal operation the display will show either the temperature or one of the following indicators:

	Room High Temperature Alarm		Air Temperature Probe T1 Failure
	Room Low Temperature Alarm		Evaporator Temperature Probe T2 Failure (Freezer Cabinets Only)
	Door Open Alarm		

Shelf Supports

The unit has slotted removable shelf supports allowing flexible shelf positioning. Ensure the trayslides/clips are fitted parallel and level.

When loading product onto the shelves ensure that it does not overflow the rear and sides as this will restrict airflow throughout the cabinet.

Lighting

Where fitted the light can be turned on or off by pressing and releasing the  button.

Load/Product Distribution and Airflow

Product should be evenly distributed throughout the cabinet or counter.

For both cabinets and counters, no more than 40kg should be placed on each shelf, basket or drawer. Product should be placed in a way that ensures air can circulate around/through it.

Cleaning and Maintenance

Important: Before cleaning, the unit should be put into standby and then the power supply should be turned off at the mains. Please do not plug or unplug the unit with wet hands. Only when cleaning has been completed and the unit is dry should the cabinet be turned back on at the mains. Using high pressure water jets or lances is prohibited; the warranty will not cover problems that result from failure to heed the above warning.

Regular Maintenance:

- > As and when required remove all product and shelving (beware of cold and or sharp components) from the unit. Clean exterior and interior surfaces with mild liquid detergent, following the directions on the pack at all times. Rinse surfaces with a damp cloth containing clean water. Never use wire wool, scouring pads/powders or high alkaline cleaning agents i.e bleaches, acids and chlorines as these may cause damage. Dry the unit thoroughly before re-using.
- > Cleaning the condenser at least once a month (it is located behind the unitcover); beware of hot and possible sharp components.
On both the cabinet and counter remove the unit cover securing screw(s) positioned underneath the front of the unitcover. Pull the cover forward and it will release from the securing pegs to allow access to the condenser coil.
Once the cover is removed please do not leave the unit unattended. Use a soft brush or vacuum cleaner to remove dirt/dust. If there are any stubborn grease deposits left on or through the condenser call your supplier to carry out a full service (this is normally chargeable). Failure to maintain the condenser may invalidate the warranty of the condensing unit and cause premature failure of the motor/compressor. **DO NOT** use a wire brush to clean the condenser. Re-fit any panels removed after cleaning.
- > Door gaskets should be inspected on a regular basis and replaced if damaged. To clean, wipe with a warm damp soapy cloth followed by a clean damp cloth. Finally thoroughly dry before closing the door.
- > Shelves, baskets and their supports should be removed to clean. The shelves are dishwasher safe; however the supports should be cleaned with warm soapy water then rinsed and dried.

Before calling your supplier please make sure that:

- a) The plug has not come out of the socket and the mains power supply is on i.e. is the controller illuminated?
- b) Check to see if the unit is in standby
- c) The fuse has not blown
- d) The cabinet is positioned correctly – cold or warm air sources are not affecting the performance
- e) The condenser is not blocked or dirty
- f) The products are placed in the unit correctly
- g) Defrost is not in progress or required

If the reason for the malfunction cannot be identified, disconnect the electrical supply to the unit and contact your supplier. When requesting a service call, please quote the model and serial number which can be found on the silver label located on the inside of the unit (starts E.....).

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