Rota-Filler 3000[®]

Instruction Manual	2
Manuel d'instructions	17
Manual de Instrucciones ..	32
Bedienungsanleitung	47

U.S. Pat. No. D510,629
Euro. Pat. App. 1 555 067
Com. Design No. 000203799-0001
China Design No. ZL20430066669.8
Other Patents Pending

Heathrow Scientific[®]

HEATHROW SCIENTIFIC, LLC
620 Lakeview Parkway • Vernon Hills, IL 60061 • USA
Phone: +1-847-816-5070 • Fax: +1- 847-816-5072

Email: info@heathrowscientific.com
www.heathrowscientific.com

Contents

Safety Precautions	3
Product Description	4
Components	5
Technical Data	6
Getting Started	
Charging the batteries.....	7
Mounting the holder	8
Using the stand.....	8
Pipetting	
Adjusting the nozzle	9
Attaching the pipette.....	9
Filling the pipette	9
Dispensing	10
Maintenance	
Cleaning the pipette adapter and replacing the filter.....	11
Checking for leaks.....	11
Recharging the batteries.....	12
Replacing the batteries	13
Trouble Shooting	14
Ordering Information	14
Return for Repair.....	15
Warranty.....	15

Safety Precautions

The Rota-Filler 3000[®] may be involved with hazardous materials. This instruction manual can not address all safety hazards. It is the responsibility of the user to consult and observe all health and safety precautions and to assess the instrument's suitability to the task.

1. Before operating the instrument, read this instruction manual.
2. Observe general laboratory safety precautions and regulations.
3. Follow all precautions from the Material Safety Data Sheet (MSDS) for any reagent you use with this instrument.
4. Never operate the instrument in a hazardous or flammable environment.
5. Do not pipette flammable solvents such as Acetone or Ether.
6. Always work in a manner which neither endangers the user, nor any other person.
7. Use the instrument only for pipetting, and within the recommended limits of its chemical resistance and mechanical properties.
8. Never use force when using this instrument.
9. When not in use, store the instrument on stand or in holder.
10. Should the instrument fail to work to specification, immediately stop pipetting. Clean and troubleshoot the instrument according to the instructions under "Trouble Shooting" (page 14) before any further use of the instrument. Contact an authorized dealer or the manufacturer if trouble persists.
11. Only use original manufacturer's accessories (AC adapter, stand, and holder) and spare parts (filters, nose piece, and pipette adapter).
12. Only original recharger can be used for recharging the NiMH batteries.
13. The recharger may only be used together with this instrument.
14. Do not immerse the recharger in any liquid nor use it with wet hands.
15. Repairs are to be performed by trained and authorized service personnel only.
16. Opening the instrument or improper use of the instrument voids the warranty. If there is a failure during the warranty period, contact Heathrow Scientific[®] for warranty service.
17. Instrument must only be used with original manufacturer's NiMH batteries.
18. Batteries must be completely discharged before disposal. Disposal should be performed in accordance with all governmental regulations.


Warning!

Improper use of the instrument or the batteries such as short-circuiting, mechanical damage, or overheating, may cause an explosion.


Product Description

The Rota-Filler 3000[®] has been designed to assist in the filling and dispensing of pipettes. The adjustable pipet nozzle rotates to three positions for maximum comfort and convenience. It accepts graduated and volumetric pipettes made of glass or plastic. Use for measuring liquids from 0.1 to 100 mL. It is not designed for use with Pasteur pipettes.

If used correctly, the pipetted liquid will only contact the pipette. The vacuum or excess pressure necessary for filling or dispensing liquid is generated by a pump. The fill or dispense speed of the pump can be set using the pump speed control. Set on high speed for use with large-volume pipettes or low speed for use with small-volume pipettes.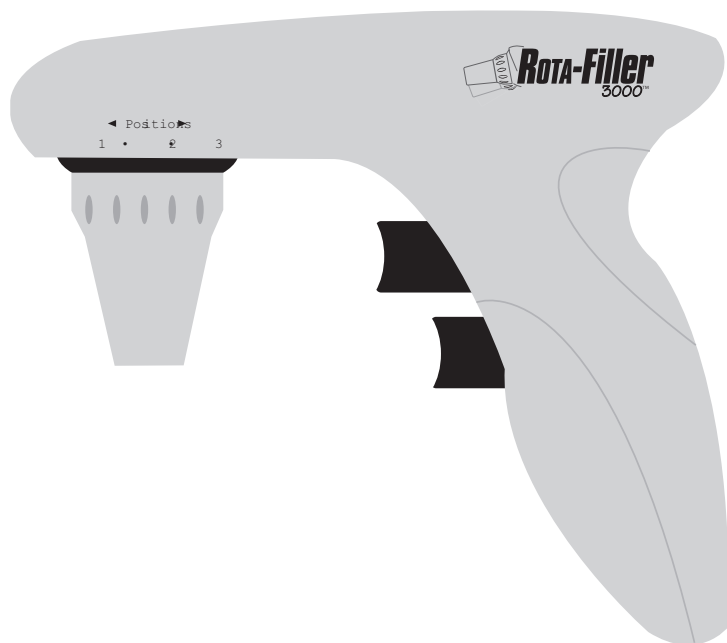

Components

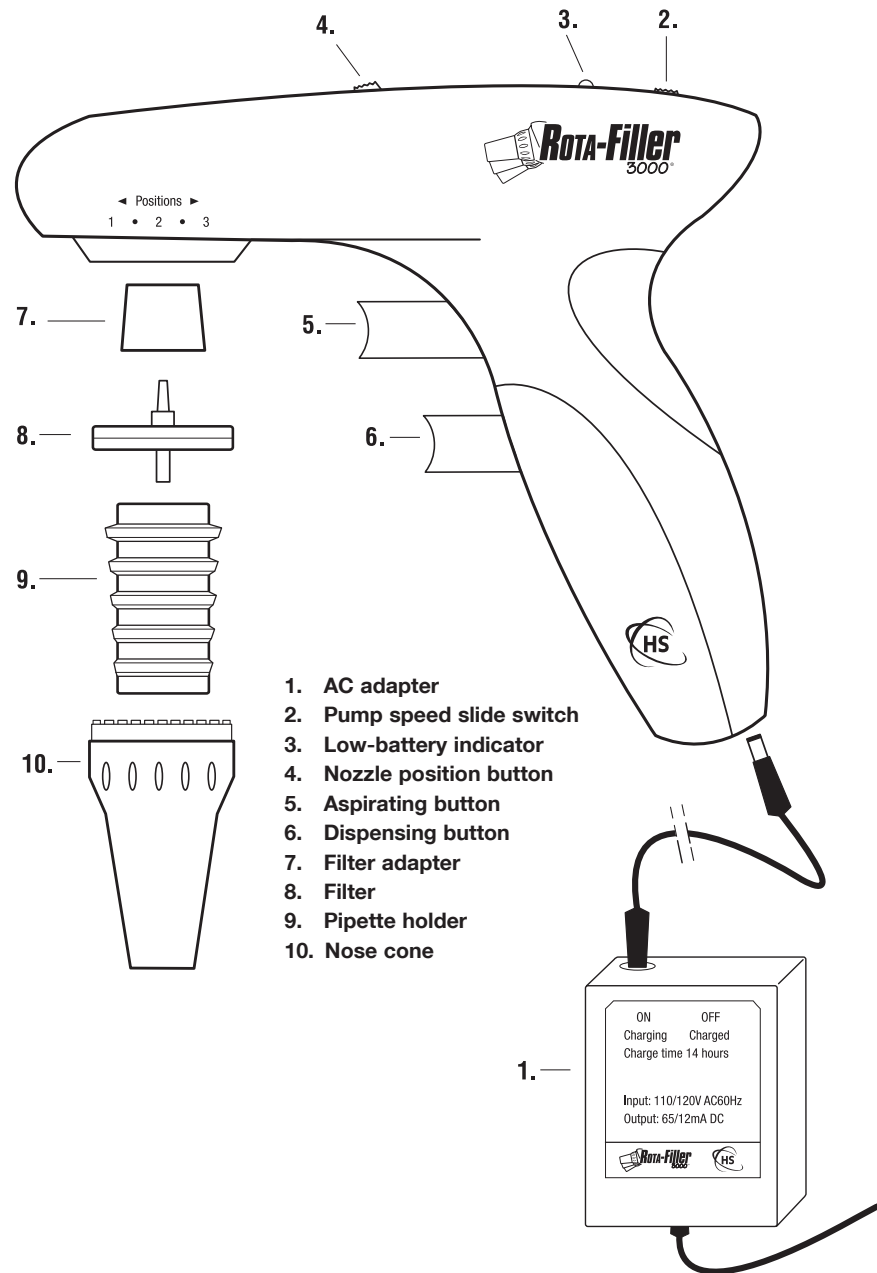Technical Data

Pipettes:	1 to 100 mL pipettes, glass or plastic
Working and recharging temperature:	50 to 95°F (10 to 35°C)
Batteries:	Two replaceable 1.2V nickel metal hydride (NiMH); approximately 8 hours operation
Charger:	65 mA/h for 14 hours
Materials	
Housing:	ABS
Nose cone:	Polypropylene
Pipette holder:	Silicone
Filter:	PTFE 25mm dia 0.45µm filter; Polypropylene housing

The Rota-Filler 3000[®] must not be used with liquids whose vapors are incompatible with silicone, ABS+Polyoxymethylene, PTFE, polypropylene, and NBR.

Warning!

This instrument must not be used or recharged in an explosive atmosphere. Highly flammable liquids (e.g. ether, acetone, and other liquids with a flash point below 0°C) must not be pipetted.

This mark is the confirmation that the unit conforms to the EU guidelines and has been tested according to the Low Voltage Directive 73/23/EEC and Electromagnetic Compatibility Directive 89/336/EEC.

Disposal Responsibilities:

The equipment you purchased may contain hazardous substances that could impact the environment. Per regulations on electronic devices in the European Community, you must use the appropriate disposal systems to avoid exposure of these substances to the environment. The disposal systems will reuse or recycle hazardous materials from your equipment responsibly.

The crossed-out wheeled bin symbol invites you to use those systems.

If you need more information on the collection, reuse and recycling systems, please contact your local or regional waste administration. You can also contact us for more information on the environmental performance of our products.

Getting Started

Remove the contents from the package, examining them carefully for breakage, defects, or shortages. Save the packaging.

Charging the Batteries

For transport safety reasons, the Rota-Filler 3000[®] is shipped with minimal battery charge. The batteries must be charged for 14 hours before initial use.

Warning!

Use original battery charger only!

The use of a different charger could damage the batteries and would void the warranty. Confirm that the voltage of the adapter matches the voltage of the electrical outlet.

Before starting to charge, confirm that the voltage or the battery adapter is identical to the main power supply at the outlet.

1. The temperature range for recharging is 50 to 95°F (10 to 35°C).
2. Plug the charger jack into the base of the Rota-Filler 3000[®].
3. Insert the plug of the adapter into a suitable wall outlet. The red indicator on the charger will light up.
4. Place the Rota-Filler 3000[®] on the benchtop stand.
5. Allow to charge 14 hours from empty to fully charged.
6. Once the Rota-Filler 3000[®] is fully charged, the adapter automatically switches to a trickle charge operation and the red indicator will turn off. If the charge is interrupted and then re-started, the adapter will start the full 14 hour charging cycle again. This will not damage the batteries.

Notes

- When fully charged, the Rota-Filler 3000[®] can be operated for approximately 8 hours.
- If using the Rota-Filler 3000[®] daily, do not charge the battery until the “low-battery indicator” LED has lit. A residual charge of approximately 10% is available when the LED has lit.
- The Rota-Filler 3000[®] can be operated during recharging.
- If the LED on the adapter does not light up when the Rota-Filler 3000[®] is connected to the adapter, the Rota-Filler 3000[®] and adapter should both be tested and if necessary, repaired or replaced.

Mounting the Holder

For storage, insert the Rota-Filler 3000® in the wall holder.

Mounting options:

1. The magnetized holder will grip a metal surface to provide support for the Rota-Filler 3000®.
2. Wall mount the holder with the included screws.

Using the Stand

For charging, place the Rota-Filler 3000® in either of the two orientations. The Rota-Filler 3000® can be charged in either position. When the Rota-Filler 3000® is placed with the handle lying down, a pipette can stay within the controller within the stand during charging. Do not place near doors or the edge of benchtop.

Pipetting


Adjusting the Nozzle

Press back the nozzle release button on the top of the Rota-Filler 3000® to adjust the pipet nozzle for maximum comfort and convenience. The wide-angle position is ideal for restricted spaces such as a fume hood. The smaller-angle position works well for tight benchtop applications.

Attaching the Pipette

Prior to attaching the pipette, inspect the pipette's top end for damage. A chipped pipette may damage the silicone pipette adapter.

Holding the pipette as near to the top as possible, carefully push it into the nose cone for a secure fit.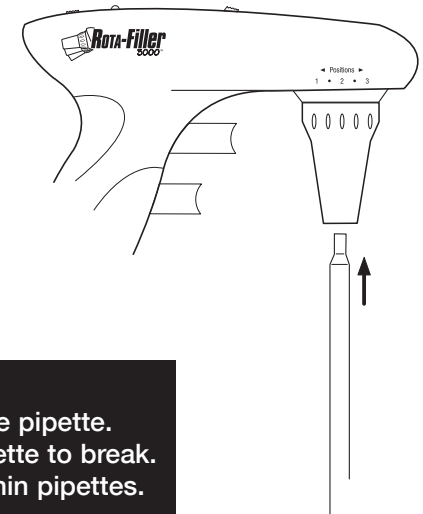

Caution!


Do not use force when inserting the pipette. Excessive force can cause the pipette to break. Use extra caution when inserting thin pipettes.

Once the pipette has been securely attached, always hold the Rota-Filler 3000® in a vertical position with the pipette tip facing down.

Filling the Pipette

Set the pump speed with the pump speed slide switch on the top of the Rota-Filler 3000®. Set on high speed for use with large-volume pipettes or low speed for use with small-volume pipettes. Select the gravity mode for “to contain” pipettes for low speed aspiration and gravity-driven delivery.

Immerse the pipette tip into the liquid. Slowly press the aspirating button (upper pipetting button). The speed of aspiration is dependent on the amount the button is depressed. Fill the pipette so that the meniscus is just above the desired volume.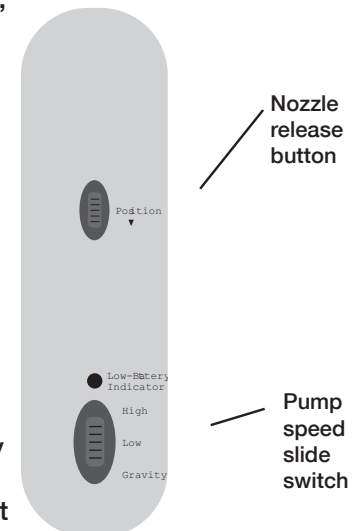

Wipe the outer surface of the pipette tip to remove residual liquid. Place the tip of the pipette against the side wall of the container. Slowly press the dispensing button (lower pipetting button) until the meniscus is exactly level with the desired volume.

Caution!


Take care to avoid aspirating the liquid into the Rota-Filler 3000[®] which can impair performance. If liquid is aspirated into the unit, clean the pipette adapter and nose cone per instructions under “Maintenance” (next page). If necessary, replace the filter.

Dispensing

Place the pipette tip at an angle against the inside of the receiving vessel. Slowly press the dispensing button (lower pipetting button). The dispensing speed is dependent upon the amount the button is depressed.

In the “Blow-Out” mode where the pump speed slide switch is set to “High” or “Low”, the dispensing button activates the pump to blow-out all liquid from the pipette.

In the “Gravity” mode, the dispensing button does not activate the pump. The liquid is dispensed by gravity. The pipette volume can be regulated using various degrees of finger pressure. This mode is for “To Contain” pipettes.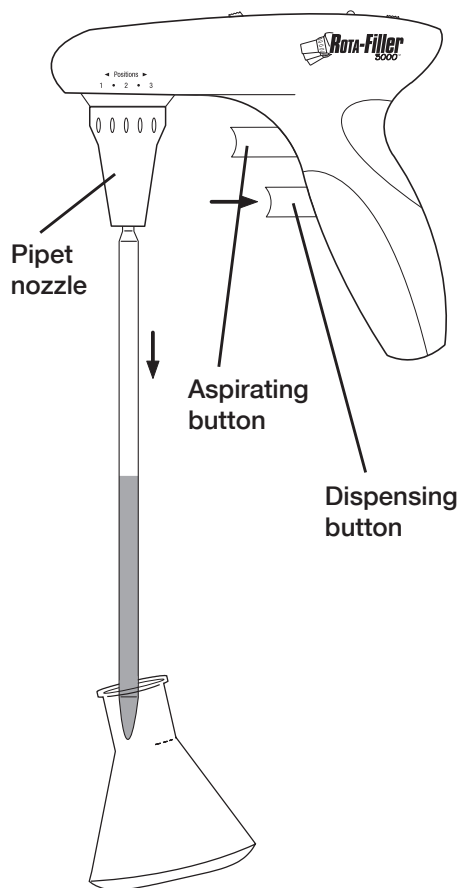

Maintenance

The Rota-Filler 3000[®] is normally maintenance-free. When necessary, the housing can be wiped occasionally using a damp cloth.

Cleaning the Pipette Adapter and Replacing the Filter

Caution!

When cleaning and exchanging the filter, wear protective gloves and safety goggles to avoid the risk of splashes and contamination.

If liquid has entered the nose cone or if the aspirating efficiency has decreased, observe the following instructions:

1. Hold the pipette over a container and remove the pipette.
2. Gently unscrew the nose cone.
3. Remove the filter.
4. Remove the pipette adapter from the nose cone by pushing it upwards.
5. Rinse the pipette adapter with demineralized water and allow to dry at room temperature.
6. With the Whatman[®] name and filter size facing down, push a new filter into the pipette adapter. The filter's smooth sided connector must be facing the pipette adapter. The filter's stepped, larger diameter connector should be facing the filter adapter. Push the filter adapter onto the filter.
7. Reassemble the Rota-Filler 3000[®] by reversing the steps above.
8. Test Rota-Filler 3000[®] for leaks, see “Checking for Leaks” (below).

Notes

- Not all membrane filters are suitable for use with the Rota-Filler 3000[®]. For proper performance, only use the manufacturer's recommended replacement filters.
- The pipette adapter, filter, and filter adapter may be steam sterilized at 121°C for 20 minutes.
- It is the user's responsibility to ensure proper sterilization.

Checking for Leaks

When the pipette has been filled, liquid should not drip from the pipette unless appropriate buttons are depressed. Should the pipette drip, disassemble the pipette adapter, filter, and filter adapter. Carefully reassemble. If necessary, replace worn parts. See “Ordering Information” (page 14).

Recharging the Batteries

One full charge of the NiMH batteries allows approximately 8 hours of continuous pipetting. Decreasing motor power and a lit “Low-Battery Indicator” LED, indicate the need for recharging.

For environmental reasons, the Rota-Filler 3000[®] is powered by rechargeable NiMH batteries, which do not contain Cadmium. These type of batteries work best when “trained”—the batteries are discharged fully before recharged for several recharge sessions.

Warning!

Use original battery charger only!

The use of a different charger could damage the batteries and would void the warranty.


Confirm that the voltage of the adapter matches the voltage of the electrical outlet.

Before starting to charge, confirm that the voltage of the battery adapter is identical to the main power supply at the outlet.

1. The temperature range for recharging is 50 to 95°F (10 to 35°C).
2. Plug the charger jack into the base of the Rota-Filler 3000[®].
3. Insert the plug of the adapter into a suitable wall outlet. The red indicator on the charger will light up.
4. Place the Rota-Filler 3000[®] on the benchtop stand.
5. Allow to charge 14 hours from empty to fully charged.
6. Once the Rota-Filler 3000[®] is fully charged, the adapter automatically switches to a trickle charge operation and the red indicator will turn off. If the charge is interrupted and then re-started, the adapter will start the full 14 hour charging cycle again. This will not damage the batteries.

Notes

- When fully charged, the Rota-Filler 3000[®] can be operated for approximately 8 hours.
- If using the Rota-Filler 3000[®] daily, do not charge the battery until the “low-battery indicator” LED has lit. A residual charge of approximately 10% is available when the LED has lit.
- The Rota-Filler 3000[®] can be operated during recharging.
- If the LED on the adapter does not light up when the Rota-Filler 3000[®] is connected to the adapter, the Rota-Filler 3000[®] and adapter should both be tested and if necessary, repaired or replaced.

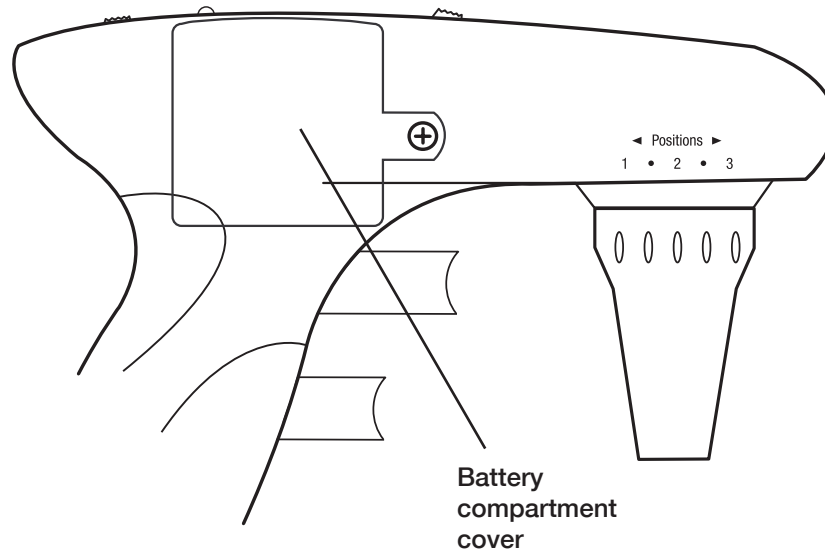Replacing the Batteries

Instrument must only be used with the original manufacturer’s NiMH batteries. Batteries must be completely discharged before disposal. Disposal should be performed in accordance with all governmental regulations.

Warning!

Improper use of the instrument or the batteries such as short-circuiting, mechanical damage, or overheating, may cause an explosion.

1. Place the Rota-Filler 3000[®] on flat surface with logo side down.
2. Use a Phillips-head screwdriver to remove the blue screw (see photo above).
3. Remove battery compartment cover.
4. Gently remove batteries.
5. Insert new batteries according to markings on plastic in the battery compartment.
6. Replace battery compartment cover.
7. Replace screw on battery compartment cover.

Warning!

Use original battery charger only!

The use of a different charger could damage the batteries and would void the warranty.

Confirm that the voltage of the adapter matches the voltage of the electrical outlet.

Trouble Shooting

Trouble	Possible Cause	Solution
Reduced aspirating efficiency	<ul style="list-style-type: none"> – Filter wetted – Batteries discharged 	<ul style="list-style-type: none"> – Replace filter – Recharge the batteries
Pipette drips	<ul style="list-style-type: none"> – Filter installed upside down – Pipette or pipette adapter damaged 	<ul style="list-style-type: none"> – Reinstall filter (see page 11) – Check pipette and pipette adapter; replace if damaged
Pipette is loose	<ul style="list-style-type: none"> – Pipette adapter is dirty or damaged 	<ul style="list-style-type: none"> – Clean pipette adapter; replace if damaged
Reduced operating time with fully-charged batteries	<ul style="list-style-type: none"> – Batteries worn – Instrument or adapter failed 	<ul style="list-style-type: none"> – Send in for repair (see page 15) – Send in for repair (see page 15)

If the above recommendations do not solve current problems, contact Heathrow Scientific® directly.

Ordering Information

Rota-Filler 3000® with 115 VAC adapter, holder, stand, and two spare filters.....	HS3000RFA
Pipette holder, silicone.....	HS3000RFP
Nose cone, polypropylene.....	HS3000RFR
Replacement filters, 0.45µm, pack of 5	HS3000RFC

Return for Repair

Important:

Transporting hazardous materials without a permit is a violation of federal law.

Heathrow Scientific® will not accept any Rota-Filler 3000® that is not appropriately cleaned and decontaminated.

In the unlikely event of repair, or where damage to the unit necessitates return, contact Heathrow Scientific® and obtain return authorization before sending your Rota-Filler 3000® for service.

Warranty

Limited Warranty

Heathrow Scientific® warrants the Rota-Filler 3000® will be free from material defects in workmanship and material for 1 year from the date of purchase.

If you believe that there is a defect in the product, you must, during the warranty period, notify Heathrow Scientific®, provide proof of purchase, and return the product to Heathrow Scientific® with a Return Authorization form. To obtain a Return Authorization form, please call 1-800-621-7193.

If Heathrow Scientific® is properly notified and, after inspection confirms that there is a defect and the warranty period has not expired, Heathrow Scientific® will repair, modify, or replace the product, at its sole option, at no charge.

OTHER THAN THIS LIMITED WARRANTY, HEATHROW SCIENTIFIC® MAKES NO WARRANTIES, EXPRESS OR IMPLIED, WITH RESPECT TO THE QUALITY OR PERFORMANCE OF THE PRODUCT, INCLUDING BUT NOT LIMITED TO THE IMPLIED WARRANTIES OF MERCHANTABILITY AND FITNESS FOR PARTICULAR PURPOSE WHICH ARE HEREBY DISCLAIMED AND EXCLUDED. HEATHROW SCIENTIFIC® WILL IN NO EVENT BE LIABLE FOR ANY LOSS OF USE, LOSS OF PROFITS, CONSEQUENTIAL, SPECIAL, EXEMPLARY OR PUNITIVE DAMAGES.

THIS WARRANTY DOES NOT COVER:

- ANY DEFECT OR DAMAGE CAUSED BY IMPROPER OR UNREASONABLE USE OF THE PRODUCT. (THE ROTA-FILLER 3000® IS DESIGNED FOR USE ONLY BY TRAINED LABORATORY TECHNICIANS. USE BY ANYONE ELSE WILL VOID THIS WARRANTY.)
- ANY PRODUCT THAT HAS BEEN, IN HEATHROW SCIENTIFIC®'S SOLE JUDGEMENT, TAMPERED WITH, ALTERED, OR REPAIRED BY ANYONE OTHER THAN HEATHROW SCIENTIFIC®.
- ANY PRODUCT THAT IS INOPERATIVE BECAUSE OF: (a) WEAR OCCASIONED BY USE, (b) NEGLIGENCE, (c) ACCIDENT, (d) INCORRECT MAINTENANCE, OR (e) USE UNDER ABNORMAL CONDITIONS OF TEMPERATURE, DIRT OR CORROSION, OR USE WITH ABRASIVE OR CORROSIVE MATERIALS.
- ACCESSORY PARTS, SUCH AS RUBBER PARTS, THAT ARE DAMAGED BY LIQUIDS OR MISUSE.

IN NO EVENT WILL HEATHROW SCIENTIFIC®'S OBLIGATION UNDER THIS WARRANTY EXCEED THE PRICE OF THE PRODUCT.

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All rights reserved.