REAX 20

Betriebsanleitung
Instruction Manual
Mode d'Emploi
Instrucciones de Empleo
Istruzioni per l'uso

D	DEUTSCH	Seite	3 - 12
E	ENGLISH	page	13 - 22
F	FRANCAISE	page	23 - 32
ES	ESPAÑOL	página	33 - 42
I	ITALIANO	pagina	43 - 52

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EN 61010-1 + EN 61010-2-051

Thank you for purchasing a Heidolph Instruments product. This item has been designed, made and inspected in compliance with DIN EN ISO 61010 for long-term and flawless operation.

The REAX 20 is an overhead mixer with a controllable speed range, designed to mix, settle out and separate soil samples. Stoppered cylinders and sediment deposit bottles between 160 and 270 mm high and a diameter up to 136 mm can be employed with the REAX 20. Due to its remarkable capability for continuous use, the REAX 20 is perfectly suitable for the analysis of water, sewage and sediment (according to DIN 38414 pt. 4: "determination of the elutriation of water").

SUMMARY

STANDARD HARDWARE & OPTIONS..... 14

GENERAL INFORMATION 14

SAFETY INFORMATION 15

SET-UP..... 16

1 **Set-up**..... 16

2 **Electric hook-up** 16

3 **Bottle installation** 16

3.1 **2000 ml bottles** 16

3.2 **1000 ml bottles** 17

3.3 **500 ml bottles** 17

OPERATION AND CONTROLS..... 18

EMERGENCY STOP 18

CLEANING & SERVICING..... 19

DISASSEMBLY & STORAGE..... 19

DISPOSAL 19

TROUBLESHOOTING 20

SPECIFICATIONS..... 20

WARRANTY, LIABILITY & COPYRIGHT 21

QUESTIONS / REPAIR WORK..... 21

CE-DECLARATION OF CONFORMITY..... 22

Important information

Advice about power cord / mains supply

Caution: mandatory action

Caution: fire- and explosion hazard

Advice about maintenance / repair

STANDARD HARDWARE & OPTIONS

	product	quantity	P/N 230/240V / 50/60Hz	P/N 115V / 50/60Hz
	REAX 20/4	1	541-20004-00	541-20004-05
	or			
	REAX 20/8	1	541-20008-00	541-20008-05
	or			
	REAX 20/12	1	541-20012-00	541-20012-05
	Instruction Manual	1	01-005-002-18	01-005-002-18

Options

	item	P/N
	Retrofitting set Reax 20 for 4 1-L bottles	549-26000-00
	Retrofitting set Reax 20 for 4 0,5-L bottles	549-27000-00
	Tensioning plate for caps Ø94 mm	11-001-001-81

GENERAL INFORMATION

Unpack your item carefully.

Inspect for damage and report such damage or missing parts to your supplier right away.

Read your instruction manual carefully. Take time to save time when working with your product. Make sure that every user has read and understood the Instruction Manual.

Please store the instruction manual in a place easily accessible to every user.

IF ALL ELSE FAILS, READ THESE INSTRUCTIONS !

A so-called EURO-plug (DIN 49441 CEE 7/VII 10/ 16 A 250 V) is standard on all of the products.

For the Continental US they feature a US-standard plug (NEMA Pub.No.WDI.1961 ASA C 73.1 . 1961 page 8 15A 125V).

For using the item in a country with deviating outlet / plug systems, we recommend to use approved adapters or to have an electrician replace the standard plug with one suiting your needs.

As shipped, the item features a protective ground wire. When replacing the original plug, make sure to reconnect this protective ground wire in the new plug !

SAFETY INFORMATION

Please comply with all safety and accident-prevention regulations as in force for laboratory work!

Use extra care when working with flammable substances; refer to safety data sheets.

Use extra care when working in the vicinity of flammable and explosive substances. Motors are of non-sparking type, the item itself however is not explosion-proof.

When connecting your item with your local power supply, please make sure your item is designed for your local voltage; go by the data plate on the item.

Please connect your item with a protective-ground outlet only.

Turn your power switch OFF whenever the item is not used, or before disconnecting the plug.

Repair work is limited to technicians approved by Heidolph Instruments.

Your item needs a solid stand.

Lab bench needs to be of rigid design, and features an anti-skid surface coat.

Before starting the item, make sure all samples are attached safely (must not move while shaking).

The unit must be placed on a stable, flat, level surface. Allow sufficient space around the unit for safe operation.

Take care that the turning rotor cannot collide with anything else on the laboratory table. Operating personnel have to be aware of the risk of accidents and observe safety rules.

**Before reactivating the Emergency Stop button turn main switch to OFF.
Else rotation will start immediately**

SET-UP

1 Set-up

The unit must be placed on a stable, flat, level surface. Allow sufficient space around the unit for safe operation.

Take care that the turning rotor cannot collide with anything else on the laboratory table. Operating personnel have to be aware of the risk of accidents and observe safety rules.

2 Electric hook-up

Connect mains plug.

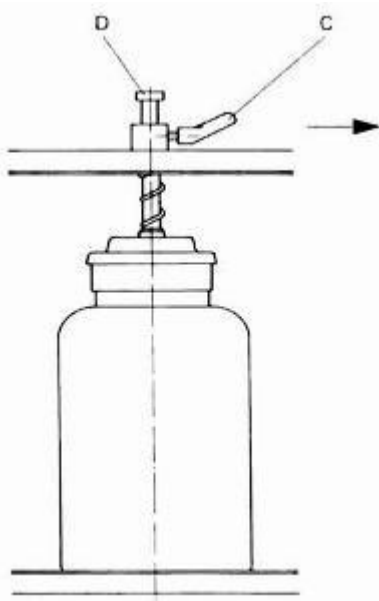
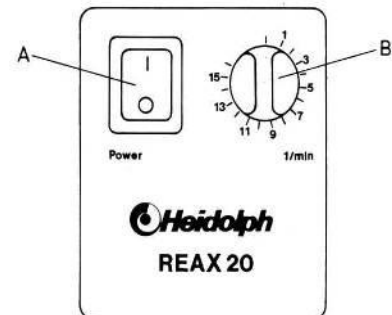
3 Bottle installation

3.1 2000 ml bottles

The standard version of the unit is designed for 2000 ml bottles. Set the rotor in the proper position in order to install the bottles.

To do this, switch (A) on and set speed control knob (B) to 1 rpm. Turn switch (A) off when levers (C) of one side are in the upright position.

Follow the instructions below and then repeat the entire process for the other side of the rotor.

Release lever (C) by turning it anticlockwise. Pull up on button (D) in order to raise the bottle holder upwards. Insert the bottle and take care that the bottle holder centers exactly above the screw cap of the bottle.

Push button (D) down in the direction of the bottle. Tighten lever (C) clockwise.

Repeat the procedure according to the number of bottles. Check the stability of the bottles and re-tighten them, as described above, if necessary.

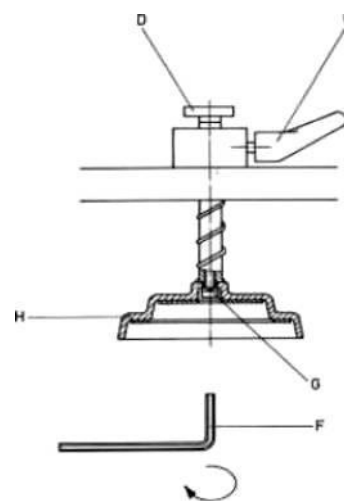
The angle of rotation of lever (C) is limited to 180°. In case this is not enough to tighten the bottle holder (H), it is possible to reposition the lever using a ratchet mechanism. Pull the lever in the direction shown in Diagram (inwards, towards the center) and hold, pull the lever in the anti-clockwise direction, release the inward pressure on the lever, and then tighten the lever in the clockwise direction. Always load the REAX 20 in a symmetrical way so as to distribute the weight equally.

3.2 1000 ml bottles

The bottle holders first have to be retrofitted with retrofitting set REAX 20 for 4 - 1 l bottles (available as an accessory, order no 549.26000.00).

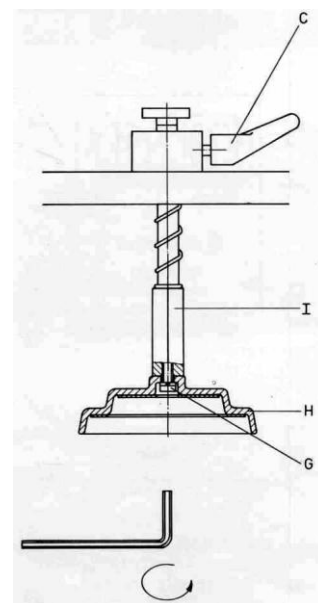
Proceed as follows:

- Tighten lever (C) clockwise.
- Loosen screw (G) with the supplied allen key.
- Remove bottle holder (H).
- Pull back the spiral spring while screwing in the extension pin (I) on the thread pin.
- Use allen key (F) to tighten the bottle holder (H) with screw (G) according to Diagram 2.
- 1000 ml bottles are inserted as described under chapter 3.1 above.

3.3 500 ml bottles

Follow the same procedure as in chapter 3.2 but use retrofitting set for 4 - 0.5 l bottles (order no.: 549.27000.00).

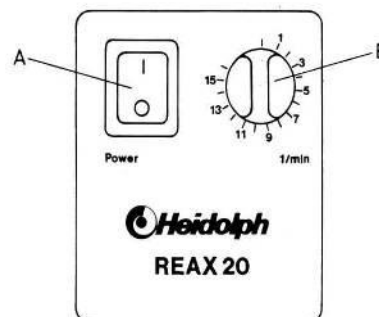
OPERATION AND CONTROLS

-  Before connecting power cord with mains outlet, make sure that:
 - your item is designed for your local voltage and frequency (data plate on item).
 - master switch is set to "0" and all controls are in "min." position (turn CCW completely);
-  Carefully close your vessels and select appropriate shaking intensity (if vessel remain open) to avoid splashes and spillage.
-  We recommend starting with low shaking speed and gradually increasing frequency to avoid accidental shaking at high speed.
-  When handling hazardous fluids, make reference to applicable safety information.


Make sure the bottles are inserted properly.

Switch on the instrument with switch (A) and the green pilot lamp in switch (A) lights. Set the required speed with speed control knob (B). The set speed is indicated on the scale.

EMERGENCY STOP

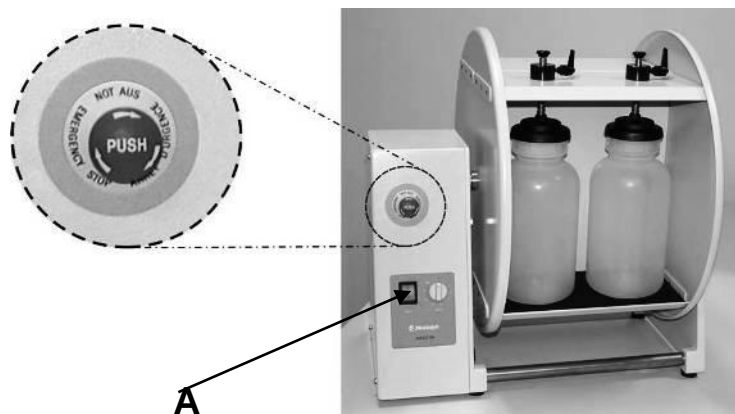-  REAX 20 is equipped with an emergency stop. In case of hazard from rotating parts press emergency stop!.
-  To reactivate, turn Emergency Stop button clockwise, and again press power switch (A). Make sure hazard is cleared.

CLEANING & SERVICING

Cleaning: wipe housing clean with a damp cloth (add some sort of mild liquid soap).

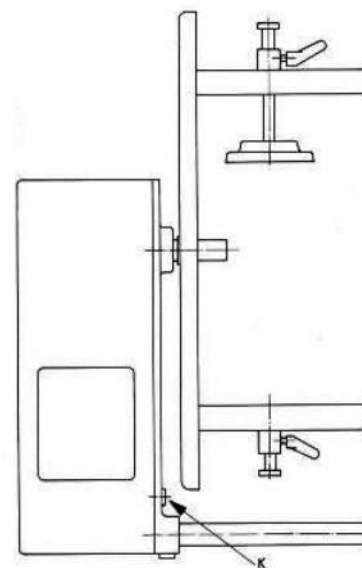
Note

To avoid damage to the surface finish, avoid using chlorine bleach, chlorine-based detergents, abrasive substances, ammonia, rags or cleaning agents containing metal particles.

The units have no parts that are subject to wear and tear and are maintenance free. A micro fuse accessible from the outside (see K) assures electric safety.

The item is maintenance-free. Repair work is limited to technicians so approved or appointed by Heidolph Instruments. Please call your local Heidolph Instruments Dealer for a Heidolph Instruments Field Representative (also refer to page 21)

DISASSEMBLY & STORAGE

DISASSEMBLY

Turn the unit off and remove the bottles.

Remove the plug from the electrical outlet.

TRANSPORT AND STORAGE:

In order to guarantee a safe transport the REAX 20 is fixed with two plastic transport straps in the original packing.

Place the unit and its parts into the original packing or a container with the necessary protection to prevent breakage during transport. Seal the original packing or container with packing tape.

Store the packed unit in a dry place.

CAUTION: When transporting this unit ensure the unit is not dropped or subject to unnecessary jolting.

DISPOSAL

For disposal, please comply with your local or national regulations. Divide into metal, plastic, etc.

Packing material to be treated as described above (material split).

TROUBLESHOOTING

Work on electric and electronic components is limited to qualified personnel.

No shaking

- check power cord
- check circuit breakers

- Thermal motor circuit breaker triggered by motor overload
Remedy:
Wait about 20 minutes, decrease load applied on shaker plate.
- Mechanical parts broken (humming motor noise) or electronic failure (no motor noise).
- Micro-fuse (K) blown, change fuse.

SPECIFICATIONS

REAX 20

Power supply	230/240 V 50/60 Hz / 115 V 50/60 Hz (optional)
Input power	280 W
Shaker frequency	1 – 16 1/min
Speed control	electronically, continuously variable,
Speed display	analog scale
Mixing movement	overhead rotation
Drive	capacitor motor, motor overheating protection with automatic reset
Capacity	Reax 20/4 - 4 bottles Reax 20/8 - 8 bottles Reax 20/12 - 12 bottles
Max. bottle diameter	136 mm
Max. bottle height	270mm
Standard bottle	a plastic wide necked bottle with a volume of 2000 ml. Along with the proper accessories, the 500 ml and 1000ml (sizes are also applicable)
Dimensions (l x b x h)	REAX 20/4: 490 x 465 x 520 mm REAX 20/8: 770 X 465 x 520mm REAX 20/12: 1050 x 465 x 520 mm
Weight	Reax 20/4 23 kg Reax 20/8 28 kg Reax 20/12 33 kg
Protection class	IP 21
ambient temperature	0°C to 40°C at 80% rel. humidity Approved for installation in gassing and conditioning cabinets (make reference to temperature limits)

WARRANTY, LIABILITY & COPYRIGHT

Warranty

Heidolph Instruments warrants that the present product shall be free from defects in material (except wear parts) and workmanship for three years from the date shipped off the manufacturer's warehouse.

Transit damage is excluded from this warranty.

To file for such warranty service, contact Heidolph Instruments (phone:++49-9122-9920-68) or your local Heidolph Instruments Dealer. If defects in material or workmanship are found, your item will be repaired or replaced at no charge.

Misuse, abuse, neglect or improper installation is not covered by this warranty.

Alterations to the present warranty need Heidolph Instruments' consent in writing.

Exclusion Clause

Heidolph Instruments cannot be held liable for damage from improper use or misuse. Remedy for consequential damage is excluded.

Copyright

Copyright in pictures and wording of the present Instruction Manual is held by Heidolph Instruments.

QUESTIONS / REPAIR WORK

If any **aspect** of installation, operation or maintenance remains unanswered in the present Manual, please contact the following address:

For repair services please call Heidolph Instruments (phone: +49 - 9122 - 9920-68) or your local, authorized Heidolph Instruments Dealer.

Note

You will receive approval for sending your defective item to the following address:

Note

If you are based in the United States of America, please contact Heidolph US:

Heidolph Instruments GmbH & Co. KG
Lab Equipment Sales
Walpersdorfer Str. 12
D-91126 Schwabach / Germany
Tel.: +49 – 9122 - 9920-68
Fax: +49 – 9122 - 9920-65
E-Mail: sales@heidolph.de

Heidolph Instruments, LLC
Lab Equipment Sales
2615 River Rd.
Cinnaminson, NJ 08077
Phone: 856-829-6160
Fax: 856-829-7639
E-Mail: heidolph@snip.net

Safety Information

When shipping items for repair that may have been contaminated by hazardous substances, please:

- advise exact *substance*
- take proper protective measures to ensure the safety of our receiving and service personnel
- mark the pack *IAW Hazardous Materials Act*

CE-DECLARATION OF CONFORMITY

We herewith declare that the present product complies with the following standards and harmonized documents:

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