REAX top

REAX control

Betriebsanleitung
Instruction Manual
Mode d'Emploi
Instrucciones de
Empleo
Istruzioni per l'uso

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EN 61010-1 + EN 61010-2-051

CONTENTS

Illustrations	5-6
Packing list	2
Brief description/General information	15
Safety instructions	16
A. Setting up	16
1. Setting up the appliance	16
2. Electrical connection.....	16
3. Assembly and accessories	16
3.1. Tray, small and tray, large	16
B. Control and operation	17
4. ON/OFF switch and setting speed.....	17
5. Shaking with tilt operation.....	17
6. Shaking in continuous operation.....	17
Routine cleaning and maintenance	17
Dismantling, Transport and storage.....	18
Disposal.....	18
Faults and how to remedy them	18
Technical Data.....	19
Warranty, Liability & Copyright	20
CE-Declaration of conformit.....	20

BRIEF DESCRIPTION

- REAX top and REAX control are the successor models to the well-known, tried and tested test tube shaker REAX 2000.
 - REAX top is equipped with speed control.
 - REAX control is equipped with speed control for a constant speed even in the low range.
- On both models, shaking is actuated by the pressure of the test tube on the tray. What is new is the practical switchover function from periodic to continuous operation.

GENERAL INFORMATION

-  Unpack the appliance and parts with care. Compare the contents with the packing list. Check for damage and report any defects or deviations from the packing list to the supplier without delay.
-  Read the operating instructions carefully and in full and ensure that everyone who operates the appliance has read the instructions carefully before use.
-  Keep the operating instructions in a place which is accessible to everyone.
-  This appliance comes fitted with a Euro plug (DIN 49441 CEE 7/VII 10/16 A 250 V) or, for North America, with a US standard plug (NEMA Pub.No.WDI-1961 ASA C 73.1-1961 page 8 15 A 125 V).
- If the appliance is to be connected up to different grid, an approved adapter must be used or the plug supplied replaced by a qualified electrician with a 3-pin plug approved for the grid in question
- The appliance is earthed upon delivery. If the original plug is replaced, make sure the appliance's protective conductor is connected to the new plug.

SAFETY INSTRUCTIONS

Observe all safety and accident prevention regulations in the laboratory!

The motor works without sparks. Nevertheless, the manufacturer cannot guarantee that the appliance is explosion-safe.

Wear safety goggles and suitable working clothes when working with the shaker.

Repairs must be carried out only by technicians authorized.

Extreme care is called for when handling toxic or other hazardous substances.

To prevent the shaker sliding off the table top as a result of the vibrations generated during operation, make sure the feet are securely attached to the table top.

A. SETTING UP

1. Setting up the appliance

- Set up the test tube shaker on a smooth and solid table top. The weight of the appliance and the 4 parabolically shaped rubber feet make for stability.
- The feet can adhere securely to the table top only if both the feet and the surface to which they are to be attached are clean. If necessary, clean the feet with a damp cloth and ethanol.

2. Electrical connection (Figure 1, 2)

- Before connecting the appliance into the mains, make sure that the mains voltage matches that required on the type plate.

- (for other plug standards, see the General Information on page 15).

- the ON/OFF switch (1) is switched off (in the “0” position).

- the speed setting knob (2) is turned to the left as far as it will go.

- Plug the shaker into the mains using the mains lead. The plug for the mains lead is at the rear of the appliance.

3. Assembly and accessories

3.1. Tray, small (accessory 23.07.06.05.01) and tray, large (accessory 549.19000.00) (Figures 1, 2, 4)

- The tray (9) is simply clipped onto and pulled off the appliance by hand. When clipping on the tray, make sure the slot in the coupling area of the tray is in line with the lug on the eccentric of the appliance.

B. CONTROL AND OPERATION

4. ON/OFF switch and setting speed (Figures 1, 2)

- The appliance is switched on at the ON/OFF switch (1). The green control lamp in the ON/OFF switch (1) indicates that the appliance is ready for operation.
- The speed (shaking intensity) is selected using the speed setting knob (2). It can be increased by turning this knob to the right.
- The REAX top has a dial (5) and the REAX control a scale (5) with precision speed display in 1/min.

5. Shaking with tilt operation (Figures 1, 2, 8)

- Before starting the shaker, select a low speed to prevent splashing when the appliance starts up.
- The operating mode is selected with the slide switch (6). To select tilt operation, turn the slide switch to the right as far as it will go. The symbol ▼ (7) to the left of the slide switch (6) indicates the operating mode.
- Shaking is actuated by the pressure of the test tube on the tray and stops again automatically as soon as the test tube is removed.

6. Shaking in continuous operation (Figures 1, 2, 7)

- Before starting the shaker, select a low speed to prevent splashing when the appliance starts up.

- The operating mode is selected with the slide switch (6). To select continuous operation, turn the slide switch to the left as far as it will go. The symbol ● (8) to the right of the slide switch (6) indicates the operating mode.
- In this position, the tray are shaken continuously.

ROUTINE CLEANING AND MAINTENANCE

- Use a damp cloth (mild soap solution) to wipe clean the housing and surfaces of the appliance.
- Important: Never use chlorine bleach or chlorine-based detergents, scouring agents, ammonia, cleaning wool or cleaning agents with metallic components to clean the appliance, as these would damage the surfaces of the appliance.
- The appliance is maintenance-free. Any repairs which may be necessary must be carried out by an authorized Heidolph technician. Please contact your local Heidolph dealer or Heidolph representative (see page 20).

- To ensure the feet adhere securely to the table top, clean the feet at regular intervals using a damp cloth and ethanol.

DISMANTLING, TRANSPORT AND STORAGE

Dismantling

1. Switch off the appliance and disconnect it from the mains (allow to cool).
2. Remove any apparatus around the shaker so that the appliance can be dismantled without difficulty. Remove the glass set and dismantle the entire appliance and all the parts in the reverse order to assembly.

Transport and storage

1. Place the appliance and the parts in the original packing or in another suitable container in order to avoid damage during transit. Seal the packing with adhesive tape.
2. Keep the appliance in a dry place.

Caution: Avoid knocks and vibrations when transporting the appliance.

DISPOSAL

Please make sure that any old or defective appliances are taken to a collection point and disposed of professionally. Collect the metal, glass and plastic parts separately.

Please also dispose of the packing materials in an environment-friendly manner (sorting).

FAULTS AND HOW TO REMEDY THEM

No shaking movement:

- Check the mains lead and power supply.
- The ON/OFF switch is not switched on (see Operation, page 17).
- The slide switch is not in the right position (see Operation, page 17).
- The speed selected is too low (increase speed).
- In tilt operation, not enough pressure on the tray (increase pressure by hand).
- In tilt operation, the reaction vessel exerts pressure on the tray at an angle (switch over to continuous operation, if necessary) (see Operation, page 17).
- The motor is defective (requires service).
- The switch is defective (requires service).
- The control electronics are defective (requires service).

TECHNICAL DATA

Type	REAX top	REAX control
Order number	541.10000.00	541.11000.00
Type of motion	circular and vibrating	circular and vibrating
Shaking frequency [1/min]	0-2400	0-2400
Orbit [mm]	5	5
Speed display	dial	scale
Control precision	uncontrolled	± 2%
Switch	for periodic and continuous	for periodic and continuous
Connected load [W]	51	51
Unit weight [kg]	2,8	2,8
Dimensions (W x H x D) [mm]	134 x 105 x 172	134 x 105 x 172
Voltage/frequency	230 V/50 Hz (others also available)	
Permissible duty factor	100%	
Overheating protection prolonged	The motor switches off after overloading with self- resetting function	
Degree of protection	IP 22	
Permissible ambient temperature	5 - 40°C at 85 % relative humidity, no condensation	

WARRANTY, LIABILITY & COPYRIGHT

Warranty

Heidolph Instruments warrants that the present product shall be free from defects in material (except wear parts) and workmanship for 3 years from the date shipped off the manufacturer's warehouse.

Transit damage is excluded from this warranty.

To obtain such warranty service, contact Heidolph Instruments (phone: +49 – 9122 - 9920- 68) or your local Heidolph Instruments Dealer. If defects in material or workmanship are found, your item will be repaired or replaced at no charge.

Misuse, abuse, neglect or improper installation are not covered by this warranty promise. Alterations to the present warranty promise need Heidolph Instruments' consent in writing.

Exclusion Clause

Heidolph Instruments cannot be held liable for damage from improper use or misuse. Remedy for consequential damage is excluded.

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Note

You will receive clearance for sending your defective item to the following address:

Heidolph Instruments GmbH & Co. KG
Lab Equipment Sales
Walpersdorfer Str. 12
D-91126 Schwabach / Germany
Tel.: +49 – 9122 - 9920-74
Fax: +49 – 9122 - 9920-65
E-Mail: sales@heidolph.de

Safety Information

When shipping items for repair that may have been contaminated by hazardous substances, please:

- advise each *substance*
- take adequate *protective* action not to harm our parts receiving and service personnel
- *mark* the pack IAW Hazardous Materials Act

QUESTIONS/REPAIR WORK

If any aspect of installation, operation or maintenance remains unanswered in the present manual, please contact us at the following address:

For equipment repairs jobs please call Heidolph Instruments (phone: +49 – 9122 - 9920-68) or your local authorized Heidolph Instruments dealer.

CE-DECLARATION OF CONFORMITY

We here with declare that the present product complies with the following standards and harmonized documents:

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