



Specifications	
External Dimensions mm (WxDxH)	650 x 600 x 1982
Internal Dimensions mm (WxDxH)	600 x 600 x 760
Working Aperture mm (WxH)	590 x 300
Face Velocity	0.5 m/sec
Airflow	320 m ³ /hr
Carbon Filter Type	Formalin
Filter Size mm	600 x400 x 115
Filter Weight	14kg
Pre-Filter	Filtrete
Internal Lighting	18 watt fluorescent
Fan	Centrifugal, Brushless & Sparkless
Water Supply	Optional
Foot Switch Control	Optional
Hand Operated Controls	Standard
Built-in 50 litre Formalin Tank	Optional
Noise Level	<57 db(A)
Electrical Supply	2 x 240v - 400watts
Construction	Head: Epoxy Powder Coated Zintec Steel with 6mm clear acrylic front and side upper panels.

Circulaire 650FDS Compact Formalin Dispensing Station

The Circulaire 650FDS provides a compact and safe solution to the dispensing of Formalin in Operating Theatres, Pathology Laboratories & Mortuaries.

It is free standing and has a footprint of less than 0.4m². It is extremely simple to use, just place your container of formalin into the base section, connect up and press the front mounted button (or optional foot switch) to dispense the formalin through the tap.

The upper cabinet provides full fume containment making sure that all formalin fumes are kept well away from the operator. The large capacity carbon filter ensures that no fumes are returned into the working area. This filter provides 3 to 4 years of life under normal use and is extremely easy to change when required. Where additional safety is required an outlet carbon filter can be fitted along with a filter saturation alarm to detect chemical breakthrough.



WolfLabs

Pricing on any accessories shown can be found by keying the part number into the search box on our website.

The specifications listed in this brochure are subject to change by the manufacturer and therefore cannot be guaranteed to be correct. If there are aspects of the specification that must be guaranteed, please provide these to our sales team so that details can be confirmed.

www.wolflabs.co.uk

Tel : 01759 301142

Fax : 01759 301143

sales@wolflabs.co.uk

Please contact us if this literature doesn't answer all your questions.