
HRD

DIGITAL HAND REFRACTOMETER

 **OPERATION MANUAL**

 **GUIDA UTENTE**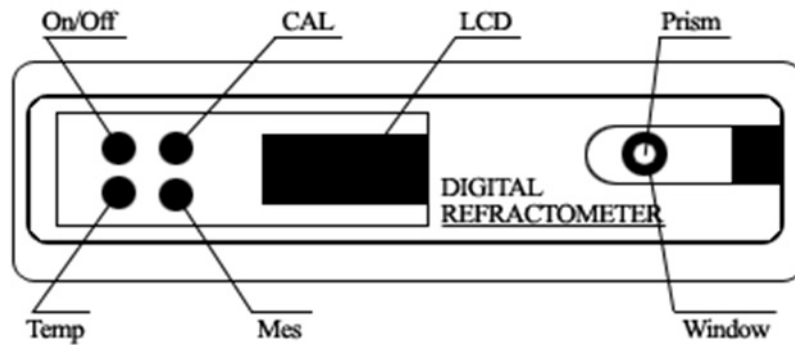

OPERATION STEPS

A. Original Point Setting (Please do not forget this step)

1. Completely clear prism window .
Drop standard solution of distilled water to fill the prism window.

2. Hold “**CAL**” button for five seconds till it twinkles “rEF” on LCD. Within 10 seconds after showing “rEF”, press **Mes** button and hold it for five seconds. The instrument start to set Original Point. After Original Point Setting is finished, “End” will be displayed on LCD.

B. Measurement

1. Completely clear prism window. Drop solution to fill the prism window.

2. Press **Mes** button

C. Conversion of Status of Scales

1. Turn it off.
2. Repeat step 1-2 to set other scale. (The chosen status of scale will be stored no matter if power is off.)
3. Synchronously hold “**CAL**” button and **Mes** button. Turn it on by pressing **On/Off** button.

OTHER OPERATION

1. Battery replacement

In the case of no battery or low battery, a low battery alert of “ on the upper left of LCD, you should install or replace the battery. Open the battery cover on the back of the instrument, install a battery or replace the old battery with a new one.

2. Power On and Off

Press the “**On/Off**” button to turn it on or off. Temperature in Centigrade will display when power on. The instrument will turn off automatically if there is no any operation in three minutes.

3. Measurement of Temperature

Temperature can be displayed at any time by press the button during measurement. The conversion of Fahrenheit and Centigrade can be made by pressing **Temp** button.

PRECAUTIONS


1. Don't expose the instrument to low, high temperature and sun light for long time, to avoid LCD losing effectiveness.
2. Because the instrument is very precise, avoid violent shocks.
3. To avoid damage, don't disassemble and assemble the instrument or change the inner circuit and parts.
4. Zero-setting should be implemented strictly according to instrument.
5. Be sure to clean the prism surface and window of stage before and after every measurement.
6. To avoid that accuracy is affected by evaporation, be sure to implement measurement immediately after dripping solution on window of stage.
7. It might cause wrong results if kept measuring under low voltage.
8. Don't use it under strong light (as sunlight, lamp etc.)
9. Don't use the instrument in humid or corrosive environment.
10. Prevent liquid entering into battery house.

Art.13 Dlsg 25 July 2005 N°151. "According to directives 2002/95/EC, 2002/96/EC and 2003/108/EC relating to the reduction in the use of hazardous substances in electrical and electronic equipment and waste disposal."

The basket symbol on equipment or on its box indicates that the product at the end of its useful life should be collected separately from other waste.

The separate collection of this equipment at the end of its lifetime is organized and managed by the producer. The user will have to contact the manufacturer and follow the rules that he adopted for end-of-life equipment collection. The collection of the equipment for recycling, treatment and environmentally compatible disposal, helps to prevent possible adverse effects on the environment and health and promotes reuse and/or recycling of materials of the equipment. Improper disposal of the product involves the application of administrative penalties as provided by the laws in force.

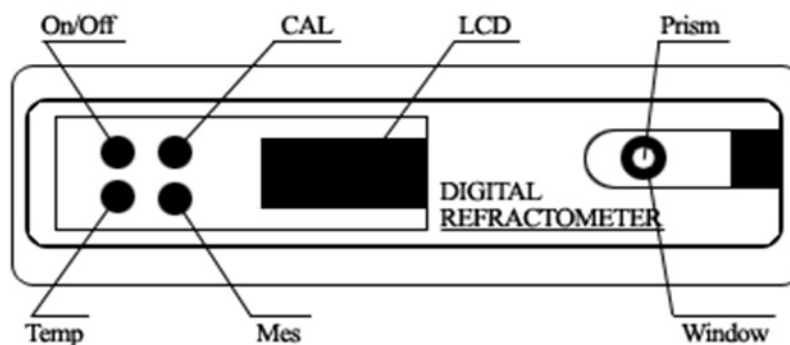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