Digital Hotplates & Stirrers


Instructions for use

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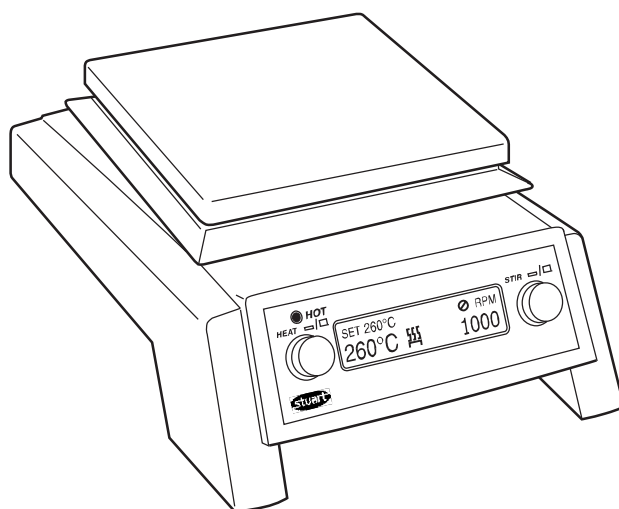

Barloworld
Scientific

English

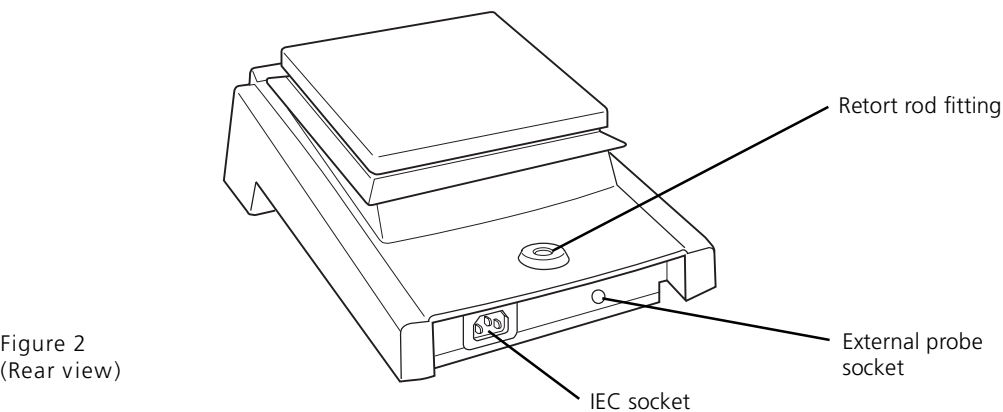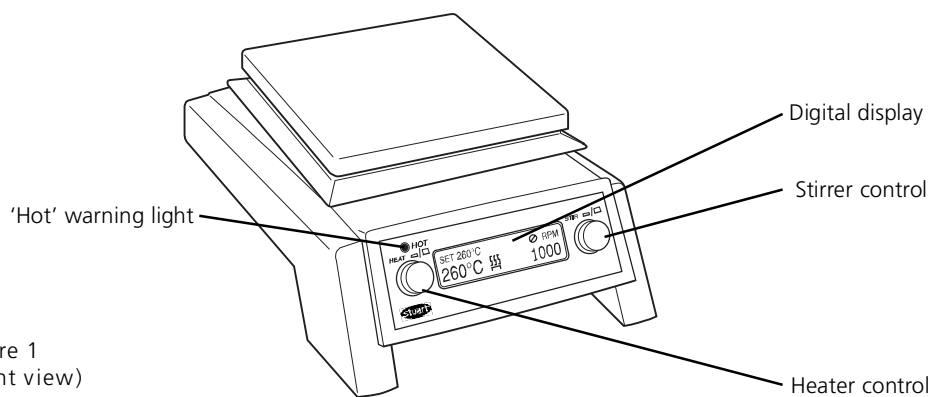Digital display

Display at switch on, heater & stirrer off, probe attached.

Display at switch on, heater & stirrer off, probe removed.

Full display with probe attached.

Full display with probe removed.

Thank you for purchasing this Barloworld Scientific product. To get the best performance from the equipment, and for your personal safety, please read these instructions carefully before use.

Barloworld Scientific hotplates and stirrers are designed to operate under the following conditions:-

- ❖ For indoor use only
- ❖ Use in a well ventilated area
- ❖ Ambient temperature range +5°C to +40°C
- ❖ Altitude to 2000m
- ❖ Relative humidity not exceeding 80%
- ❖ Mains supply fluctuations not exceeding 10% of nominal
- ❖ Overvoltage category II IEC60364-4-443
- ❖ Pollution degree 2 IEC664
- ❖ Use with a minimum distance all round of 200mm from walls or other items

If the equipment is not used in the manner described in this manual the protection provided by the equipment may be impaired.

Electrical Installation

THIS EQUIPMENT MUST BE EARTHED

Before connection please ensure that the line supply corresponds to that stated on the rating label.

Power requirements:

Model		
CD162	230V, 50/60Hz,~	550W
SD160	230V, 50/60Hz,~	700W
SD162	230V, 50/60Hz,~	750W
CD162/LV	100-120V, 50/60Hz,~	550W
SD160/LV	100-120V, 50/60Hz,~	700W
SD162/LV	100-120V, 50/60Hz,~	750W

There is an IEC socket at the rear of the instrument for connection to the mains supply, (see figure 2). The unit is supplied with two mains leads fitted with IEC plugs for connection to the instrument. One has a U.K. 3 pin plug and the other has a 2 pin "Shuko" plug for connection to the mains. Choose the lead appropriate for your electrical installation and discard the other.

Should neither lead be suitable, take the lead with the U.K. plug and replace the plug with a suitable alternative. See the enclosed instruction

sheet for advice on how to carry out this procedure.

Should the mains lead require replacement a cable of 1mm² of harmonised code H05W-F connected to an IEC 320 plug should be used.

IF IN DOUBT CONSULT A QUALIFIED ELECTRICIAN

The mains lead should be connected to the instrument BEFORE connection to the mains supply.

Operation

Use as a Hotplate

CD162, SD160 and SD162

Ensure that the external temperature probe is NOT connected to the instrument.

The control knob on the left hand side of the display controls the heating function.

To switch on the heating function briefly press and release the knob. Repeat to turn off the heating function.

When switched on the display will show actual plate temperature at the bottom of the screen and the set temperature directly above. The heating icon will show , indicating that the displayed temperatures are those of the plate surface.

The plate temperature can be selected and controlled to a resolution of 1°C.

To increase the set temperature turn the knob clockwise until the required setting is shown in the display. To decrease the setting turn the knob anti-clockwise.

The unit will now begin to heat until the set temperature is reached and then will hold the plate at this value. The actual plate temperature is continuously displayed at the bottom of the screen. When the set temperature is reached an alarm will sound. When the plate surface is too hot to touch (> 50°C) a red warning light on the front panel will begin to flash. This will continue to flash while the plate temperature is above 50°C even after the heating function is turned off.

WARNING: The top surface of the instrument may be HOT. In free air a

surface temperature of 450°C can be achieved on the ceramic top CD162 model and 325°C on the metal top SD160 and SD162 models. Do not leave heaters switched on when not in use.

Use as a Stirrer

The control knob on the right hand side of the front panel controls the stirrer speed.

To switch on the stirring function briefly press and release the knob. Repeat to turn off the stirring function. When switched on the icon on the display will revolve and the display will show the set speed.

Turning the knob clockwise increases speed from 100rpm to 1300rpm. Turning the knob anti-clockwise decreases the speed.

Use as a Temperature Controller using the External Probe SD162, CD162 only

These units are supplied with a temperature probe to allow direct control of liquid temperature. NOTE: Before attaching probe, mains supply must be turned off.

To use the probe first insert the connector into the socket on the rear of the instrument (see fig 2) and place the probe into the vessel containing the liquid. Switch on the heating function by briefly depressing the control knob. The heating icon in the display will show  indicating that the displayed temperatures are those of the liquid rather than the plate surface.

When switched on the display will show actual liquid temperature at the bottom of the screen and the set temperature directly above.

The liquid temperature can be selected and controlled to a resolution of 1°C.

To increase the set temperature turn the knob clockwise until the required setting is shown in the display. To decrease the setting turn the knob anti-clockwise.

The unit will now begin to heat until the set temperature is reached and then will hold the liquid at this value. The actual liquid temperature is continuously displayed at the bottom of the screen.

Note that the maximum liquid temperature is limited to 200°C.

Guidance Notes

1. For optimum temperature control please ensure the contents of the beaker are stirred adequately.
2. Keep the unit out of strong drafts.
3. Allow the temperature adequate time to stabilise.

Safety Advice

- ❖ Do not use the instrument to heat inflammable liquids.
- ❖ Never lift or carry the instrument until it has been switched off and allowed to cool for at least 30 minutes. The hot warning light will give guidance.
- ❖ The unit should be carried using both hands with the fingers under the side edges.
- ❖ Never move or carry the unit with containers on the top plate or while still connected to the mains supply.
- ❖ There is a danger of liquid spillage if containers are over-filled and stirred at high speed. Always build stirrer speed slowly and never stir more rapidly than necessary.
- ❖ NEVER place a cold glass vessel onto a hotplate which is already hot.

Cleaning and Care

Before attempting cleaning:-

Ensure that the top is cool

Disconnect from the mains electricity supply.

The metal casing should be cleaned using a damp cloth and a mild detergent solution.

Ceramic top units:

A damp cloth will normally remove most types of contamination. For more difficult stains a domestic cream cleanser is recommended.

Cleaning is made easier if spillages are attended to promptly. In any case, spillages of alkali, phosphoric acid and hydrofluoric acid MUST be removed immediately as these chemicals can attack and damage the glass ceramic. Ensure that the appropriate safety precautions are observed.

During cleaning and general operation take care not to scratch the surface of the top plate as this could result in subsequent thermal breakage.

WARNING: A ceramic top which is scratched, chipped, chemically etched or damaged must not be used.

Metal top units:

The metal top plate should be cleaned using a damp cloth and a mild detergent solution. Cleaning is made easier if spillages are attended to promptly. In any case, spillages of acids and alkalis MUST be removed immediately as these chemicals can attack and damage the the surface of the aluminium alloy. Ensure that the appropriate safety precautions are observed.

Accessory

A rod is available to support apparatus used with the instrument. The instrument is equipped with a fitting on the rear to accept the support rod - catalogue number SR1, (see figure 2). To fit the rod to your instrument, first isolate from the mains supply and allow to cool. Then simply screw the threaded end of the support rod into the fitting on the rear of the instrument.

Fault finding

In the event of malfunction the units can self diagnose certain fault conditions. In this case the display will show a message in the form ERROR where is a 4 digit code which identifies the fault.

If the error message appears followed by any other code it indicates an internal component failure. Turn the unit off, isolate from the mains electricity supply and wait two minutes before reconnecting and switching back on. If error codes other than those listed above continually display contact the Technical Service Department of Barloworld Scientific Ltd for advice.

Certain faults can be easily remedied by the user and are identified as follows:-

Error code	Fault condition	Remedy
000C	Probe has become disconnected while in use	Reconnect probe
0001	Temperature probe out of sample	Replace probe in sample container
0002	Probe failed open circuit	Renew probe
0004	Probe failed short circuit	Renew probe
0008	Probe temperature above 200°C	Turn unit off and allow to cool
0010	Over temperature protection activated	Turn unit off and allow to cool
0200	Internal case temperature too high	Turn unit off and allow to cool. Ensure the unit is well ventilated before attempting to operate again.

Servicing and Repair

This product range does not require any routine servicing.

Note: There are no internal user replaceable parts.

In the event of product failure it is recommended that any repair is only undertaken by suitably qualified personnel. For advice or to receive a service manual please contact the Technical Service Department of Barloworld Scientific Ltd.

Only spare parts supplied by Barloworld Scientific or its agent should be used. Fitting of non-approved parts may affect the performance of the safety features of the instrument.

Note: The magnetic stirrer drive utilises strong magnets.

If in doubt, please contact the Technical Service Department of
Barloworld Scientific Ltd.
Stone, Staffordshire ST15 0SA
United Kingdom
Tel: +44 (0) 1785 812121
Fax: +44 (0) 1785 813748
e-mail equipment@barloworld-scientific.com
www.barloworld-scientific.com

Spares

The following spare parts are available from your laboratory supplier. For a comprehensive parts list please contact the Technical Service Department of Barloworld Scientific Ltd quoting model and serial number.

Description	Cat. No.
Control knob	SD160/DCK
Rubber feet, set of 4	SB160/RF
Retort rod,	SR1
Probe, stainless steel	CD162/1
Probe, PTFE	CD162/2

Warranty

Barloworld Scientific Ltd warrants this instrument to be free from defects in material and workmanship, when used under normal laboratory conditions, for a period of three (3) years. In the event of a justified claim Barloworld Scientific will replace any defective component free of charge.

This warranty does NOT apply if,

1. A ceramic top has broken due to mechanical impact, scratching, chipping or chemical etching.
2. Any repair has been made or attempted other than by Barloworld Scientific or its agent.

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Stone, Staffordshire ST15 0SA
United Kingdom

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Notes

These products meet the relevant EC harmonised standards for radio frequency interference and may be expected not to interfere with, or be affected by, other equipment with similar qualifications. We cannot be sure that other equipment used in their vicinity will meet these standards

and we cannot guarantee that interference will not occur in practice. Where there is a possibility that injury, damage or loss might occur if equipment malfunctions due to radio frequency interference, or for general advice before use, please contact the Technical Service Department of Barloworld Scientific Ltd.

Declaration of Conformity

Digital Hotplate Stirrer

Model SD160, SD162 & CD162

These products comply with the requirements of the EU Directives listed below:

89/336/EEC Electromagnetic Compatibility Directive amended by 93/68/EEC.
73/23/EEC Low Voltage Directive amended by 93/68/EEC

Compliance with the requirements of these Directives is claimed by meeting the following standards:


EN 61326: 1997 + Amendments A1: 1998, EN 61000-4-2: 1995, EN 61000-4-4: 1995, EN 61000-4-5: 1995, EN 61000-4-6: 1996, EN 61000-4-9: 1993, EN 61000-4-11: 1994, EN 61000-3-2: 1995, EN 61000-3-3: 1995, EN 61010-1: 1993 + Amendments A1: 1992 & A2: 1995, EN 61010-2-010: 1994.

Compliance Certificates and Full Reports.

Ref: RETS0265/1, RETS0265/2 and RETS0265/3

From "Epsilon Technical Services" an Independent Accredited Test House showing compliance to the above Standards, are available on request.

CE Mark Affixed '02.

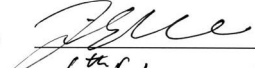Signed:  (Mr D. E. Hicks)

Date: 6th February 2006

Authority: Technical/Development Manager

For

Barloworld Scientific Ltd - Stone - Staffs - ST15 0SA - UK
Tel: +44 (0) 1785 812121 - Fax +44 (0) 1785 813748

INSPECTION REPORT

MODEL _____


ELECTRICAL SAFETY

- | | |
|---------------------|-------------------------------------|
| 1. Earth continuity | ☒ |
| 2. Insulation | ☒ |
| 3. Flash test | ☒ |

FUNCTIONAL

- | | |
|---------------------------|-------------------------------------|
| 1. Indicators | ☒ |
| 2. Heater/stirrer control | ☒ |
| 3. Visual acceptance | ☒ |

QUALITY CONTROL INSPECTOR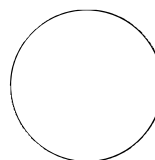

**Barloworld
Scientific**

Barloworld Scientific France SAS
ZI du Rocher Vert - BP 79
77793 Nemours Cedex
France
Tel: +33 1 64 45 13 13
Fax: +33 1 64 45 13 00
e-mail: bibby@bibby-sterilin.fr

Barloworld Scientific Italia Srl
Via Alcide de Gasperi 56
20077 Riozzo di Cerro al Lambro
Milano Italia
Tel: +39 (0)2 98230679
Fax: +39 (0)2 98230211
e-mail: marketing@barloworld-scientific.it
www.barloworld-scientific.it

Barloworld Scientific Ltd
Beacon Road Stone
Staffordshire ST15 0SA
United Kingdom
Tel: +44 (0)1785 812121
Fax: +44 (0)1785 813748
www.barloworld-scientific.com

Barloworld Scientific US Ltd
350 Commerce Drive
Rochester NY 14623
USA
Tel: (800) 828-6595
Fax: (585) 334-0241
www.dynalabcorp.com

Afora S.A.
Calle Aribau 240
08006 Barcelona Spain
Tel: +343 93-306 98 00
Fax: +343 93-306 98 23
e-mail: marketing@afora.com
www.afora.com