

DigiGenius Quick Start Guide



This guide explains basic operations for first-time use, how to use the direct print function and how to install the software and download images to a computer. For more information about usage precautions and detailed instructions, please refer to the other related manuals (page 2).

You must install the software bundled on the Canon Digital Camera Solution Disk before connecting the camera to the computer.

Please refer to the Troubleshooting section of the Software Starter Guide if you have trouble installing the software or downloading images.

Before using the DigiGenius to capture gel images it is advisable to familiarize yourself with the camera and its controls. Please review this complete manual before starting work.

Content

For detailed instructions, see the other manuals below.



Camera User
Guide

Preparing the camera	3
Shooting	4
Reviewing an image	5
Selecting menus and settings	6
List of menu options	7

Setting up the DigiGenius — for UV illuminated DNA gels	15
--	----

Setting up the DigiGenius — for white light illuminated protein gels ...	18
---	----

Printing images by connecting the camera directly to a printer	21
---	----

Printer User Guide
(included with your
Canon printer)


Downloading images to a computer (Windows)	24
---	----

Downloading images to a computer (Macintosh)	31
---	----


Printing to a thermal printer	36
-------------------------------------	----

Transfer of images to GeneTools for analysis	37
---	----

Preparing the Camera

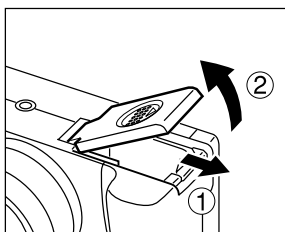


1 Slide the battery cover in the direction of the arrows.

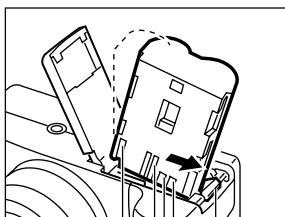

Terminals
Battery lock


2 Press the battery lock while inserting the Battery Pack BP-511 and close the battery cover.

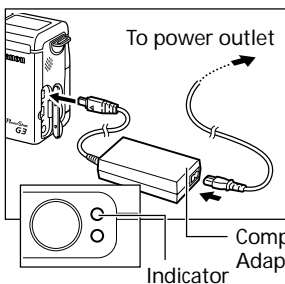


Compact Power Adapter CA-560
Indicator

3 Charge the battery.

- The battery is approximately 90% charged (adequate charge for use) when the orange indicator beside the viewfinder stops blinking and lights steadily.
- After charging, detach the compact power adapter's DC plug from the camera.

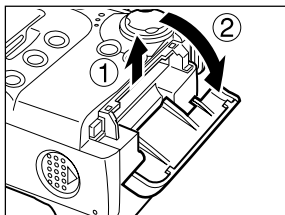

4 Slide the **CF** card slot cover in the direction of the arrow and lift it open.

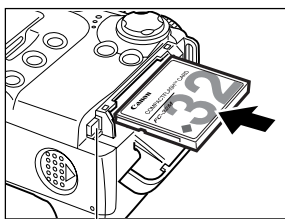


CF card eject button


5 Install the CF card and close the **CF** card slot cover.


- Insert the CF card with the label facing out and push the card in until the CF card eject button fully extends.

Shooting

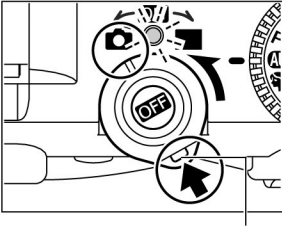


1 Remove the lens cap and turn the power on.

- Open the LCD monitor and turn the mode lever to  (shooting) while pressing the release lever.

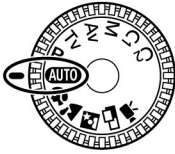


2 Turn the mode dial to **AUTO**.

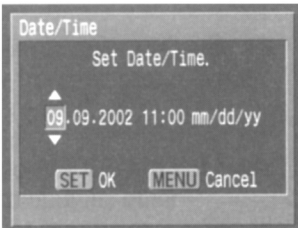


3 Set the date and time. (First time the camera is powered on)


- The Set Date/Time menu will appear on the LCD monitor.

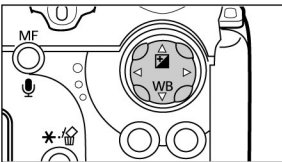

- Select the items you wish to set using the ◀ or ▶ arrow on the omni selector and set the value using the ▲ or ▼ arrow on the omni selector.

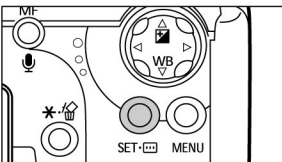


- To put the settings into effect, press the **SET** button.

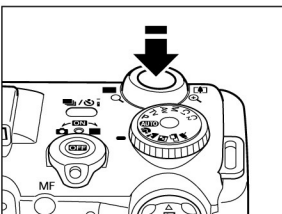

4 Focus.

- Aim the camera and lightly press the shutter button halfway. Two beeps will sound when the autofocus is set.

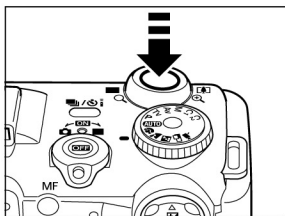


5 Shoot.

- Press the shutter button all the way. You will hear the sound of the shutter close when the shot is complete.
- The recorded image appears in the LCD monitor for approximately 2 seconds.


Reviewing an Image

Displays the recorded images on the LCD monitor.

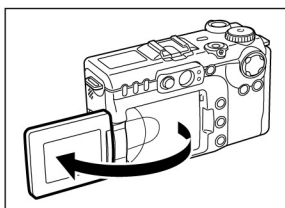


1 Open the LCD monitor.

- The LCD monitor can be freely opened to various position.

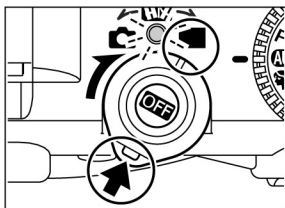


2 Turn the mode lever to .

- The camera powers up in Replay mode.
- The power/mode indicator lights green.
- The last recorded image appears.

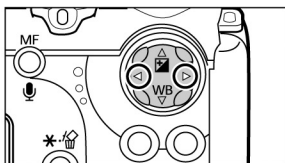

3 Move between images using the or arrow on the omni selector.

To Immediately Delete the Displayed Image

1. Press the  button while the image is displayed.
2. Confirm that [Erase] is selected and press the **SET** button.



Selecting Menus and Settings


- 1 Press the **MENU** button.
 - In Shooting mode, the [RECORD] (Rec.) menu appears.
 - In Replay mode, the [REPLAY] (Replay) menu appears.



- 2 Switch between menu tabs using the ◀ or ▶ arrow on the omni selector.



- 3 Select a menu item using the ▲ or ▼ arrow on the omni selector.

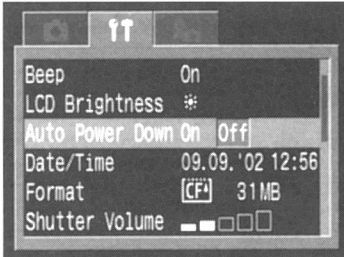

- 4 Select setting contents using the ◀ or ▶ arrow on the omni selector.
 - Press the **SET** button for items followed by an ellipsis (...) then select a setting. Press the **SET** button again to confirm the setting.

- 5 Press the **MENU** button to close the menu.







List of Menu Options





Rec. Menu

Some items may not appear in certain Shooting modes.


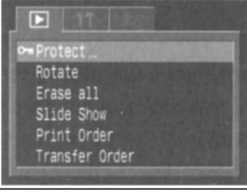

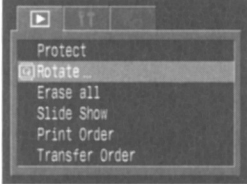
*Default setting


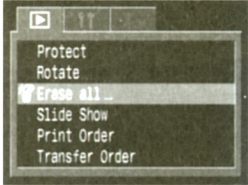

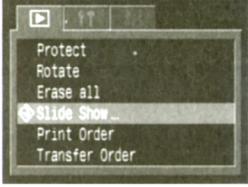

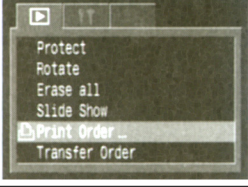

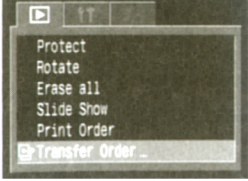
Item	Menu Screen	Settings Content
Flash Sync		Sets the timing at which the flash fires. <ul style="list-style-type: none"> • 1st-curtain* • 2nd-curtain
Slow Synchro		Sets the slow synchro. <ul style="list-style-type: none"> • On • Off*
Flash Adjust		Sets whether or not the flash is adjusted automatically. <ul style="list-style-type: none"> • Auto* • Manual
Red-eye		Sets whether or not the red-eye reduction lamp fires. <ul style="list-style-type: none"> • On* • Off
Cont. Shooting		Sets the continuous shooting option. <ul style="list-style-type: none"> •  (Standard)* •  (High speed)
Self-timer		Sets the duration before the self-timer releases the shutter once the shutter button is pressed. <ul style="list-style-type: none"> • 2 seconds • 10 seconds*

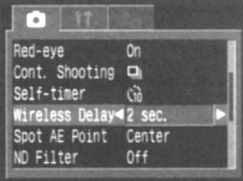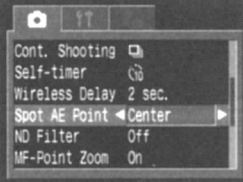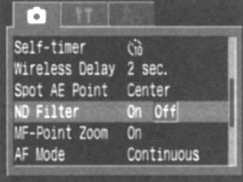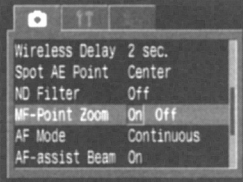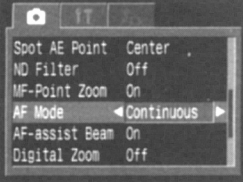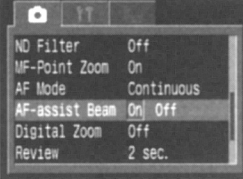Item	Menu Screen	Settings Content
Wireless Delay		Sets the duration before the self-timer releases the shutter once the remote control shutter button is pressed. <ul style="list-style-type: none"> • 0 seconds • 2 seconds* • 10 seconds
Spot AE Point		Sets the Spot AE Point option, Center (the AE point is fixed to the center) or AF Point (the AE point corresponds to the AF point). <ul style="list-style-type: none"> • Center* • AF Point
ND Filter		Sets the ND filter. <ul style="list-style-type: none"> • On • Off*
MF-Point Zoom		Sets whether or not the focusing point is magnified when using the manual focus. <ul style="list-style-type: none"> • On* • Off
AF Mode		Sets the frequency of autofocus activation. When the LCD monitor is off, this setting locks in single mode. <ul style="list-style-type: none"> • Continuous* • Single
AF-assist Beam		Sets the output of the AF-assist beam according to shooting conditions when using the auto focus. <ul style="list-style-type: none"> • On* • Off

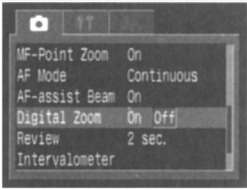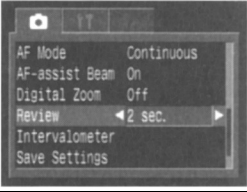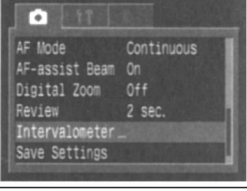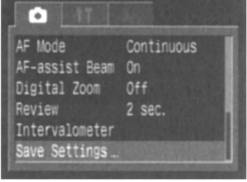



Item	Menu Screen	Settings Content
Digital Zoom		Sets whether or not the digital zoom combines with the optical zoom to zoom images. <ul style="list-style-type: none"> • On • Off*
Review		Sets the duration images display on the LCD monitor after the shutter button is released. <ul style="list-style-type: none"> • Off • 2 sec.* to 10 sec.
Intervalometer		Shoots automatically at set intervals.
Save Settings		Saves settings selected in the shooting menu and the function menu to the mode dial's C1 or C2 mode.




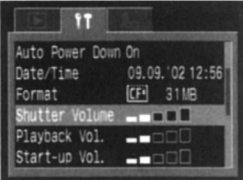
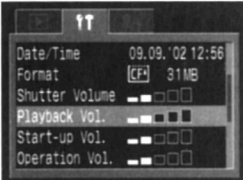
Replay Menu













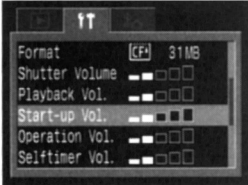






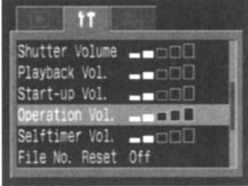




Item	Menu Screen	Settings Content
 Protect		Protects images from accidental erasure.
 Rotate		Rotates images 90 or 270 degrees clockwise in the display.




Item	Menu Screen	Settings Content
 Erase all		Erases all images from a CF card (excluding protected images).
 Slide Show		Plays images in an automated slide show.
 Print Order		Sets the images that will be printed on a DPOF-compatible printer or at a photo lab service, the number of copies and other parameters.
 Transfer Order		Sets the images that will be transferred with your e-mail software.

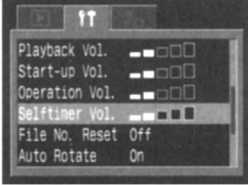Set up Menu






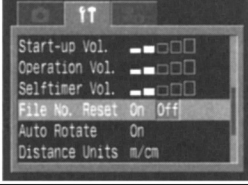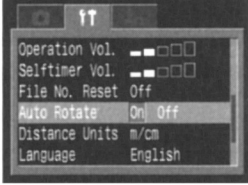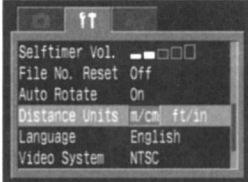


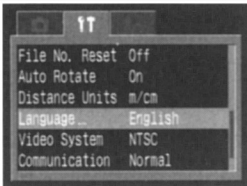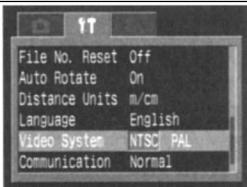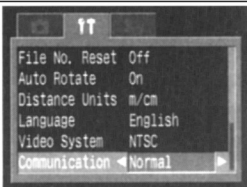

Item	Menu Screen	Settings Content
Beep		Sets whether or not the beep sounds when the shutter button is pressed or when a menu operation is performed. <ul style="list-style-type: none"> • On* • Off (Still beeps to warn of an error even when set to Off)
LCD Brightness		Sets the brightness level of the LCD monitor. <ul style="list-style-type: none"> •  (Normal)* •  (Bright)

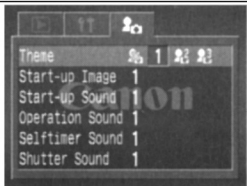



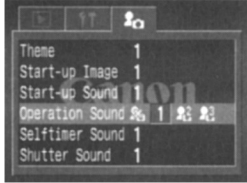


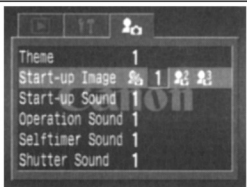
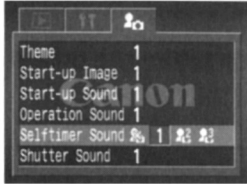



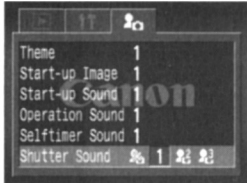


Item	Menu Screen	Settings Content
Auto Power Down		<p>Sets whether or not the camera automatically powers down after a specified period of time elapses without something being pressed.</p> <ul style="list-style-type: none"> • On* • Off
Date/Time		<p>Sets the date, time and date format.</p>
Format		<p>Formats (initializes) a CF card.</p>
Shutter Volume		<p>Adjusts the sound volume when the shutter button is pressed all the way. There is no shutter sound in Movie mode.</p> <ul style="list-style-type: none"> • □□□□ (Off) • ■□□□ (1) • ■■□□ (2)* • ■■■□□ (3) • ■■■■□□ (4) • ■■■■■ (5)
Playback Vol. (Replay)		<p>Adjusts the sound volume when a sound annotation or a movie is played.</p> <ul style="list-style-type: none"> • □□□□ (Off) • ■□□□ (1) • ■■□□ (2)* • ■■■□□ (3) • ■■■■□□ (4) • ■■■■■ (5)

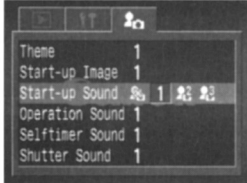Item	Menu Screen	Settings Content
Start-up Vol.		<p>Adjusts the sound volume when the camera powers up.</p> <ul style="list-style-type: none"> •  (Off) •  (1) •  (2)* •  (3) •  (4) •  (5)
Operation Vol.		<p>Adjusts the sound volume of the Operation Sound.</p> <ul style="list-style-type: none"> •  (Off) •  (1) •  (2)* •  (3) •  (4) •  (5)
Selftimer Vol.		<p>Adjusts the sound volume of the Selftimer Sound.</p> <ul style="list-style-type: none"> •  (Off) •  (1) •  (2)* •  (3) •  (4) •  (5)
File No. Reset		<p>Sets how file numbers are assigned to images when new CF cards are inserted.</p> <ul style="list-style-type: none"> • On • Off*
Auto Rotate		<p>Sets whether or not images shot with the camera on end are automatically rotated in the display.</p> <ul style="list-style-type: none"> • On* • Off
Distance Units		<p>Sets the format of the MF indicator's displayed distance units.</p> <ul style="list-style-type: none"> • m / cm* • ft / in

Item	Menu Screen	Settings Content
Language		<p>Sets the language used in the menus and messages on the LCD monitor. The menu may appear differently in some regions.</p> <ul style="list-style-type: none"> • English* • Deutsch • Français • Nederlands • Dansk • Suomi • Italiano • Norsk • Svenska • Español • Chinese • Japanese <p>You can change the language during image replay by holding the SET button and pressing the JUMP button. (Except when the camera is connected to a printer)</p>
Video System		<p>Sets the video output signal standard.</p> <ul style="list-style-type: none"> • NTSC • PAL
Communication		<p>Sets the camera and computer communication settings.</p> <ul style="list-style-type: none"> • Normal* • PTP

My Camera Menu

Item	Menu Screen	Settings Content
Theme		<p>Selects a common theme for each My Camera settings item.</p> <ul style="list-style-type: none"> •  (Off) •  * •  • 
Start-up Image		<p>Sets the start-up image when you turn on the camera.</p> <ul style="list-style-type: none"> •  (Off) •  * •  • 

Item	Menu Screen	Settings Content
Start-up Sound		<p>Sets the start-up sound when you turn on the camera.</p> <ul style="list-style-type: none"> •  (Off) • 1 * •  • 
Operation Sound		<p>Sets the sound when the multicontroller or any button other than the shutter button is operated.</p> <ul style="list-style-type: none"> •  (Off) • 1 * •  • 
Selftimer Sound		<p>Sets the sound informing you that the photo will be taken in 2 seconds.</p> <ul style="list-style-type: none"> •  (Off) • 1 * •  • 
Shutter Sound		<p>Sets the shutter sound when you depress the shutter button. There is no shutter sound while a movie is being shot.</p> <ul style="list-style-type: none"> •  (Off) • 1 * •  • 