Prime and PrimeG

OPERATOR'S MANUAL

CONTENTS

	PAGE
INTRODUCTION	3
BEFORE USE	4
Unpacking	4
SAFETY INFORMATION	5
English	5
Français	7
Deutsch	9
Italiano	11
Español	13
CONTACT INFORMATION	15
PRIME SPECIFICATION	16
INSTALLATION	18
GENERAL SOFTWARE FEATURES	20
Introduction	20
Template programs	20
Incremented/decremented time and temperature	21
Gradient (PrimeG only)	21
Oligonucleotide T_m calculator	22
PROGRAMMING TEXT, TEMPERATURE AND TIME	23
Text entry	23
Temperature entry	23
Time entry	27
INSTRUMENT SETTINGS AND DEFAULTS	30
User preferences	30
Instrument settings	31
Program defaults	33
Service settings	34
Gradient calculator	35
Oligonucleotide T_m calculator	35
CREATE A NEW PROGRAM	37
Name the program	37
Insert/delete a stage	38
Insert/delete a step	39
Insert/delete a pause	39
Program options	39


	PAGE
Settings	43
View	43
Delete	44
Up and down arrows	44
Save	44
Cancel	44
RUN A PROGRAM	45
Running a program from the shortcut list	45
The pre-run screen	45
The run screen	46
Searching for a program not in the shortcut list	47
Running a program not on the shortcut list	49
EDITING A PROGRAM	50
Editing saved programs	50
Editing templates	50
Editing password-protected programs	50
CURRENT STATUS	51
VIEW FILES	52
File searching	52
Viewing a program file	54
Viewing a temperature log file	54
Copying files	55
Deleting files	57
UPDATING THE PRIME SOFTWARE	58
REMOVING AND EXCHANGING BLOCKS	59
UPGRADING FROM PRIME TO PRIMEG	59
FAQs	60
TECHNICAL SUPPORT AND SERVICING	61
ADDITIONAL INFORMATION	62
User Maintenance	62
Fault Finding	62
Fuses	62
Insulation Testing	62
ACCESSORIES	63
REPLACEMENT PARTS	63

INTRODUCTION

The Prime is a stand-alone, flexible and versatile thermal cycler designed to accurately control the temperatures of samples between 4°C and 100°C. It has many scientific applications including PCR¹, reverse transcription, ligation and sequencing. It has user-interchangeable blocks to allow for most commonly used consumables including 0.2ml x 96 well plates, 0.2ml tubes, 0.5ml tubes and 384 well plates. The heated lid is fully adjustable to accommodate a wide range of different consumables.

The Prime is programmed via a 5.7" colour VGA LCD touch screen using simple and intuitive software. Programs and run progress can be viewed as a temperature/step graphical display, allowing the user to readily check the status or position within a run. Files can be saved with user names and/or keywords which allows for rapid searching and retrieval. A USB port on the front of the unit allows transfer of files to and from a USB memory stick for viewing in the Techne Workbench PC software and for the easy update of instrument software.

The Prime can easily be upgraded to the PrimeG which incorporates a block gradient function and gradient calculator, allowing for simple and rapid optimisation of reaction conditions.

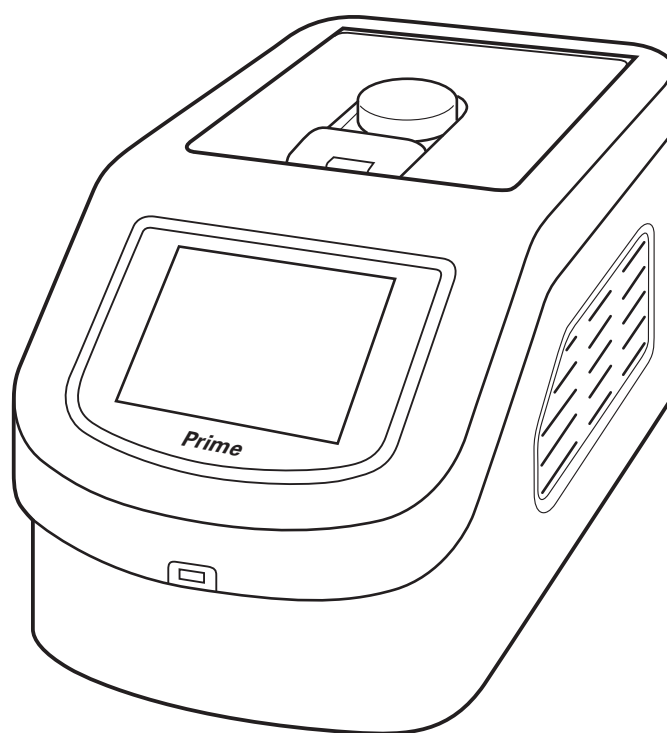¹Purchase of this instrument conveys a limited non-transferable immunity from suit for the purchaser's own internal research and development and applied fields other than human in-vitro diagnostics under non-real-time thermal cycler patents of Applied Biosystems LLC.

BEFORE USE

Before using the Prime please ensure you have read this manual carefully. If there is any doubt relating to the proper use of this equipment, the staff at Bibby Scientific Ltd. or your supplier will be happy to assist you.

UNPACKING

When unpacking the unit please ensure that the following have been removed from the packaging:

- Prime with block
- Mains cables (UK, EU and US)
- Operator's manual
- Warranty card

The user is advised to keep the original packaging in case the instrument ever needs to be returned for service or repair. Bibby Scientific Ltd. accepts no responsibility for damage incurred unless the unit is correctly packed and transported in its original packaging.

SAFETY INFORMATION

Please read all the information in this manual before using the Prime.

WARNING

HIGH TEMPERATURES ARE DANGEROUS: they can cause serious burns to operators and ignite combustible material. Users should be aware of the following potential hazards:

- USE CARE AND WEAR PROTECTIVE GLOVES TO PROTECT HANDS
- DO NOT use combustible substances near hot objects
- DO NOT operate the instrument in the vicinity of inflammable liquids or gases
- DO NOT place any liquid directly into the instrument.

OPERATOR SAFETY

All operators of Techne equipment must have available the relevant literature needed to ensure their safety. It is important that only suitably trained personnel operate this equipment, in accordance with the instructions contained in this manual and with general safety standards and procedures. If the equipment is used in a manner not specified by Bibby Scientific Ltd. the protection provided by the equipment to the operator may be impaired.

All Techne instruments are designed to conform to international safety requirements and are fitted with an over-temperature cut-out. If a safety problem should be encountered, switch off the unit at the mains socket and remove the plug from the electricity supply.

INSTALLATION

The instrument should be carried using both hands. Never move or carry the instrument when in use or connected to the mains electricity supply.

1. All Techne instruments are supplied with a power cable; this may be integral or plug-in.

2.  Before connecting the instrument to the mains electricity supply, check the voltage against the rating plate (located on the back of the unit). **Please note that the unit must be earthed to ensure proper electrical safety.** Connect the mains cable to a suitable plug according to the table below.

Connections	220/240V, 50/60Hz	110V/120V, 50/60Hz
Live	Brown	Black
Neutral	Blue	White
Earth	Green/yellow	Green

3. UK ONLY: The fused plug supplied with the mains cable is fitted with a 5 amp fuse to protect the instrument and the operator.
4. The units are rated to operate at 100-240V, 50/60Hz.
5. Place the unit on a suitable flat bench or in a fume cupboard if required, ensuring that the air inlet vents on the underside are free from obstruction.
6. Plug the mains cable into the socket on the back of the instrument.
7. Switch on the instrument using the switch located on the front of the unit.

Replacement cable

Should the mains lead need replacement, a cable of 1mm² of harmonized code H05VV-F connected to an IEC320 plug should be used. **IF IN DOUBT CONSULT A QUALIFIED ELECTRICIAN.**

ENVIRONMENTAL CONDITIONS

The Prime is designed operate under the following conditions:

- Indoor use
- Ambient temperature range +5°C to +40°C
- Altitude to 2000m
- Relative humidity not exceeding 80%
- Mains supply fluctuations not exceeding 10%
- Over voltage category II IEC 60364-4-443
- Pollution degree 2

Note: The control specifications are quoted at an ambient temperature of 20°C. The specification may deteriorate outside an ambient temperature of between 10°C and 30°C.

The instrument has been tested for radio frequency interference and is certified under EN61326.

GUARANTEE

The instrument is guaranteed against any defect in material or workmanship for the period specified on the enclosed guarantee card. This period is effective from the date of purchase; within this period all defective parts will be replaced free of charge provided that the defect is not the result of an accident, misuse or negligence. Servicing under this guarantee should be obtained from the supplier of the instrument.

This manual has been prepared for the convenience of Techne's customers and nothing in this manual shall be taken as a warranty, condition or representation concerning the description, merchantability, fitness for purpose or otherwise of the unit or components.

Notwithstanding the description and specification(s) of the instruments contained in the operator's manual, Techne reserves the right to make such changes as it sees fit to the instruments or to any of the components.

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CONTACT INFORMATION

For technical, sales or servicing information, contact your local Techne dealer or:

UK

Bibby Scientific Ltd.
Beacon Road
Stone
Staffordshire
ST15 0SA
UK
Tel: +44 (0)1785 812121
Fax: +44 (0)1785 810405
E-mail: sales@bibby-scientific.com
www.techne.com

France

Bibby Scientific Limited
Bâtiment Le Deltaparc Parc Silic PN2
7 rue du Canal
BP 55437 VILLEPINTE
95944 ROISSY Charles de Gaulle
France
Tel: +33(0)148 63 78 03
Fax: +33(0)148 63 78 01
E-mail: ventes@bibby-scientific.com
www.bibby-scientific.com

Middle East

Bibby Scientific Middle East Ltd.
PO Box 27887, Engomi 2433
Nicosia
Cyprus
Tel: +357 22 660 423
Fax: +357 22 660 424
e-mail: sales@bibby-scientificme.com

North and South America

Bibby Scientific US Inc. t/a Techne Inc.
3 Terri Lane, Suite 10
Burlington, N.J. 08016
USA
Toll Free (in NA): 800-225-9243
Tel: +1 609 589 2560
Fax: +1 609 589 2571
E-mail: labproducts@techneusa.com
www.techneusa.com

Italy

Bibby Scientific Italia Srl
Via Alcide de Gasperi 56
20070 Riozzo di Cerro al Lambro (MI)
Italia
Tel: +39 (0)2 98230679
Fax: +39 (0)2 98230211
E-mail: marketing@bibby-scientific.it
www.bibby-scientific.it

Asia

Bibby Scientific - Singapore
Prudential Tower, Level 26
30 Cecil Street
Singapore 049712
Tel: +65 6631 2976
Fax: +44 (0) 1785 810405
e-mail: info@bibby-scientific.com
www.bibby-scientific.com

PRIME SPECIFICATION

TEMPERATURE

Block temperature range	4°C to 100°C
Block uniformity (at 55°C)	<±0.3°C
Temperature accuracy (at 55°C)	<±0.25°C
Gradient availability (PrimeG only)	Between 30°C and 80°C
Maximum gradient (PrimeG only)	29°C (19°C with the 384-well block)
Minimum gradient (PrimeG only)	1°C
Column uniformity with a 10°C gradient (PrimeG only)	<±0.3°C
Pre-run sample cooling	Yes, 4°C
Temperature set point adjustment	0.1°C

HEATING/COOLING RATE

Maximum heating rate	3.4°C/s
Minimum heating rate	0.1°C/s
Maximum cooling rate	1.0°C/s
Minimum cooling rate	0.1°C/s

HEATED LID

Selectable heated lid temperature	35 to 115°C or off
Lid uniformity	±5°C
Pre-heat lid	Yes
Warm up time from ambient	<2 min
Over-temperature cut-out	Yes
Regulated heated lid pressure	Manually adjustable

The heated lid is only operational if the block temperature is set above 35°C

PROGRAMMING

Program interface	5.7" colour VGA LCD touch screen
Maximum number of programs stored	1000
Maximum number of stages per program	10
Maximum number of steps	10
Maximum number of cycles per stage	99
Programmable ramp rate	Yes, 0.1°C/s steps
Maximum hold time	4h 59m 59s
Minimum hold time	1s
Gradient calculator (PrimeG only)	Yes
Incremented/decremented temperature	Yes
Incremented/decremented time	Yes
Pause facility	Yes
Program naming	Alpha numeric plus symbols
Password protection	Yes
Keyword selection	Yes
Run completion time	Yes
End of program alarm	Yes
Auto resume on power failure	Yes, always or never
Oligonucleotide T _m calculator	Yes, based on the Nearest-Neighbour method ¹
Software updates	Free of charge from www.techne.com

¹Breslauer, K.J.; Frank, R.; Blocker, H. and Marky, L.A. (1986) Proc. Natl. Acad. Sci. USA 83, pp 3746-3750.

COMMUNICATION PORTS

USB port on front of unit	Connection for USB memory stick
USB port on rear of unit	Connection for servicing and calibration only

DIMENSIONS

Height (mm)	240
Width (mm)	240
Length (mm)	420
Weight (kg)	9.4

POWER


Voltage	100-240V, 50/60Hz
Power	425W

INSTALLATION

See also the **Safety Information** sections on pages 5 to 14.


FRONT VIEW

1. Lid adjustment knob.
2. Lid release latch.
3. 5.7" colour VGA LCD touch screen display.
4. USB port for memory stick.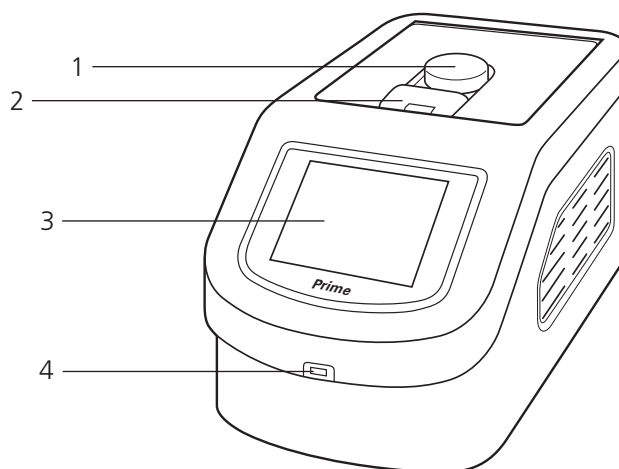


REAR VIEW

1. Mains cable inlet.
2. On/Off rocker switch
3. USB port for service and calibration only.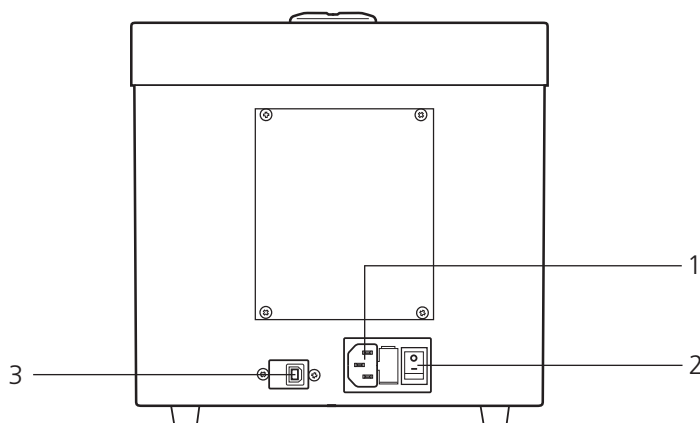


INSTALLATION


1. Place the unit on a suitable flat bench, ensuring that the air inlet vents on the sides are free from obstruction.
2. Insert the block into the unit ensuring the connectors on the block and unit are aligned.
3. Slide the quick release bars on the block towards the closed padlock symbols to lock the block in place.
4. Plug the mains cable into the socket on the back of the instrument.
5. Connect to the mains electricity supply with the plug provided or one wired correctly for the supply. Switch the power ON using the switch located on the rear of the unit. The front display will then light up.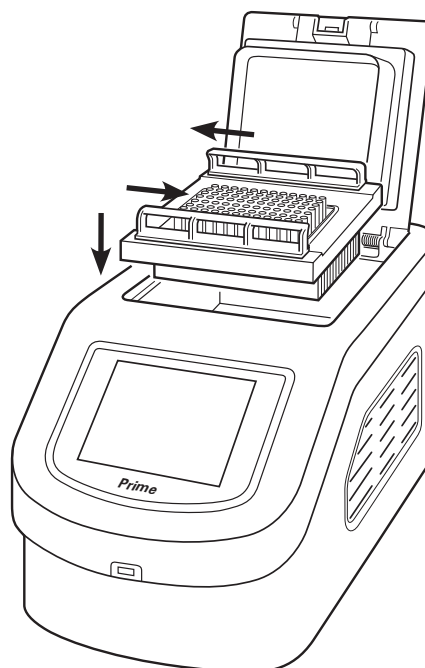

6. Release the heated lid by lifting the latch forward and use it to lift the lid. To close the lid, lower the lid onto the block and press the latch to secure in place.
7. Place the samples in the block. If individual tubes or strip tubes are being used, space these out evenly across the block to equalise the pressure from the heated lid.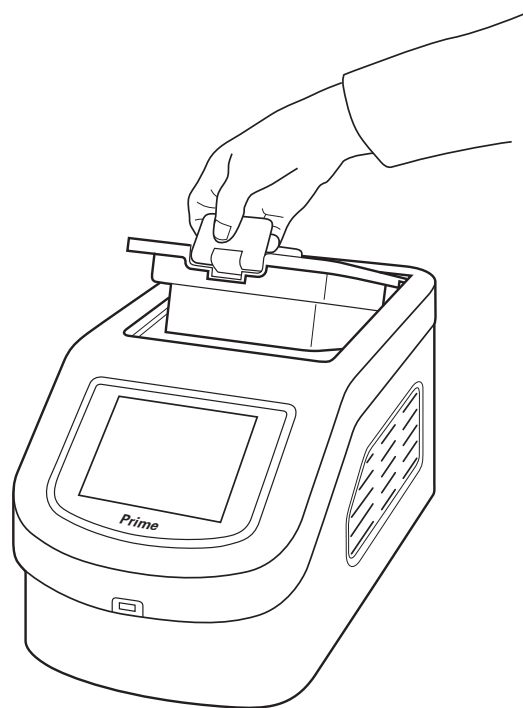

8. The heated lid has a rotating knob to adjust the lid pressure on the samples, allowing for a variety of consumables to be used. To adjust the pressure:
 - a. Place the samples in the block and place the lid in the down position.
 - b. Adjust the orange rotating knob anticlockwise to give the least pressure then close the lid fully.
 - c. Gently rotate the knob clockwise until you can feel pressure beginning to be applied.
 - d. Rotate the knob a further quarter of a turn and this will give the correct pressure. Do not over-tighten.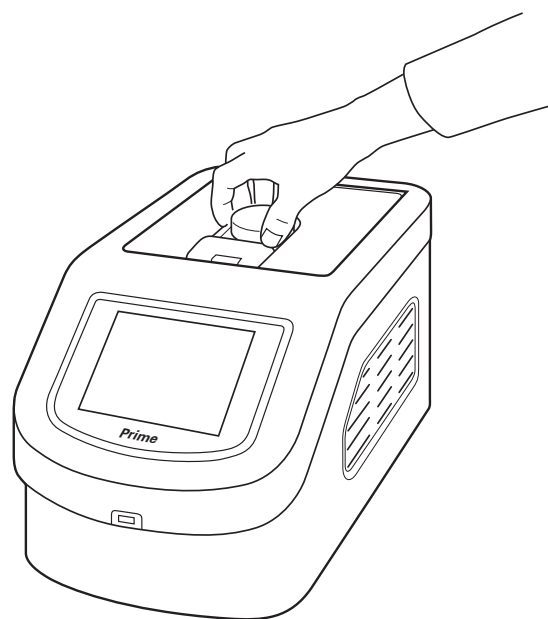


9. Once the pressure has been set it does not need to be adjusted unless a different block or consumable is used.

GENERAL SOFTWARE FEATURES

INTRODUCTION

The Prime software allows the user to quickly and easily create thermal cycling programs using a simple and clear touch screen format. The software is structured into five modules which are accessed from the Home Screen. These allow the user to:

- Run an existing program, selecting from templates or programs saved on the unit.
- Create a new program.
- Set instrument and program defaults and preferences. Includes calculator tools.
- View the current status of the unit.
- View and search for files stored on the unit.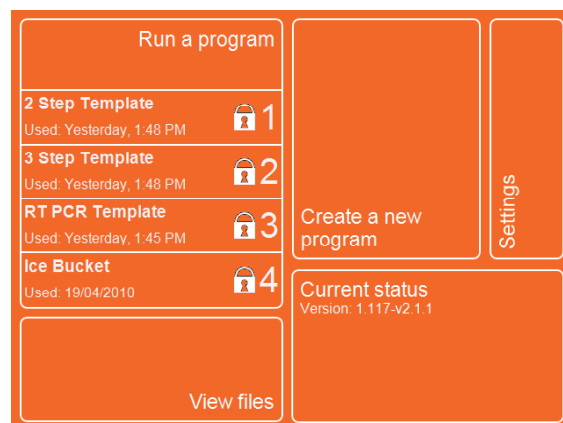

TEMPLATE PROGRAMS

To help with programming, a number of program templates are installed on the unit which are available to copy and edit or can be run directly without changes. Details of the installed program templates are given below.

Instrument and program defaults

Parameter	2 Step Template	3 Step Template	RT-PCR Template
Heated Lid	105°C	105°C	105°C
Heated lid before program	On	On	On
Sample cooling	On	On	On
Pause before program	Off	Off	Off
Initial denaturation	94°C, 05m00s	94°C, 05m00s	94°C, 05m00s
Hot start	Off	Off	Off
Final extension	72°C, 05m00s	72°C, 05m00s	72°C, 07m00s
Final hold	10°C	10°C	10°C

Thermal cycling conditions

	2 Step Template	3 Step Template	RT-PCR Template
Stage 1			
Number of cycles	30	30	1
Step 1	94.0°C, 00m30s	94.0°C, 00m30s	45.0°C, 40m00s
Step 2	60.0°C, 01m00s	55.0°C, 00m30s	94.0°C, 05m00s
Step 3		72.0°C, 00m30s	
Stage 2			
Number of cycles			40
Step 1			94.0°C, 01m00s
Step 2			55.0°C, 00m50s
Step 3			72.0°C, 01m00s


The Ice Bucket and Ligation programs are simple temperature holds, with the former having an infinite hold at 10°C and the latter at 15°C. In both of these programs the heated lid is switched off and the initial denaturation, hot start and final extension functions disabled.

INCREMENTED/DECREMENTED TIME AND TEMPERATURE

The software allows the user to either increment or decrement the time and/or temperature of a step within a cycling stage. These features are used with applications such as touchdown PCR where the annealing temperature is gradually decreased during the cycling process and long PCR, where due to the size of the product, long annealing/extension times are required and the extension time is increased by 15 to 20s per cycle during the final stages of the reaction.

GRADIENT (PRIMEG ONLY)

The gradient feature of the PrimeG can be useful in optimising the annealing conditions for primers. A gradient can be set around a temperature in any step of a program. The set temperature is the temperature in the middle columns of the block and the gradient is the variation at the two extremes of the block; the left hand columns being the coolest and the right hand columns the hottest.

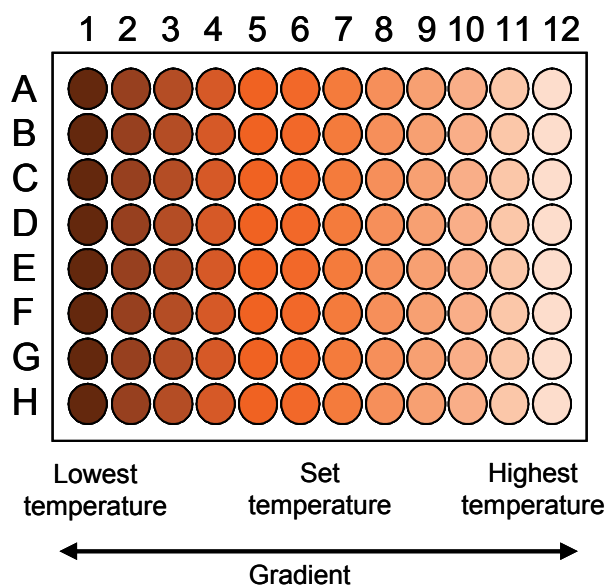The maximum temperature gradient range which can be set is 29°C and the minimum is 1°C; the lowest temperature is 30°C and the hottest temperature is 80°C. Examples are given in the table below:

Lowest temperature (°C)	Set temperature (°C)	Highest temperature (°C)	Gradient (°C)
50	55	60	10
40.5	55	69.5	29
50.5	65	79.5	29

A gradient calculator is included in the **Settings** module so the user can obtain an accurate estimation of the temperature of each column during the gradient stage. The temperatures for any particular gradient are calculated from an algorithm stored in the block firmware.

OLIGONUCLEOTIDE T_m CALCULATOR

One of the most important parameters in a PCR is the annealing temperature of the reaction. This can be determined theoretically using various software programmes to calculate the melting temperature (T_m) of the primers.

The T_m is defined as the temperature at which half the DNA strands are single stranded and half are double stranded when base paired to a complementary strand. Using annealing temperatures that are well below the T_m of the primers can result in mismatching, false priming and may lead to primer-dimer artefacts. Annealing temperatures that are significantly higher than the primer T_m may result in reduced priming or prevent priming altogether.

The Prime uses the Nearest-Neighbour method for calculating the T_m of an oligonucleotide¹.

The T_m is calculated using the following equation:

$$T_m = \{(\Delta H^\circ \times 1000) / (A + \Delta S^\circ + R \ln(C/4))\} - 273.15 + 16.6 \times \log[Na^+]$$

Where:

- ΔH° = The sum of the Nearest-Neighbour enthalpy changes (Kcal/mol)
- A = Helix initiation correction constant
- ΔS° = The sum of the Nearest-Neighbour entropy changes (cal K⁻¹mol⁻¹)
- R = The Gas Constant (1.99 cal K⁻¹mol⁻¹)
- C = Concentration of the oligo (M)
- [Na+] = Concentration of monovalent ions (M)

¹Breslauer, K.J.; Frank, R.; Blocker, H. and Marky, L.A. (1986) Proc. Natl. Acad. Sci. USA 83, pp 3746-3750.

The Oligonucleotide T_m calculator can be found in the **Settings** module where the user can input the oligonucleotide sequence, its concentration in the reaction and the salt concentration of the reaction mix.

PROGRAMMING TEXT, TEMPERATURE AND TIME

This section gives instructions for the basic programming of text, temperature and time which are required in later sections.

TEXT ENTRY

When giving a name to a program or stage, or when entering a keyword or a password, a text entry screen will be opened.

Touch individual characters to create the name.

- For words with all upper case letters touch **CAPS**.
- Touch **SHIFT** if only the first letter is required in upper case.
- For numbers touch **123**.
- For commonly used symbols touch **SYM**.
- Touch the **space bar** (not labelled) for spaces between characters.
- To clear the typed name, touch **CLR**.
- To delete the last character touch **DEL**.

Once the name is complete, touch **OK**.


TEMPERATURE ENTRY

When touching a temperature parameter a temperature entry screen will open. There are two forms of temperature entry screen:

- A **simple** form for adjusting the temperature of the instrument parameters such as the heated lid and program parameters such as initial denaturation and final extension.
- An **advanced** form when programming step temperatures within a cycling stage that includes the options for introducing an incremented/decremented step temperature, for example in touchdown PCR, or a temperature gradient.

Simple temperature entry

- Enter the temperature, type in the required value. The decimal point can be used for 0.1°C resolution.
- To clear the entry press **DEL**.
- If an entered temperature is invalid, a prompt will indicate the acceptable range.
- Once the temperature is correct, touch **OK** to return to the previous screen.
- Touch **Cancel** to return to the previous screen if no changes are required.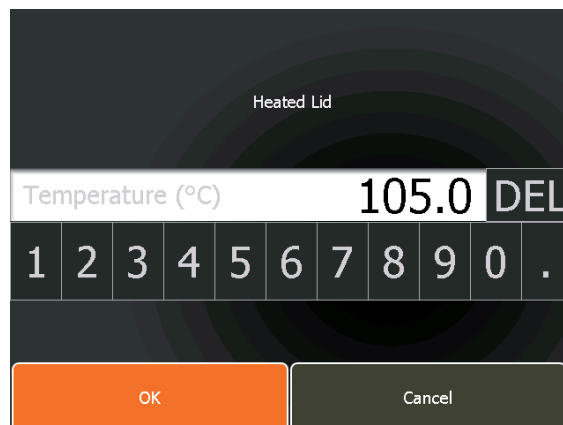

Advanced temperature entry

- First ensure the number of cycles for the stage has been set then enter the required hold temperature as described above.

Note: Increase/Decrease Per Cycle will be inactive unless the number of cycles set is greater than 1.

To increase the step temperature per cycle:

- For an **increase** in temperature per cycle, touch the button below **Increase Per Cycle** to turn this function on. Next touch the temperature entry button below this.

Temperature (°C)										55	DEL
1	2	3	4	5	6	7	8	9	0	.	
Increase Per Cycle					Decrease Per Cycle						
On					Off						
+0.0°C					+0.0°C						
OK					Cancel						

A new screen will open which is similar to the simple temperature entry screen. The maximum increase in temperature per cycle (based on the number of cycles and the temperature limits of the unit) will be shown.

- Enter the required value.
- Touch **DEL** to clear the entry.
- If an entered incremented temperature is invalid, a prompt will indicate the acceptable range.
- Touch **OK** to accept and return to the previous screen.

Increase Per Cycle Maximum increase per cycle is 1.6°C											
Up to 1 decimal place										0.5	DEL
1	2	3	4	5	6	7	8	9	0	.	
OK					Cancel						

The increase in temperature per cycle will now be displayed.

- Touch **OK** again to return to the programming screen.
- Touch **Cancel** to return to the programming screen if no changes are required.

Temperature (°C)										55	DEL
1	2	3	4	5	6	7	8	9	0	.	
Increase Per Cycle					Decrease Per Cycle						
On					Off						
+0.5°C					+0.0°C						
OK					Cancel						

The temperature value will now appear orange in the programming screen indicating that it contains a modified function.

Test #1

Program Options

Settings

Stage 1

35 cycles

View

Delete

Step 1

Temperature

95.0°C

Time

00:01:00

Ramp Rate

Maximum

Step 2

Temperature

55.0°C

Time

00:01:00

Ramp Rate

Maximum

Insert a new step

Save

Cancel

To decrease the step temperature per cycle:

- For a **decrease** in temperature per cycle, touch the button below **Decrease Per Cycle** to turn this function on. Next touch the temperature entry button below this.

Temperature (°C)

55 DEL

1 2 3 4 5 6 7 8 9 0 .

Increase Per Cycle

Decrease Per Cycle

Off

On

+0.0°C

+0.0°C

OK

Cancel

A new screen will open which is similar to the simple temperature entry screen. The maximum decrease in temperature per cycle (based on the number of cycles and the temperature limits of the unit) will be shown.

- Enter the required value.
- Touch **DEL** to clear the entry.
- If an entered decremented temperature is invalid, a prompt will indicate the acceptable range.
- Touch **OK** to return to the previous screen.

Decrease Per Cycle

Maximum decrease per cycle is 1.9°C

Up to 1 decimal place

1.5 DEL

1 2 3 4 5 6 7 8 9 0 .

OK

Cancel

The decrease in temperature per cycle will now be displayed.

- Touch **OK** again to return to the programming screen.
- Touch **Cancel** to return to the programming screen if no changes are required.

Temperature (°C)

55 DEL

1 2 3 4 5 6 7 8 9 0 .

Increase Per Cycle

Decrease Per Cycle

Off

On

+0.0°C

-1.5°C

OK

Cancel

The temperature value will now appear orange in the programming screen indicating that it contains a modified function.

Note: It is not possible to both increase and decrease the temperature of the same step.

To enter a temperature gradient on a step (PrimeG only):

On the PrimeG an additional **Temperature Gradient** button is available.

- Touch the button below **Temperature Gradient** to turn this function on. Next touch the temperature entry button below this.

A new screen will open which is similar to the simple temperature entry screen. The maximum temperature gradient (based on the set hold temperature and the temperature limits of the unit) will be shown.

- Enter the required value.
- Touch **DEL** to clear the entry.
- If an entered gradient temperature is invalid, a prompt will indicate the acceptable range.

- Touch **OK** to return to the previous screen. The temperature gradient for the step will now be displayed.
- Touch **OK** again to return to the programming screen.
- Touch **Cancel** to return to the programming screen if no changes are required.

The temperature value will now appear orange in the programming screen indicating that it contains a modified function.

The screenshot shows the 'Test #1' programming screen. At the top, there are tabs for 'Test #1', 'Program Options', and 'Settings'. Below this, 'Stage 1' is selected, showing '35 cycles'. The screen is divided into two steps. Step 1 has a Temperature of 95.0°C, Time of 00:01:00, and Ramp Rate of Maximum. Step 2 has a Temperature of 55.0°C (highlighted in orange), Time of 00:00:30, and Ramp Rate of Maximum. On the right side, there are buttons for 'View', 'Delete', up and down arrows, 'Save', and 'Cancel'. At the bottom, there is a button labeled 'Insert a new step'.

TIME ENTRY

When touching a time parameter a time entry screen will open. There are two forms of time entry screen:

- A **simple** form for adjusting the time of the program parameters such as initial denaturation and final extension.
- An **advanced** form when programming step hold times within a cycling stage that includes the options for introducing incremented/decremented time, for example in long PCR.

Simple time entry

- Enter the time using the + or – buttons to increase or decrease the number of hours, minutes or seconds.
- Using the – button with the value set at 0 will scroll to the maximum permitted value.
- Touching **Hide Hours** will remove the hour time selector and just display selectors for minutes and seconds.
- Once the time is correct, touch **OK** to return to the previous screen.
- Touch **Cancel** to return to the previous screen if no changes are required.

The screenshot shows the 'Simple time entry' screen. At the top, there is a button labeled 'Hide Hours'. Below it, the text 'Final Extension Time' is displayed twice. The screen features three input fields: 'Hours' with a value of 0, 'Minutes' with a value of 5, and 'Seconds' with a value of 0. Each field has minus and plus buttons for adjustment. At the bottom, there are two buttons: 'OK' (orange) and 'Cancel' (grey).

Advanced time entry

- First ensure the number of cycles for the stage has been set then enter the required hold time as described above.

Note: **Increase/Decrease Per Cycle** will be inactive unless the number of cycles set is greater than 1.

To increase the step hold time per cycle:

- For an **increase** in hold time per cycle, touch the button below **Increase Per Cycle** to turn this function on. Next touch the time entry button below this.

A new screen will open which is similar to the simple time entry screen. The maximum increase in hold time per cycle (based on the number of cycles and the time limits of the unit) will be shown.

- Enter the required value using the + and - buttons.
- If an entered incremented time is invalid, a prompt will indicate the acceptable range.
- Touch **OK** to return to the previous screen.

The increase in hold time per cycle will now be displayed.

- Touch **OK** again to return to the programming screen.
- Touch **Cancel** to return to the programming screen if no changes are required.

The time value will now appear orange in the programming screen indicating that it contains a modified function.

To decrease the step hold time per cycle:

- For a **decrease** in hold time per cycle, touch the button below **Decrease Per Cycle** to turn this function on. Next touch the time entry button below this.

A new screen will open which is similar to the simple time entry screen. The maximum decrease in hold time per cycle (based on the number of cycles and the time limits of the unit) will be shown.

- Enter the required value.
- If an entered decremented time is invalid, a prompt will indicate the acceptable range.
- Touch **OK** to return to the previous screen.

The decrease in hold time per cycle will now be displayed.

- Touch **OK** again to return to the programming screen.
- Touch **Cancel** to return to the programming screen if no changes are required.


The time value will now appear orange in the programming screen indicating that it contains a modified function.

Note: It is not possible to both increase and decrease the time of the same step.

INSTRUMENT SETTINGS AND DEFAULTS

Touch the **Settings** button to access the instrument Settings menu.

- Touch any of the buttons to view and edit the settings.
- Touch **More Settings** to access further options.
- Once all the settings and defaults have been set as required, touch **Save** to save the changes and return to the Home Screen.
- If you do not wish to change any settings, touch **Cancel** to return to the Home Screen.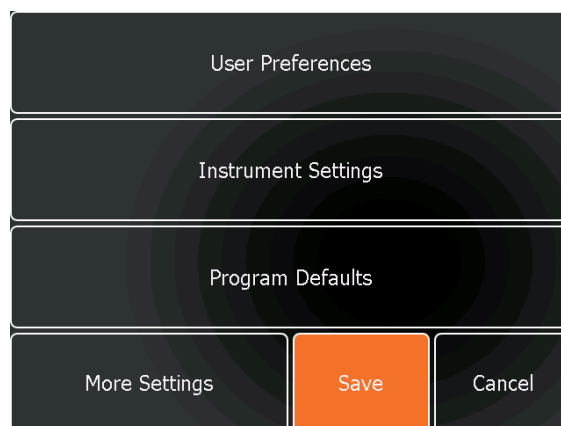

USER PREFERENCES

Language

- Touch the language button to set the required language. The options are English, French, Spanish, Italian and German.


Temperature Display

- Touch the temperature button to set the required temperature units. The options are °C and °F.

Home Screen File List Sort

This sets the user preference for listing the program shortcuts on the Home Screen.


- Touch the button to set the required sorting preference. The options are recently used first or most used first.
- Touch **More Settings** for further user preferences.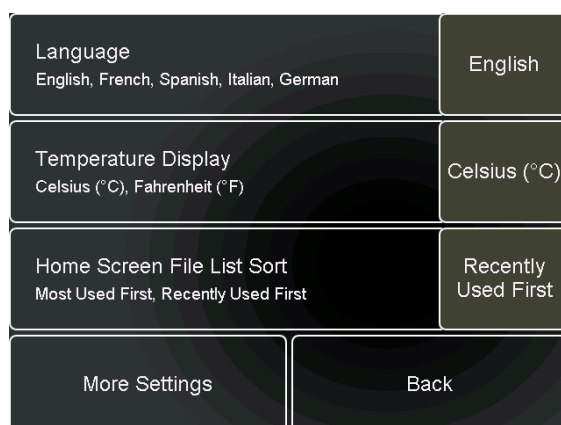

Select Multiple Items in File Manager

This allows either multiple files or only single files to be selected for copying and deleting when using the File Manager in the View files module.

- Touch to select **On** or **Off**.
- Touch **Back** to return to the Settings menu.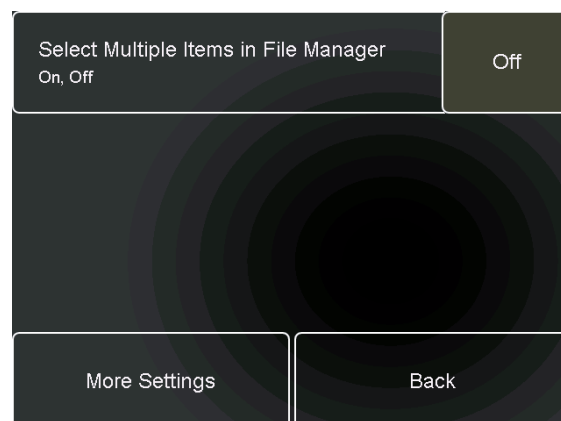

INSTRUMENT SETTINGS

Temperature Log Recording

To record the temperature logs (block temperature profile) of each run, touch the button to set this to Always. If temperature logs are not required, set to Never.

Auto Resume On Power Loss

This determines whether the unit will restart after a power failure. Setting Auto Resume to Always will always allow the unit to re-start irrespective of the length of the power failure.

Date and Time Settings

Use these settings to set the current date, time and date format on the unit.

- When set, touch **Back** to return to the Settings menu.

Temperature Log Recording Always, Never	Always
Auto-Resume On Power Loss Always, Never	Always
Date and Time Settings	
More Settings	Back

Date Current date is 15/12/2011	Set Date
Time Current time is 13:59	Set Time
Date Format dd/mm/yy (15/12/11), mm/dd/yy (12/15/11)	dd/mm/yy
More Settings	Back

Date

- Touch **Set Date** in order to enter the current date.
- Use the + and – keys to set the date values.
- When the correct date has been set, touch **OK** to return to the previous screen.
- Touch **Cancel** to return to the previous screen if no changes are required.

Enter a new date

+		+		+	
15	/	12	/	11	
-		-		-	

OK Cancel

Time

- Touch **Time** in order to enter the current time using the 24 hour clock.
- Use the + and – keys to set the time values.
- When the correct time has been set, touch **OK** to return to the previous screen.
- Touch **Cancel** to return to the previous screen if no changes are required.

Enter a new time

+		+		+	
14	:	27	:	25	
-		-		-	

OK Cancel

Date Format

Touch the date button to set the required date format.

The options are dd/mm/yy or mm/dd/yy.

- Touch **More Settings** for further date and time settings or **Back** to return to the Settings menu.

Administrator Settings

This allows an administrator to remove passwords from password-protected programs, change the Administrator password and restore the unit to its factory settings.

- Touch **None** and enter the administrator password.

The default password is 1234.

- Touch **OK** to enter the Administrator Settings menu.
- Touch **Cancel** to return to the menu without making any changes.

Administrator Password Required.
The administrator password is required to access administrator settings.

Password

The password is correct

OK Cancel

Remove Passwords

- From the Administrator menu touch **Remove Passwords**.
- Touch the file with the password to be removed.
- Touch **Remove Password**.

The password will now be removed and the program can be deleted if required.

- Touch **Close** to return to the menu.

Sort By Program Name (ABC) Search By Keyword Search By User Select All Remove Password

Program 2 Step Template Keywords:None selected User:None selected Used:Today, 10:02 PM	Program 3 Step Template Keywords:None selected User:None selected Used:Never used	Delete View Close
Program Ice Bucket Keywords:None selected User:None selected Used:Never used	Program Ligation 15°C Keywords:None selected User:None selected Used:Never used	
Program RT PCR Template Keywords:None selected User:None selected Used:Never used	Program Test #1 Keywords:None selected User:None selected Used:Never used	
Program Test #2 Keywords:None selected User:None selected Used:Never used		

Change Administrator Password

This allows the administrator to change the default password (1234).

- From the Administrator menu touch **Change Administrator Password**.
- Touch **None** entered and enter the new password.
- Touch the second **None** entered to confirm the new password.

Note: It is recommended that the administrator keeps a record of the new password.

- Touch **OK** to return to the menu or **Cancel** to return without making any changes.

Change Administrator Password
To change the administrator password, the new password must be entered twice below.

Password None entered


Confirm password None entered

OK Cancel

Restore Factory Settings

All changes to the instrument defaults and settings can be removed and the unit returned to its factory settings.

- From the Administrator menu touch **Restore Factory Settings**.
- Touch the **None** button and enter the administrator password.
- Touch **Restore Settings** and the unit will be reset to factory defaults.
- Touch **Cancel** to return to the Settings menu without making any changes.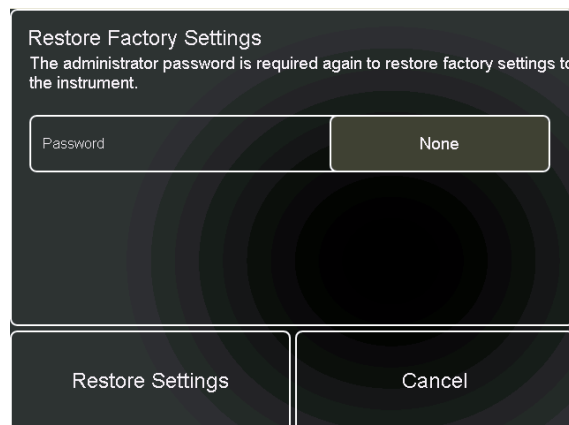


PROGRAM DEFAULTS

The program default settings will apply to every new program created on the Prime. However individual programs can be edited as required.

Heated Lid

Use this to set the heated lid temperature to approximately 10°C higher than the hottest block temperature in the program.

- Touch **On/Off** to toggle between settings.
- Touch the temperature button to set the required lid temperature (35°C to 115°C).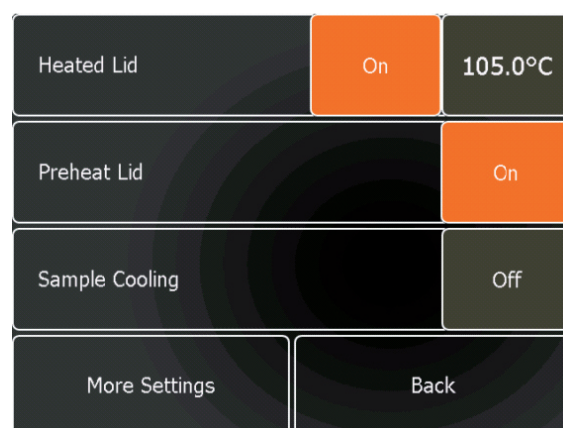

Preheat lid

Set to **On** to pre-heat the lid to the set temperature before the thermal cycling program begins.

- Touch **On/Off** to toggle between settings.

Sample Cooling

Set to **On** to cool the block to 4°C while the lid preheats.

- Touch **On/Off** to toggle between settings.
- Touch **More Settings** for further program defaults or **Back** to return to the Settings menu.

Pause

This will cause the unit to pause before the program starts at the ambient temperature of the block unless Sample Cooling has been set to On. In this case the unit will pause with the block held at 4°C and the program will not commence until the unit is manually un-paused.

- Touch **On/Off** to toggle between settings.

Initial Denaturation

Use this to denature the sample and/or activate the enzyme before the thermal cycling program begins.

- Touch **On/Off** to toggle between settings.
- Touch the time button to set the required hold time.
- Touch the temperature button to set the required temperature.

Hot Start

This will cause the unit to pause at the set temperature. The program will not continue until the unit is manually un-paused.

- Touch **On/Off** to toggle between settings.
- Touch the temperature button to set the required temperature.
- Touch **More Settings** for further program defaults or **Back** to return to the Settings menu.

Pause		Off	
Initial Denaturation	On	00:05:00	94.0°C
Hot Start		Off	95.0°C
More Settings		Back	

Final Extension

Use this to add a final extension step to the program to ensure complete amplification of the products.

- Touch **On/Off** to toggle between settings.
- Touch the time button to set the required hold time.
- Touch the temperature button to set the required temperature.

Final Hold

This will hold the block at the set temperature at the end of a program until the user manually stops the unit.

- Touch **On/Off** to toggle between settings.
- Touch the temperature button to set the required temperature.
- Touch **More Settings** to return to the top of the Program Defaults menu or **Back** to return to the Settings menu.

Final Extension	On	00:05:00	72.0°C
Final Hold	On		10.0°C
More Settings		Back	

SERVICE SETTINGS

The service settings are for the use of authorised service engineers only. A service engineer's password is required to access these settings.

GRADIENT CALCULATOR (PRIMEG ONLY)

The Gradient Calculator can be used to calculate the temperature for each column of the block for a user-specified temperature and gradient range.

- Touch the **Target Temperature** button and enter the required temperature.
- Touch **OK**.
- Touch the **Temperature Gradient** button and enter the required gradient range followed by **OK**.
- Touch **OK**.

The calculated temperatures for each of the columns of the block will now be displayed.

Note: With the 384-well block, the columns are paired as it is not possible to individually control 24 columns. Therefore only 12 column temperatures are available across this block.

Target Temperature Enter the temperature used for the gradient	55°C
Temperature Gradient Enter the temperature gradient used	10°C
50.7 51.2 52.3 53.3 54.3 54.9 55.5 56.2 56.9 58.1 59.2 59.7	
Back	

OLIGO T_m CALCULATOR

The oligo T_m calculator can be used to estimate the T_m of the PCR primers and help in selecting an appropriate annealing temperature for the thermal cycling program.

- Touch **Enter The Oligo Sequence** in order to input the sequence of the primer.

Enter The Oligo Sequence	
Primer Concentration Enter the concentration of the primer in μM	0 μM
Salt Concentration Enter the salt concentration in mM	0mM
Calculate	Reset Back

- Touch the **A C G T** buttons to enter the primer sequence.
- Touch **DEL** to delete the last character.
- Touch **CLR** to delete the whole sequence.
- Use the arrow keys to scroll through the sequence.
- Touch **OK** to return to the calculator screen.

DNA Base Entry Bases					
5'	ATTGTAATTCGTAC				3'
CLR				DEL	
A	C	G	T	<	>
OK			Cancel		

- Enter the primer concentration in the PCR mix.
- Enter the monovalent salt concentration of the PCR mix.
- Touch **Calculate** to calculate the T_m of the primer.

ACGGCAATTGTAATTCGTAC	
Primer Concentration Enter the concentration of the primer in μM	0.5 μM
Salt Concentration Enter the salt concentration in mM	50mM
Calculate	Reset Back

- To check the T_m of another primer with the same primer and salt concentrations, touch **Back** and enter the new primer sequence.
- Touch **Reset** to clear the entries.
- Touch **Back** to return to the Settings menu.


T _m Calculation 50mM (salt) and 0.5 μM (primer) ACGGCAATTGTAATTCGTAC	
58.94°C	
Calculate	Reset Back

- Once all the settings and defaults have been set as required, touch **Save** to save the changes and return to the Home Screen.
- If you do not wish to change any settings, touch **Cancel** to return to the Home Screen.

CREATE A NEW PROGRAM

From the Home Screen, touch the **Create A New Program** button. This will open a screen where you can build a program by adding stages and steps and define the temperatures and hold times.

The function of each of the buttons on this screen is described below.

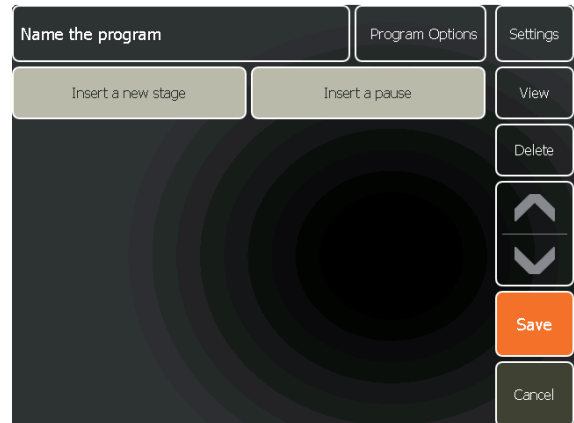
NAME THE PROGRAM

- Touch this button to give the program a name.
- Type in the required name using the text entry screen.
- Once the name is complete, touch **OK** to return to the programming screen.
- Touch **Cancel** to return to the programming screen if no changes are required.

When saving the program, if the chosen name has already been used for another program, the message box shown opposite will appear.

- Touch **OK** to return to the Programming screen.
- Touch the program name and enter a different name.
- Once the name is complete, touch **OK** to return to the programming screen.
- Touch **Save** to save the program.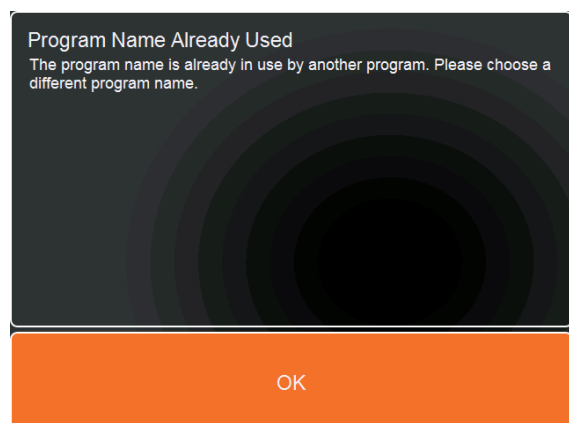

INSERT/DELETE A STAGE

A **stage** is a section of a program that consists of a number of individual steps that can be cycled (repeated) a defined number of times.

Insert a new stage

- To insert a stage in a program, touch **Insert a new stage**. A stage with one step will be inserted.

Stage 1

- Touch this to give the stage a name (optional). Enter the name using the text entry screen.

1 cycle

- Touch this to edit the number of cycles (repeats) for the stage. Enter the required number of cycles then touch **OK**.

Temperature

- Touch the temperature button to edit the temperature of the step and add any increment/decrement or gradient features.

Time

- Touch the time button to edit the hold time of the step and to add any increment/decrement time features.

Ramp rate

- Touch the ramp rate setting to adjust the ramp rate.

The range is 0.1°C/s (minimum) to 3.4°C/s maximum for heating and 1.0°C/s for cooling. The value can be entered to 1 decimal place. Ramp rate will affect the heating or cooling rate from the previous step to the current step.

- Once set, touch **OK** to return to the Programming Screen.

Delete a stage

- To **delete** a stage, touch the stage area to highlight it and then touch **Delete**.

INSERT/DELETE A STEP

Insert a new step

- Touch **Insert a new step** to insert the next step in the current stage.
- Edit the new step temperature, hold time and ramp rate as required.

Delete a step

- To delete a step, touch the step area to highlight it and then touch **Delete**.

Test #1			Program Options	Settings
Amplification			35 cycles	View
Temperature	Time	Ramp Rate	Step 1	Delete
95.0°C	00:00:30	Maximum		↑
Temperature	Time	Ramp Rate	Step 2	↓
60.0°C	00:00:40	Maximum		Save
Insert a new step				Cancel

INSERT/DELETE A PAUSE

Insert a pause

- To insert a pause between stages, touch **Insert a pause**.

The hold temperature will be that of the previous step. The unit will not continue with the remainder of the program until it is manually un-paused by the user.

Delete a pause

- To delete a pause, touch the pause button to highlight it and then touch **Delete**.

			Settings
Temperature	Time	Ramp Rate	Step 2
60.0°C	00:00:40	Maximum	View
Temperature	Time	Ramp Rate	Step 3
72.0°C	00:01:00	Maximum	Delete
Insert a new step			↑
Pause			↓
Insert a new stage			Save
Insert a pause			Cancel

Note: To insert a pause between two previously programmed stages, it is necessary to first delete the later stages to where the pause is required.

PROGRAM OPTIONS

- Touch this to view the program defaults as set up in the Settings module.

Any of these parameters can be edited by touching the appropriate button and will be saved for the current program only.


- Touch **Program** to return to the programming screen.

Pause	Off	Program
Initial Denaturation	On	94.0°C
Hot Start	Off	95.0°C
Final Extension	On	72.0°C
Final Hold	On	10.0°C
User	None selected	Save
Keywords	None selected	Cancel
Password	None	


User

User names can be entered and linked to the program to help in searching and identifying programs.

- To enter a user name, touch the **None selected** button next to User.
- User names that have already been created will then be listed.
- To create a new user name, touch the **Create a new user** button.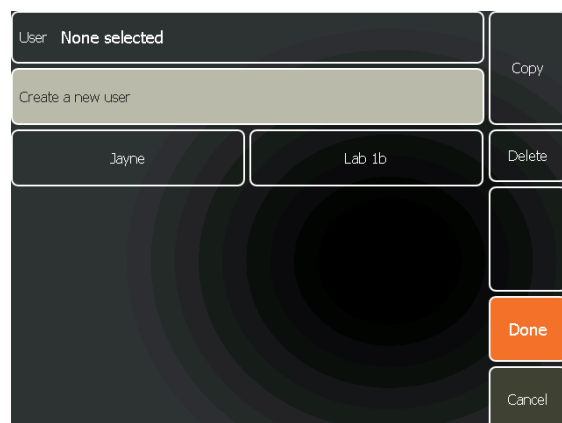

- Type in the user name and touch **OK**.

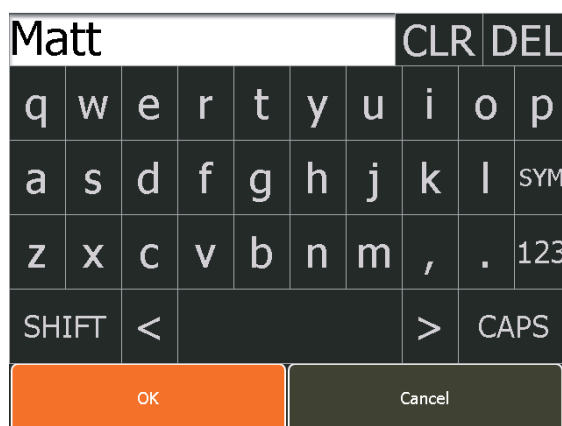
The new user name will be added to the list.

- To select a user name, touch the button to highlight it. Touch it again to de-select it.
- To delete a user name from the list, select the user name and touch **Delete**.
- Touch **Done** to return to Program Options.

The user name will now appear in the Program Options screen.


- Touch the **Program** button to return the programming screen.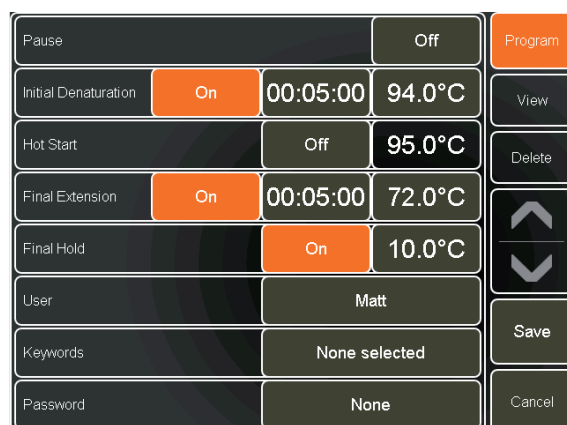

Keywords

Keywords can be entered and linked to the program to help in searching for related programs.

- To enter a keyword, touch the **None selected** button next to Keywords.
- Other keywords that have already been created will then be listed.
- To create a new keyword, touch the **Create a new keyword** button.

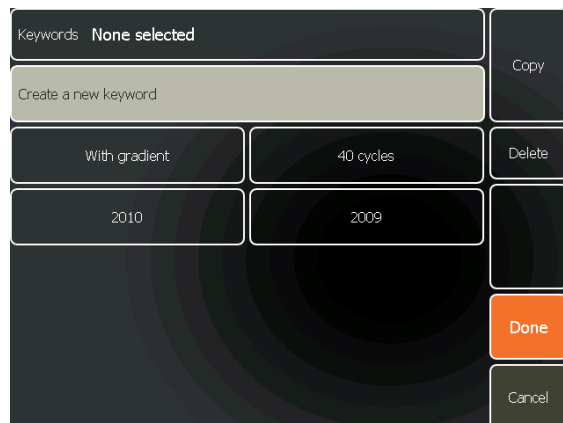
- Type in the keyword and touch **OK**.


The new key word will be added to the list.

- To select a keyword, touch the button to highlight it. Touch it again to de-select it. More than one keyword can be selected
- To delete a keyword from the list, select the keyword(s) and touch **Delete**.
- Touch **Done** to return to Program Options.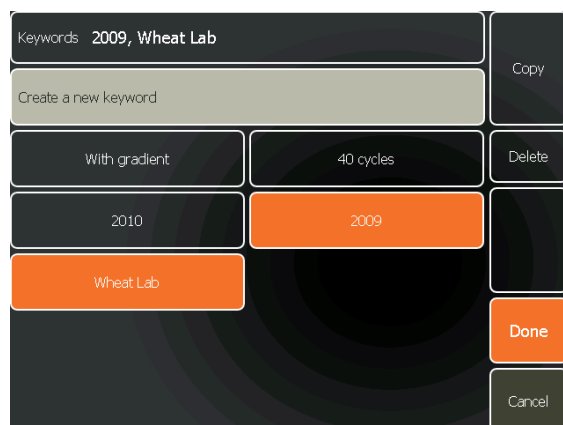

The keyword(s) will now appear on the Program Options screen.


- Touch the **Program** button to return the programming screen.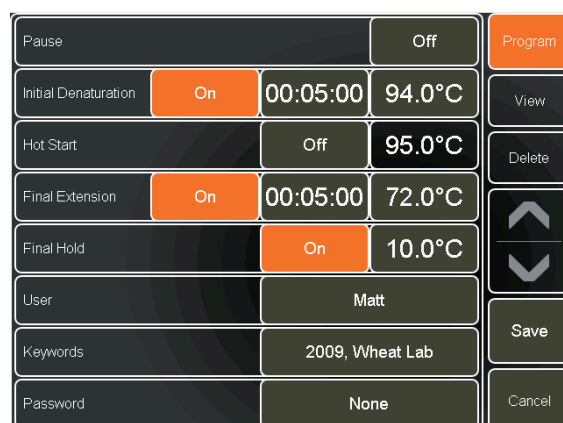


Passwords

A password can be created to protect the program from being edited and deleted by other users. The password will be requested if a user attempts to edit or delete the program.

- To create a password, touch the **None** button next to Password.
- Enter the required password then touch **OK**. The password is case-sensitive.
- Touch **Cancel** to return to Program Options without entering a password.
- Touch the **Program** button to return to the Programming screen.

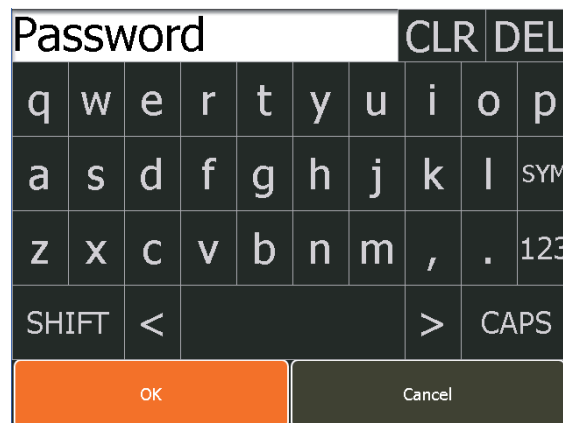
The Password creation screen features a title bar with 'Password', 'CLR', and 'DEL' buttons. Below is a QWERTY keyboard layout with a 'SYM' button for symbols. At the bottom, there are 'SHIFT', '<', '>', and 'CAPS' buttons, followed by an orange 'OK' button and a grey 'Cancel' button.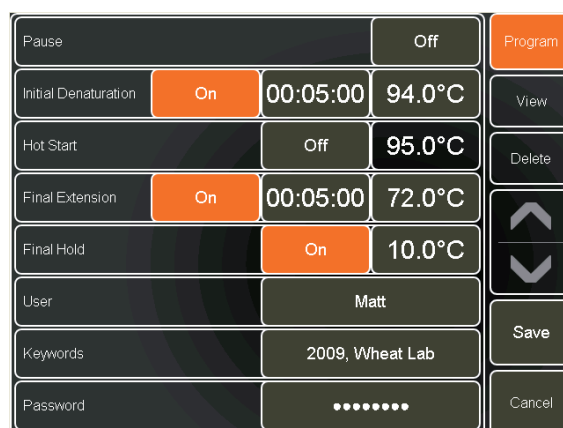


The Program Options screen displays various settings in a table-like format. On the right side, there are buttons for 'Program', 'View', 'Delete', navigation arrows, 'Save', and 'Cancel'. The settings include: Pause (Off), Initial Denaturation (On, 00:05:00, 94.0°C), Hot Start (Off, 95.0°C), Final Extension (On, 00:05:00, 72.0°C), Final Hold (On, 10.0°C), User (Matt), Keywords (2009, Wheat Lab), and Password (represented by dots).

- To change or remove the password, from the Program Options screen, touch the button next to **Password**.
- On the Password Editing screen, touch the **None** button and enter the password.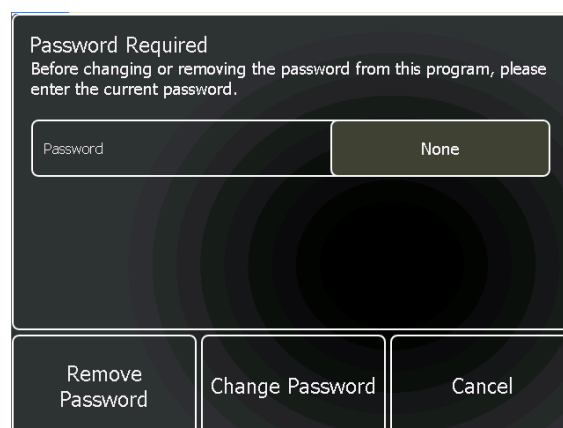

The Password Required screen has a title 'Password Required' and a message: 'Before changing or removing the password from this program, please enter the current password.' It shows a 'Password' field with a 'None' button next to it. At the bottom are three buttons: 'Remove Password', 'Change Password', and 'Cancel'.

- To remove the password, touch **Remove Password**.
- To change the password, touch **Change Password** and type in the new password, followed by **OK**.
- Touch **Cancel** to return to the Program Options screen if no change is required.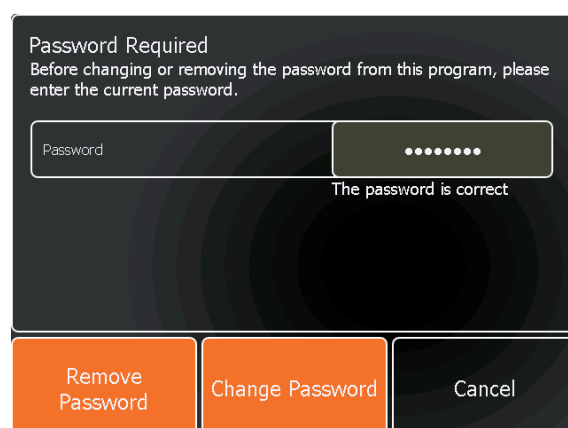


This screen is identical to the previous one, but the 'Password' field now contains dots and the text 'The password is correct' is displayed below it. The buttons at the bottom remain 'Remove Password', 'Change Password', and 'Cancel'.

SETTINGS

- Touch the **Settings** button to view the instrument defaults as set up in the instrument Settings module.

Any of these parameters can be edited by touching the appropriate button and will be saved for the current program only.

- Touch **Program** to return to the programming screen.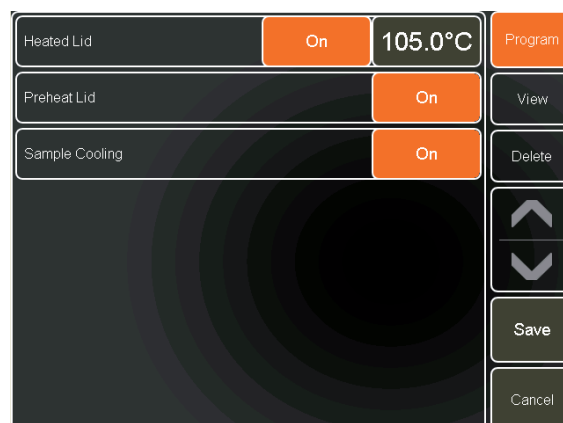


VIEW

- Touch **View** to see a graphical representation of the program.
- If the program has more stages and steps than are displayed on the screen, scroll through the program by dragging a finger across the touch screen.

The centre of the screen shows each temperature step and hold time (not to scale). Stages are indicated by a white bar spanning the steps within the stage.

The bottom of the screen shows the name of the program and the expected end time if it was started immediately.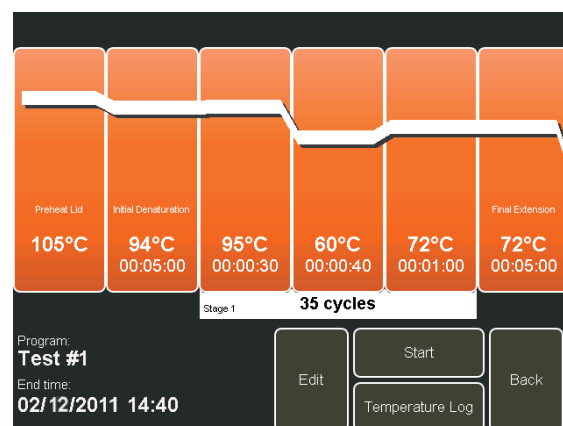

Note: The program is not saved at this point.


- Touch **Start** to start the program.
- Touch **Edit** to edit the program.
- Touch **Back** to return to the Programming Screen.

Note: The temperature log will remain blank until the program has been started.

If a step has a gradient programmed into it, the bar will be shown thicker and indicate the upper and lower temperatures of the gradient range.

If a step has an incremented or decremented temperature, this will be indicated by a white spot next to the temperature.

If a step has an incremented or decremented time, this will be indicated by a white spot next to the time.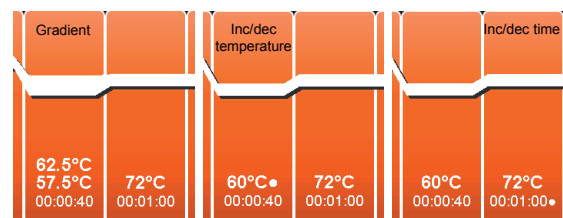

DELETE

This is used to delete an entry. First touch the required segment followed by **Delete**.

UP AND DOWN ARROWS

Use these buttons to scroll through a program.

SAVE

Touch this button to save the program.

CANCEL

Touch this button to cancel changes to the program and return to the Home Screen.


RUN A PROGRAM

From the Home Screen, touch one of the shortcuts to run one of the listed programs, or touch **Run a program** to search and select from other programs and templates saved on the unit. Selected programs can be run directly or edited before running.

RUNNING A PROGRAM FROM THE SHORTCUT LIST

On the Home Screen, the Run a program area displays shortcuts to four programs. User Preferences in the Settings module can be used to select whether this displays the four most **recently** used programs or the four most **frequently** used programs.

- To run one of the programs in the shortcut list, touch the button to open the Pre-run Screen.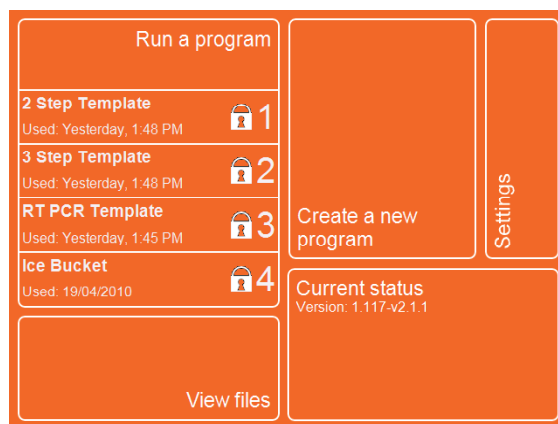

THE PRE-RUN SCREEN


The Pre-run Screen shows a graphical representation of the program allowing for a rapid check of the temperature parameters.

The centre of the screen shows each temperature step and hold time (not to scale). Stages are indicated by a white bar spanning the steps within the stage.

The bottom of the screen shows the program name, the date and the expected end time if it was started immediately.

- Touch **Start** to start the program.
- Touch **Edit** to edit the program.
- Touch **Back** to return to the previous screen.

Note: The temperature log will remain blank until the program has been started.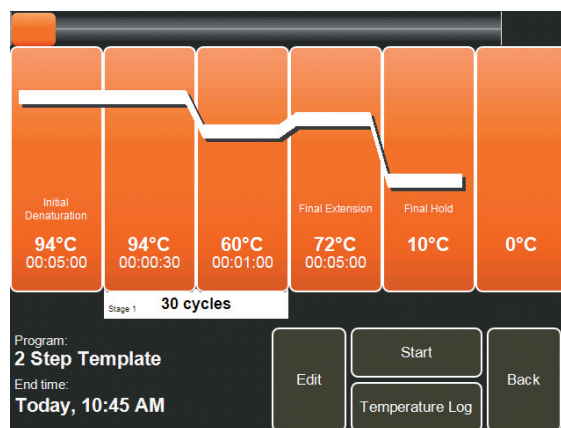

THE RUN SCREEN

The Run Screen will be displayed after a program has been started. It shows the instrument status and the progress of a run.

Once the program has been started, the timers in the top right hand corner of the screen will indicate how long the program has been running and the total run time. The progress bar at the top of the screen also indicates how far the run has progressed.

Adaptive ramp rate

The first time a program is run, the unit estimates the total run time based on the ramp rates set in the program. After the first cycle has completed the actual time taken to complete the cycle is then used to re-calculate the total run time. Subsequent runs of the same program will therefore have a more accurate estimate of total run time. The estimate also takes into account the ambient room temperature; therefore run times will also depend on ambient temperature.

The status of the block is indicated by the step bars which display the activity of the block (i.e. ramping to temperature), paused or the count down time of the step.

Progress through the program is indicated by the stage bar(s) which show at which cycle of the program the unit is currently at.

- Touch **Stop** to stop the program. A prompt will ask you to confirm the stop.
- Touch **Temperature log** to view the block temperatures. The previous two minutes of temperature readings can be viewed by scrolling across the screen using a finger.



Note: The complete log is written and saved on the unit only if Temperature Log Recording is selected as Always in the Instrument Settings.

Note: Touch **Program View** to return to the program status.

- Touch **Pause** to pause the program.
- If paused, touch **Continue** to un-pause the program.
- Touch **Back** to return to the Home Screen.

Note: It is possible to access all modules of the software while the unit is running a program.


Once the run has finished, Program Finished will be displayed instead of End time.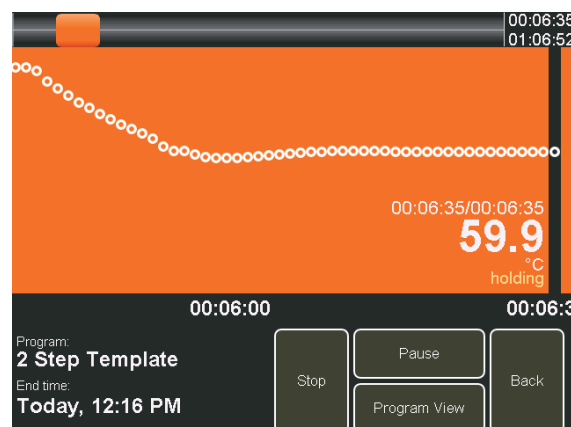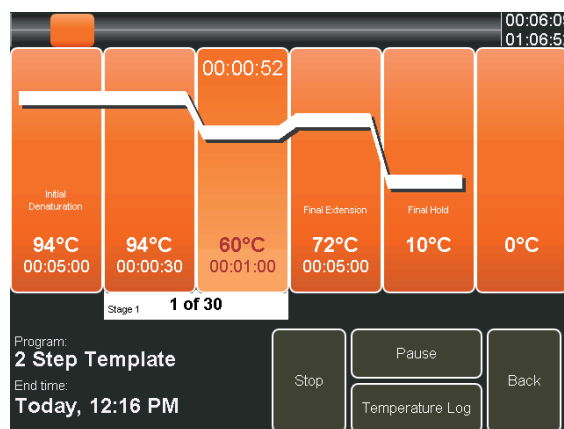

Viewing a running program

From the Home Screen:

- Touch **Current status**.
- Touch the area for the instrument.
- Touch **View**.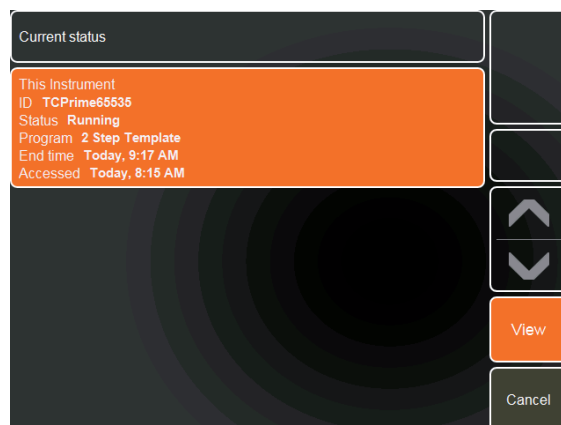


SEARCHING FOR A PROGRAM NOT IN THE SHORTCUT LIST

There are four ways to search for a program if it is not in the shortcut list:

- Search from the complete list of saved programs.
- Search by keyword.
- Search by user.
- Search by both keyword and user.

Search from a list

- To search for a program saved on the unit, touch **Run a Program** to open the Program Selection Screen.
- Use the + and – buttons to scroll up and down the list of stored programs.
- Touch **Sort By Program Name** to list the programs in order. The options are:
 - Program Name (ABC)/(ZYX)
 - Recently Used (123)/(321)
 - Most Used (123)/(321)
 - User (ABC)/(ZYX)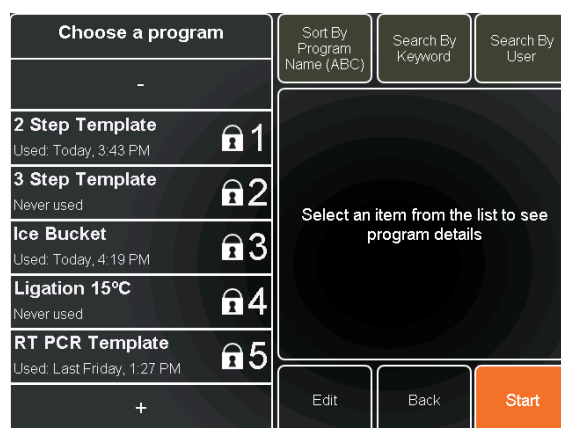


Search the programs by keyword

- Touch **Search by Keyword** and this will open the list of keywords stored on the unit.
- Select one or more keywords by touching the buttons to highlight them. Touch again to de-select.
 - When all keywords have been selected, touch **Done** to return to the Program selection screen.

Only the programs with the selected keyword(s) saved with them will be listed.

- To clear the keyword search, touch **Clear Keyword Search**.

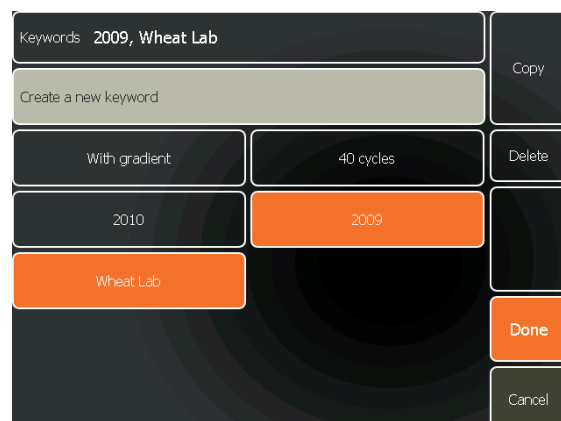Search the programs by user

- Touch **Search by User** and this will open the list of user names stored on the unit.
- Select the required user name by touching the appropriate button to highlight it. Touch again to de-select.
- When the user name has been selected, touch **Done** to return to the Program selection screen.

Only the programs with the selected user name saved with them will be listed.

- To clear the user name search, touch **Clear User Search**.


Search the programs by keyword and user

- Touch **Search by Keyword** and this will open the list of keywords stored on the unit.
- Select one or more keywords by touching the buttons to highlight them. Touch again to de-select.
- When all keywords have been selected, touch **Done** to return to the Program selection screen.

Only the programs with the selected keyword(s) saved with them will be listed.

- Touch **Search by User** and this will open the list of user names saved with previously keyword-searched programs.
- Select the required user name by touching the appropriate button to highlight it. Touch again to de-select.
- When the user name has been selected, touch **Done** to return to the Program selection screen.

Only the programs with both the selected user name and key word(s) saved with them will be listed.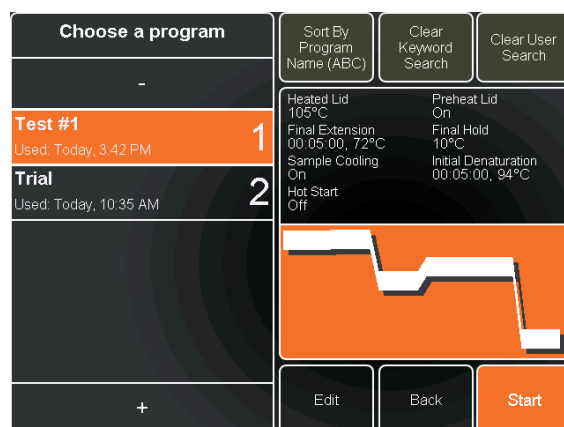


RUNNING A PROGRAM NOT ON THE SHORTCUT LIST

- Once the required program has been identified, touch it to highlight it.

The Program Selector Screen will show a summary of the program, the instrument settings and a graphical representation of the thermal cycling program.

- Touch **Edit** to edit the program if required.
- Touch **Back** to return to the Home Screen.
- Touch **Start** to proceed to the Pre-run Screen.
- From the Pre-run Screen, touch **Start** to run the program.

See **THE RUN SCREEN** on page 46 for further information on instrument status and progress during a run.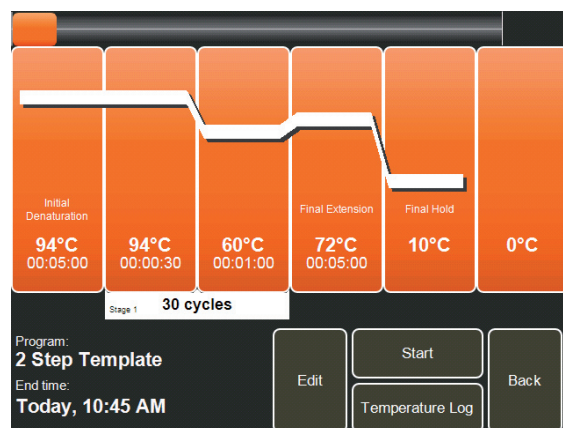

EDITING A PROGRAM

EDITING SAVED PROGRAMS

First select the required program by searching the **Run a program** module.

- To edit a selected program, touch the **Edit** button.

Any of the Program Options, Settings or thermal cycling parameters can be edited.

- Once all the required changes have been made, touch **Save**.

Note: this will overwrite the existing program.

To copy a program for editing, please see section **VIEW FILES**.

Test #1			Program Options	Settings
Amplification			35 cycles	View
Temperature	Time	Ramp Rate	Step 1	Delete
95.0°C	00:00:30	Maximum		
Temperature	Time	Ramp Rate	Step 2	
60.0°C	00:00:40	Maximum		
Temperature	Time	Ramp Rate	Step 3	
72.0°C	00:01:00	Maximum		
				Save
				Cancel

EDITING TEMPLATES

Editing templates can be a quick method of creating a program by just changing a few key parameters. However it is not possible to save changes to the templates installed on the thermal cycler.

If the edited program is a template, touching **Save** will open the message box shown opposite.

- Touch the **None** button next to New Program Name and enter a name for the edited program.
- Touch **Save As New Program** to save the changes.

The re-named program will now be listed in the program list.

This Program Is A Template
A template is a program which comes installed on the thermal cycler, and cannot be changed. You can save the changes made to the template...

New Program Name None

Save As New Program Cancel

EDITING PASSWORD PROTECTED PROGRAMS

If the edited program is password protected, touching **Save** will open the message box shown opposite. To save the changes either:

- Touch the **None** button next to Password and enter the correct password followed by **Save**.

Note: this will overwrite the existing program.

or

- Touch the **None** button next to New Program Name and enter a new name for the edited program, followed by **Save as a New Program**.

The re-named program will now be listed as a separate program in the program list.

This Program Is Locked
You can save any changes made by entering the password, or save the new program details as a new program with a different name.

Password None

New Program Name None


Save Save As New Program Cancel

CURRENT STATUS

The **Current Status** area of the Home Screen provides status information on the Prime and is used to access the Run Screen.

On the Home Screen, the Current status area shows the versions of software and firmware loaded on the unit.


- Touch the **Current status** button to view the status of the Prime.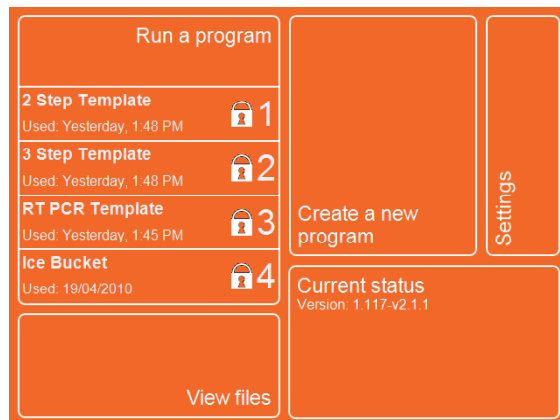

The Prime status will be shown.

If a unit is running a program, as shown by the status, the Run Screen can be viewed. Runs can then be paused or stopped as required.

- Touch the area to select the unit.
- Touch **View**.
- Touch **Cancel** to return to the Home Screen.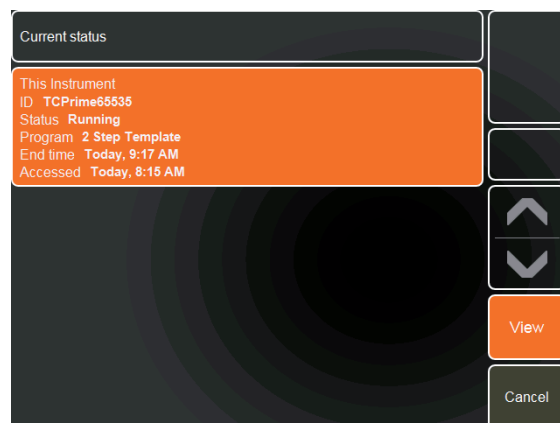


VIEW FILES

The **View files** module opens the **File Manager** and allows access to all the program and temperature log files. From here it is possible to copy program files to the unit for editing. Files can also be deleted or copied to a USB memory stick for transfer to a PC or storage.

FILE SEARCHING


There are four ways to search for a file:

- Search program files or temperature logs from the complete list of saved files.
- Search by keyword.
- Search by user.
- Search by keyword and user.
- To search for a file saved on the unit, touch **View Files** to open the File Manager screen.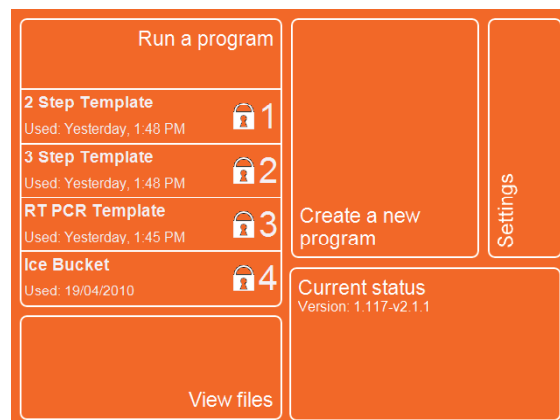

Search from a list

- Touch **Programs And Logs** to display either programs only, temperature logs only or both.
- Use the up and down arrow buttons to scroll through the list of saved files.
- Touch **Sort By Program Name** to list the programs in order. The options are:
 - Program Name (ABC)/(ZYX)
 - Recently Used (ABC)/(ZYX)
 - Most Used (ABC)/(ZYX)
 - User (ABC)/(ZYX)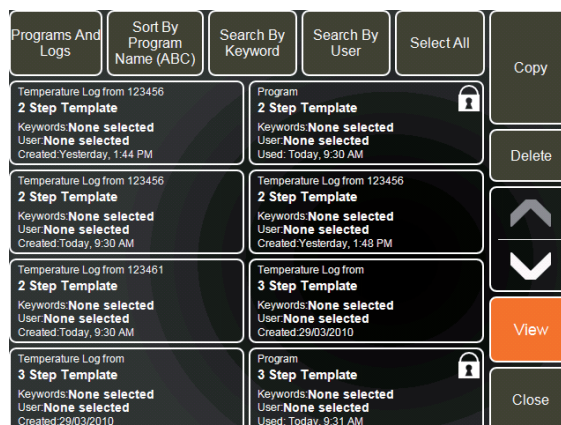


Search the files by keyword

- Touch **Search By Keyword** and this will open the list of keywords stored on the unit.
- Select one or more keywords by touching the buttons to highlight them. Touch again to de-select.
- When all keywords have been selected, touch **Done** to return to the File Manager screen.

Only the files with the selected keyword(s) saved with them will be listed.

- To clear the keyword search, touch **Clear Keyword Search**.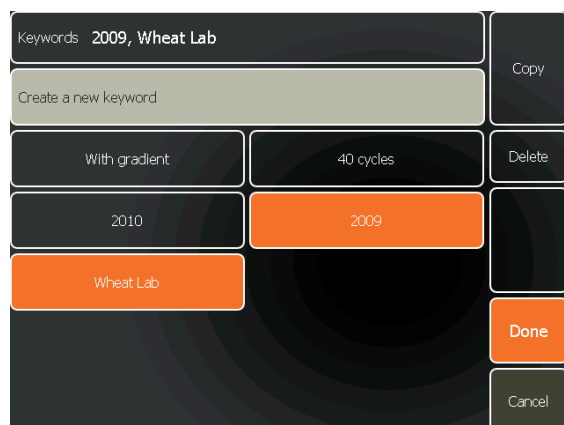


Search the files by user

- Touch **Search By User** and this will open the list of user names stored on the unit.
- Select the required user name by touching the appropriate button to highlight it. Touch again to de-select.
- When the user name has been selected, touch **Done** to return to the File Manager screen.

Only the programs with the selected user name saved with them will be listed.

- To clear the user name search, touch **Clear User Search**.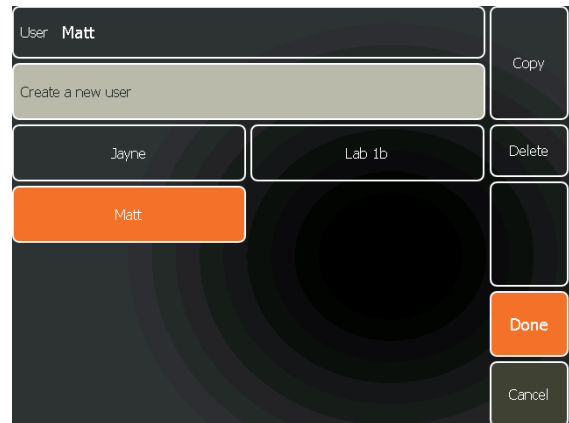


Search the files by keyword and user

- Touch **Search By Keyword** and this will open the list of keywords stored on the unit.
- Select one or more keywords by touching the buttons to highlight them. Touch again to de-select.
- When all keywords have been selected, touch **Done** to return to the Program selection screen.

Only the programs with the selected keyword(s) saved with them will be listed.


- Touch **Search By User** and this will open the list of user names saved with previously keyword-searched files.
- Select the required user name by touching the appropriate button to highlight it. Touch again to de-select.
- When the user name has been selected, touch **Done** to return to the File Manager screen.

Only the files with both the selected user name and keywords saved with them will be listed.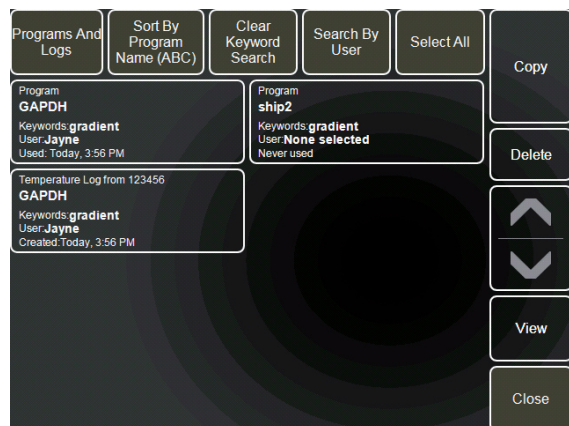


VIEWING A PROGRAM FILE

- Touch **Programs And Logs** to display Only Programs.
- Search for the required program file and touch it to highlight it.
- Touch **View** to open the temperature profile (Pre-run Screen) of the program.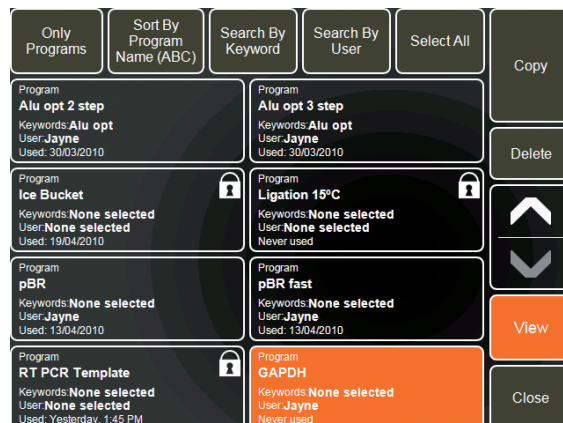

- Touch **Start** to start the program.
- Touch **Edit** to edit the program.
- Touch **Back** to return to the previous screen.

Note: The temperature log will remain blank until the program has been started.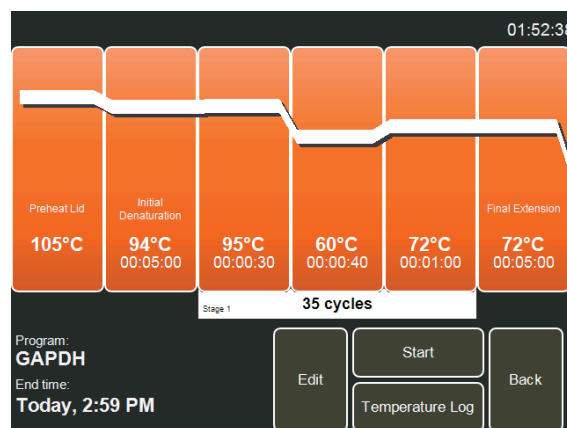


VIEWING A TEMPERATURE LOG FILE

It is not possible to view a temperature log file on the Prime. It must first be copied to a USB memory stick then viewed on a PC using the Techné Workbench software, which can be downloaded free of charge from www.techné.com

- See the directions given in the next section for copying files to a USB memory stick.
- Copy the files to a PC and open the Workbench software.
- Use **Settings** to select the path for the saved files or click on **Open** to browse for the saved files.
- Select the required temperature log and click **Open**.

Note: Since the .temperaturelogdata file is a text file, it can also be opened in applications such as Excel for further analysis.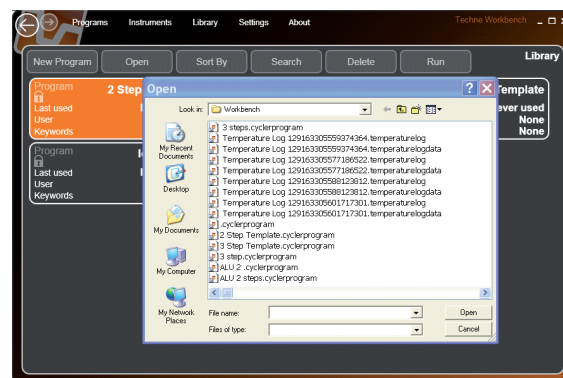


When opening a temperature log file, the Temperature Log Monitor screen will be displayed. To view the complete temperature log, drag the mouse across the screen to scroll through the data points.

The view will show a graphical representation of the actual temperature changes, second by second, throughout thermal cycling program.

Clicking on the **Program View** button of this screen will display the program that generated the temperature log.

For further information on using the Techne Workbench software, see the separate user manual which can be downloaded from www.techne.com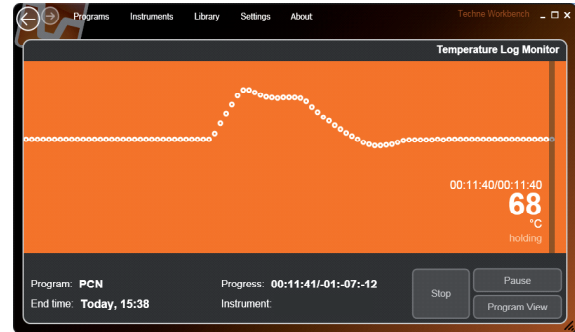

COPYING FILES

Files can either be copied onto the unit for editing, or to and from a USB memory stick.

To copy files to the unit:

- Select one or more files for copying by touching the appropriate file button.



Note: Multiple file selection is optional and can be set up in the User Preferences section of the Settings Module.

or

- Touch **Select All** to select all files.
- Touch **Copy**.

A message box will confirm that the files have been copied.

The copied file(s) will now be highlighted and listed as Copy of... and can be edited as required.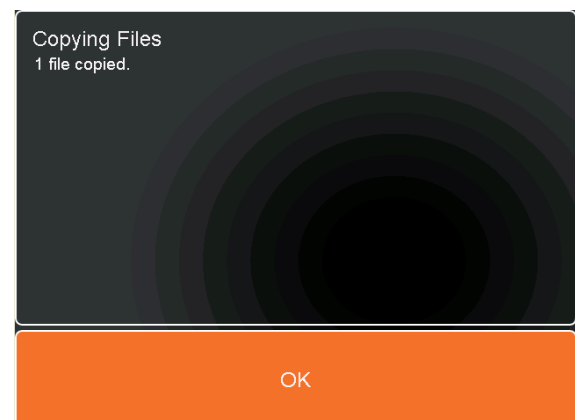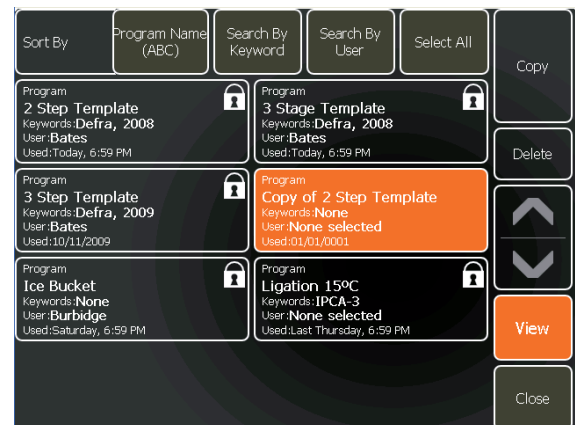

To copy files to a USB memory stick:

- Insert a USB memory stick into the port on the front of the unit.

A message screen will appear confirming that the USB memory stick has been inserted.

- Touch **Cancel** to return to the File Manager Screen.
- Select one or more files for copying by touching the appropriate file button.



or

- Touch **Select All** to select all files.
- Touch **Copy**.

A message screen will appear giving the option of saving the files to the unit or to the USB memory stick.

- Touch the upper button to save the file(s) to the USB memory stick.
- Touch **Cancel** to return to the File Manager screen.

The file(s) will be saved on the USB memory stick in a folder called FileSystem as a .cyclerprogram file.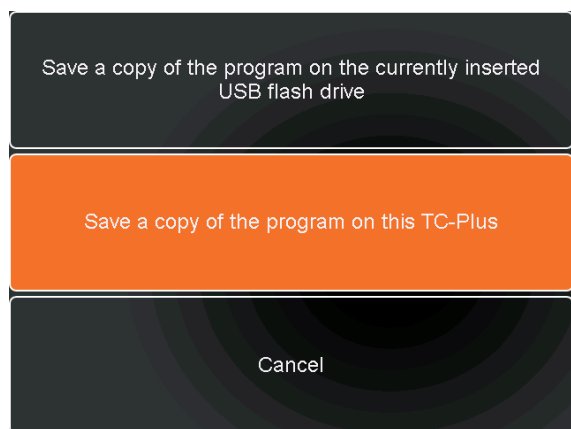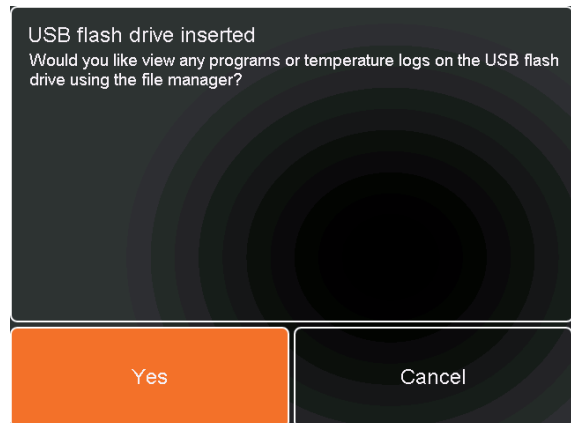

To copy files from a USB memory stick:

- Insert a USB memory stick into the port on the front of the unit.

A message screen will appear confirming that the USB memory stick has been inserted.

- Touch **Yes** to view the programs on the USB memory stick in the File Manager.

A USB button will be visible in the top right corner indicating that the files are being read from the USB memory stick.


- Select one or more files for copying by touching the appropriate file button.

or

- Touch **Select All** to select all files.
- Touch **Copy**.

A message screen will appear giving the option of saving the files to the unit or to the USB memory stick.

- Touch the upper button to save the file(s) to the Prime unit.
- Touch **Cancel** to return to the File Manager screen.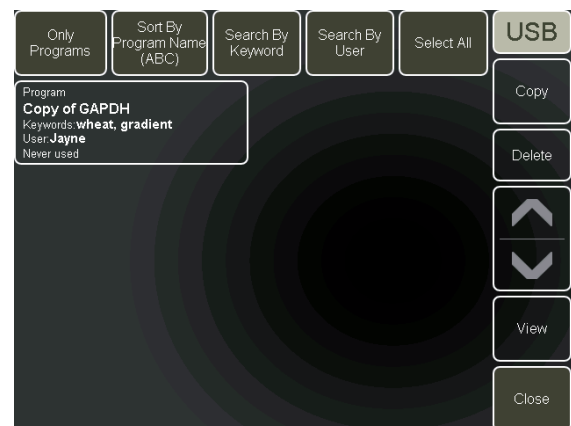

DELETING FILES

Unused or old files can be permanently deleted from the unit or a USB memory stick.

- Select one or more files for deleting by touching the appropriate file button.



Note: Multiple file selection is optional and can be set up in the User Preferences section of the Settings Module.

or

- Touch **Select All** to select all files.
- Touch **Delete**. A message box will appear asking for confirmation of delete.
- Touch **Yes** to confirm the delete.
- Touch **Cancel** to abort the delete.

If the delete is confirmed, a message box will confirm that the files have been deleted.

- Touch **OK**.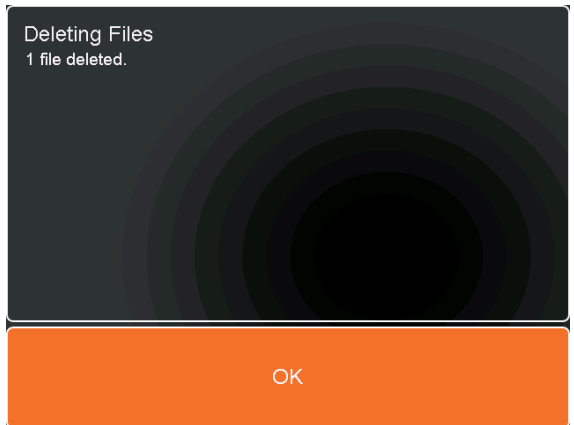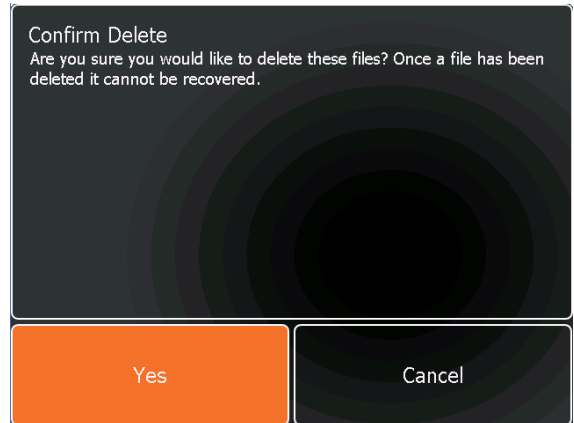

UPDATING THE PRIME SOFTWARE

The Prime software can be updated by downloading the latest version from the Techne website www.techne.com. You will need a USB memory stick on which to save the update.

Note that the update tool requires at least Windows XP SP2 and the Microsoft .Net Framework 3.5 SP1, which, if needed, can be downloaded from:

<http://www.microsoft.com/downloads/details.aspx?familyid=AB99342F-5D1A-413D-8319-81DA479AB0D7&displaylang=en>

- From the website, click on **Support** followed by **Software Downloads**.
- Complete the short registration form and click on **Submit**.

You will now have access to all software and manual downloads.


- Select the **Prime Update** and choose the option to run the file.
- Insert a USB memory stick into the PC and click to select.
- Click on **Prepare the USB memory stick**.

This will copy the update to the USB memory stick.

- Turn the Prime off and insert the USB stick into the front USB port.
- Turn the Prime on and follow the instructions on the screen.

The update will take several minutes. When finished, the home screen will be displayed.


- Remove the USB memory stick from the Prime.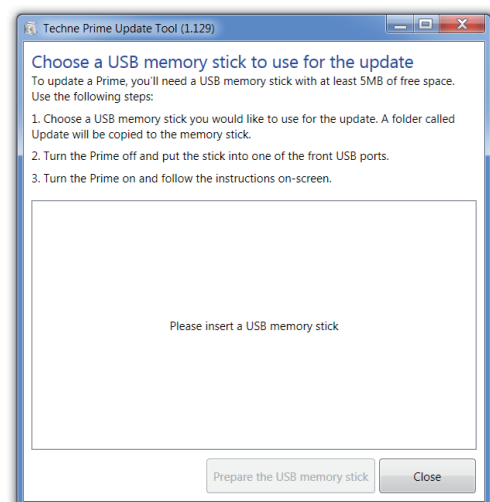

REMOVING AND EXCHANGING BLOCKS

The thermal cycling block can be removed and exchanged for one of a different format or for detailed cleaning.

1. Before removing the block, disconnect from the power supply and allow it to cool to below 50°C.
2. Slide the quick release handles in the direction of the arrows indicated and lift the block assembly out of the unit.
3. Never lift or carry the block by one quick release handle, always use both handles or support the block from underneath.
4. When re-fitting the block, or fitting a new one, slide the quick release handles in the opposite direction to lock the assembly into the unit.
5. Open or closed padlock symbols are printed on the quick release handles for easy reference.

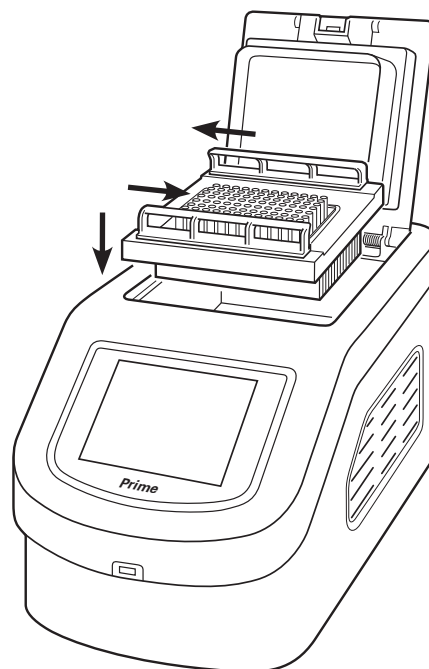UPGRADING FROM PRIME TO PRIMEG

The standard Prime thermal cycler can be easily and quickly upgraded to the gradient PrimeG version by purchasing the PRIMEX/USB upgrade. The upgrade consists of a custom USB stick which will **only** work in a specified unit based on the unit serial number.

1. Turn on the Prime then insert the upgrade USB stick into the port on the front of the unit.
2. A screen will appear indicating the upgrade to the unit. The upgrade will be performed automatically.
3. Remove the USB stick once the upgrade is complete.

FAQS

Q1 *How do I adjust the pressure of the heated lid for my tubes or plates?*

A1 For thermal cyclers with lid adjustment knobs, rotate the knob anticlockwise until there is no pressure on the consumable, then close the lid and latch it. To obtain the correct pressure gently rotate the knob clockwise until it is possible to just feel the pressure being applied. Finally, rotate the knob a further quarter of a turn; the lid is now at the correct pressure for use with the consumable. Ensure that all consumables used in the block are of the same height and are spread evenly across the block. Insert empty “dummy” tubes if necessary to spread the pressure of the heated lid evenly.

Q2 *What is the “Pause” function at the start of the program used for?*

A2 Some users prefer to preheat the heated lid before placing the samples into the unit. The pause feature is used to pause the unit after the 2 minute heated lid preheat step.

Q3 *What is “Sample Cooling”?*

A3 Sample cooling can be applied during the lid pre-heat phase. This maintains the block at 4°C while the lid is heating to temperature and avoids any increase in sample temperature due to the heated lid before the thermal cycling program begins.

Q4 *What is “Hot start”?*

A4 The Hot Start step is used to pause the machine at a specific temperature, typically around 70°C, after the initial template denaturation. The reason is to allow the manual addition of unmodified **Taq** DNA polymerase which may lose activity if added during the initial denaturation step. Heat-activated **Taq** or Hot Start enzymes do not require this step.

Q5 *What is the incremented time and temperature function for?*

A5 Incremented/decremented time and temperature are used to increase or decrease either the time or temperature incrementally over the number of cycles in a stage. Incrementation of extension time is used with ‘Long PCR’ which is when large template fragments are to be amplified (e.g. 27kb lambda DNA, 40kb genomic DNA). Decremental temperature is used for protocols such as ‘Touchdown PCR’ where the first cycle starts with a high annealing temperature and over the number of cycles in the stage there is a gradual decrease in the temperature. This ensures that only the specific product is amplified.

Q6 *What is a gradient thermal cycler?*

A6 Gradient blocks enable a particular temperature step in a protocol to be programmed so that the temperature varies across the block. By specifying a temperature and the gradient to be applied, each column will achieve a different temperature. The set temperature is in the middle of the block, the lowest on the left and the highest on the right. The gradient is the total difference across the block, for example if the set temperature was 60°C and the gradient 10°C then the temperatures would range from approximately 55°C to 65°C from left to right.

Q7 *Why is a temperature gradient required?*

A7 The annealing temperature of the specific primers used in DNA amplification often requires optimisation. Instead of running multiple experiments with different annealing temperatures the gradient thermal cycler (PrimeG) allows the testing of up to 12 different annealing temperatures simultaneously.

TECHNICAL SUPPORT AND SERVICING

If you require further technical or application assistance please contact Techne at:

E-mail: technehelp@bibby-scientific.com

Phone: +44 (0)1785 810433

Fax: +44 (0)1785 810405

For servicing information please contact:

Service Department

Bibby Scientific Ltd.

Beacon Road

Stone

Staffordshire

ST15 0SA

E-mail: service@bibby-scientific.com

Phone: +44 (0)1785 810475

Fax: +44 (0)1785 810405

We are continually striving to improve our products and software. If you have any comments and suggestions on how we can do things better please send them to us at: techne@bibby-scientific.com

ADDITIONAL INFORMATION

USER MAINTENANCE

Before cleaning your unit, disconnect it from the power supply. The outer case of the Prime may be cleaned with a cloth dipped in water or ethanol (hexane or 50% methanol can also be used). No part of the case or cover should be immersed in the solvents. Do not use aggressive solvents such as acetone or abrasive cleaners. The block may be wiped with water, ethanol or propan-2-ol and may be decontaminated by wiping with 2% Neutracon or 1% bleach solution.

Before using any cleaning or decontamination method except those recommended here, the responsible person should check with Techne that the proposed method will not damage the equipment.

FAULT FINDING

Note that this equipment should only be dismantled by properly trained personnel. Removing the outer cover exposes potentially lethal mains voltages. There are no user serviceable parts within this equipment.

Should you have any problems with your Prime which cannot be easily remedied, you should contact your supplier and return the unit if necessary. Please include details of the fault observed and remember to return the unit in its original packing. Bibby Scientific Ltd. accepts no responsibility for damage to units which are not properly packed for shipping: if in doubt, contact your supplier, giving the full serial number of the unit.

FUSES

If the display on the front panel is not lit, one of the two fuses may have blown. Check that there is no external cause, such as a faulty plug or lead. Check both fuses and replace the faulty fuse with a new one of the correct value (fuse values are given on the label next to the power inlet). Note that fuses should only be replaced by a qualified electrician.

The holder for the two fuses is built into the mains input socket. First remove the power cable and then gently prise the fuse drawer open with a flat-bladed screwdriver or similar tool. Each fuse can be removed by using the screwdriver as a lever.

Exchange the faulty fuse in the fuse holder for a working fuse of the correct value. Finally, replace the fuse drawer in the fuse compartment and push the drawer shut. Fuses which blow repeatedly indicate a serious fault and you should return the unit to your supplier for repair.

INSULATION TESTING

This equipment is fitted with RFI suppression circuitry. Any check of the electrical insulation by means of high voltage dielectric testing (for example as in BS EN 61010-1) must be carried out using only a DC voltage.

This unit contains semiconductor components which may be damaged by electric field effects.

ACCESSORIES

The following accessories can be obtained from Bibby Scientific Ltd. or your Techne distributor:

Product code	Description
5PRIME/02/B	Prime block, 96 x 0.2ml
5PRIME/05/B	Prime block, 60 x 0.5ml
5PRIME/384/B	Prime block, 384 well plates
5PRIMEX/USB	Prime base gradient upgrade

REPLACEMENT PARTS

The following replacement parts can be obtained from Bibby Scientific Ltd. or your Techne distributor:

Product code	Description
HH179(S)	UK 230V mains lead with plug
HH180(S)	European mains lead with plug
FCABLEUS	US mains lead with plug

UK

Bibby Scientific Ltd.
Beacon Road, Stone
Staffordshire ST15 0SA
United Kingdom
Tel: +44 (0)1785 812121
Fax: +44 (0)1785 810405
e-mail: sales@bibby-scientific.com
www.bibby-scientific.com

France

Bibby Scientific Limited
Bâtiment Le Deltaparc Parc Silic PN2
7 rue du Canal
BP 55437 VILLEPINTE
95944 ROISSY Charles de Gaulle
France
Tel: +33(0)148 63 78 03
Fax: +33(0)148 63 78 01
e-mail: ventes@bibby-scientific.com
www.bibby-scientific.com

North & South America

Bibby Scientific US Inc.
t/a Techne Inc.
3 Terri Lane, Suite 10
Burlington, NJ 08016 USA
Toll Free (in NA): 800-225-9243
Tel: +1 609 589 2560
Fax: +1 609-589-2571
e-mail: labproducts@techneusa.com
www.techneusa.com

Italy

Bibby Scientific Italia Srl
Via Alcide de Gasperi 56
20077 Riozzo di Cerro al Lambro
Milano Italia
Tel: +39 (0)2 98230679
Fax: +39 (0)2 98230211
e-mail: marketing@bibby-scientific.it
www.bibby-scientific.it

Middle East

Bibby Scientific Middle East Ltd.
PO Box 27842, Engomi 2433
Nicosia
Cyprus
Tel: +357 22 660 423
Fax: +357 22 660 424
e-mail: sales@bibby-scientificme.com

Asia

Bibby Scientific - Singapore
Prudential Tower, Level 26
30 Cecil Street
Singapore 049712
Tel: +65 6631 2976
Fax: +44 (0)1785 810405
e-mail: info@bibby-scientific.com
www.bibby-scientific.com